

Technical Bulletin

Organisms Inactivated by Gamma Irradiation at 25 - 35 kGy

(Summarized from in-depth validation performed by SAFC Biosciences)

Organism	Туре	Size	Log Reduction
E.coli Escherichia coli	Gram Negative rod	1.5 μm - 4.0 μm	> 6
B.pumilus Bacillus pumilus	Gram Positive rod Spore former	0.6 μm - 2.5 μm	> 6
C.albicans Candida albicans	Fungus (dimorphic)	2 µm - 3 µm	> 6
A.laidlawii Acholeplasma laidlawii	Mycoplasma	0.3 μm - 0.8 μm	> 8
JH Strauss	Bacteriophage	24 nm - 26 nm	1
BVD Bovine Viral Diarrhea (cytopathic)	Togavirus (NADL - strain) ssRNA enveloped	30 nm - 60 nm	6
PI₃ Parainfluenza Type 3	Paramyxovirus ssRNA enveloped	100 nm - 250 nm	6
IBR Infectious Bovine Rhinotracheitis	Herpes dsDNA enveloped	100 nm	5
BTV Bluetongue Virus	Orbivirus dsRNA	60 nm - 80 nm	> 3
FELV Feline Leukemia Virus	Retrovirus RNA	100 nm - 120 nm	> 3
MVM Minute Virus of Mice	Parvovirus ssDNA nonenveloped	25 nm	2
PPV Porcine Parvovirus	Parvovirus ssDNA nonenveloped	18 nm - 25 nm	1

United States

 SAFC Biosciences, Inc.

 13804 W. 107th Street

 Lenexa, Kansas 66215

 USA

 Phone
 +1 913-469-5580

 Toll free-USA
 1 800-255-6032

 Fax
 +1 913-469-5584

 E-mail
 info-na@sial.com

Europe SAFC Biosciences Ltd. Smeaton Road, West Portway Andover, Hampshire SP10 3LF UNITED KINCDOM Phone +44 (0)1264-333311 Fax +44 (0)1264-332412 E-mail info-eu@sial.com

 SAFC Biosciences Pty. Ltd.

 18-20 Export Drive

 Brooklyn, Victoria 3025

 AUSTRALIA

 Phone
 +61 (0)3-9362-4500

 Toll free-AUS
 1 800-200-404

 Fax
 +61 (0)3-9315-1656

 E-mail
 info-ap@sial.com

Asia Pacific

www.safcbiosciences.com

Warranty, Limitation of Remedies

Varranty, Limitation of Remeales SAFC Biosciences warrants to the purchaser for a period of one year from date of delivery that this product conforms to its specifications. Other terms and conditions of this warranty are contained in SAFC Biosciences' written warranty, a copy of which is available upon request. ALL OTHER WARRANTIES, EXPRESSED OR IMPLIED, INCLUDING THE IMPLIED WARRANTY OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE, ARE EXCLUDED. In no case will SAFC Biosciences be liable for any special, incidental, or consequential damages arising out of this product or the use of this product by the customer or any third party based upon breach of warranty, breach of contract, negligence, strict tort, or any other legal theory. SAFC Biosciences expressly disclaims any warranty against claims by any third party by way of infringement or the like. THIS PRODUCT IS INTENDED FOR PURPOSES DESCRIBED ONLY AND IS NOT INTENDED FOR ANY HUMAN OR THERAPELITY. ANY HUMAN OR THERAPEUTIC USE.

Additional Terms and Conditions are contained in the product Catalog, a copy of which is available upon request.

© 2006 SAFC Biosciences, Inc.

Issued April 2006 T022 0103 1205

www.safcbiosciences.com

United States

SAFC Biosciences, Inc. 13804 W. 107th Street Lenexa, Kansas 66215 USA Phone +1 913-469-5580 1 800-255-6032 +1 913-469-5584 Toll free-USA Fax E-mail info-na@sial.com

Europe

SAFC Biosciences Ltd. Smeaton Road, West Portway Andover, Hampshire SP10 3LF UNITED KINGDOM Phone +44 (0)1264-333311 +44 (0)1264-332412 info-eu@sial.com Fax E-mail

Asia Pacific

SAFC Biosciences Pty. Ltd. 18-20 Export Drive Brooklyn, Victoria 3025 AUSTRALIA +61 (0)3-9362-4500 Phone 1 800-200-404 +61 (0)3-9315-1656 info-ap@sial.com Toll free-AUS Fax E-mail