

1.00801.0001

# Label Set

## Adhesive labels 80 x 110 mm

### for self labelling

Contents:

20 adhesive labels 80 x 110 mm

168 adhesive pictograms

40 adhesive signal words

Package insert with H, P and EUH statements

The labels and pictograms of this package offer you the option to label your own individual mixtures and preparations in analogy with the GHS regulations:

The upper white field of the basic label provides a space for the designation of the contents. The two following lines serve the purpose of clearly stating the ingredients with a hazard potential and of listing all the components. The four pictogram fields can be used to affix the rhomboid pictograms included with the package. The pictograms must be affixed standing on one corner.





The signal words "Caution" and "Danger" point out the severity of the respective hazard. The signal word "Danger" is for use in connection with serious hazard categories, while the word "Caution" is used for less serious hazard categories. Please use the enclosed self-adhesive signal word labels for the empty field in the red box in the lower left space of the basic label. In the case that local regulations require other signal words in the respective language, these can be entered directly in this field. A selection of the official wordings in different languages is also enclosed.

The lower right space is used for the entry of further details as per the CLP guidelines and DIN EN ISO.

The enclosed leaflet contains the currently valid hazard statements, supplemental hazard information, precautionary statements (H, EUH, P) with the corresponding codes.

If you have any questions regarding classification and labeling, please contact us at MSDS-EU@merckgroup.com.

Specimen completed and affixed label:

<b>MOBILE PHASE A</b>			
Enthält: / Contains: <b>Acetonitril (60%), Water (10%)</b>			
<b>TFA (0.1%)</b>			
			
<b>Gefahr / Danger</b> <b>Peligro</b>	Hergestellt am / Prepared on <b>2021/03/22</b>	mind. verwend. bis / min. shelf life <b>2021/03/26</b>	
	<b>A. Schneider, Lab 001, Tel. 1234</b>		
	Hergestellt von / Prepared by		

We would like to point out that you are fully responsible for complying with existing laws and regulations. The above / following information does not release you from your sole responsibility to classify, label and pack your preparations and mixtures according to the specifications of the CLP/EU-GHS (Global Harmonized System) directive. By supplying this information we neither assume liability for the accuracy of such information nor for the accuracy or legal conformity of the classification, labeling or packaging.

#### Signal words

BG	FR	NL
Опасно	Danger	Gevaar
Внимание	Attention	Waarschuwing
ES	GA	PL
Peligro	Contúirt	Niebezpieczeństwo
Atención	Rabhadh	Uwaga
CS	HR	PT
Nebezpečí	Opasnost	Perigo
Varování	Upozorenje	Atenção
DA	IT	RO
Fare	Pericolo	Pericol
Advarsel	Attenzione	Atenție
DE	LV	SK
Gefahr	Bīstami	Nebezpečnost
Achtung	Uzmanību	Pozor
ET	LT	SL
Ettevaatus	Pavojinga	Nevarno
Hoiatus	Atsargai	Pozor
EL	HU	FI
Κίνδυνος	Veszély	Vaara
Προσοχή	Figyelem	Varoitus
EN	MT	SV
Danger	Periklu	Fara
Warning	Twissija	Varning

Version April 2021

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

© 2021 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved. The vibrant M is a trademark of Merck KGaA, Darmstadt, Germany. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly available resources.

Merck KGaA, 64271 Darmstadt, Germany, Tel. +49(0)6151 72-2440  
EMD Millipore Corporation, 400 Summit Drive, Burlington MA 01803, USA,  
Tel. +1-978-715-4321

Sigma-Aldrich Canada Co. or Millipore (Canada) Ltd. 2149 Winston Park,  
Dr. Oakville, Ontario, L6H 6J8, Phone: +1 800-565-1400

