

74807 Blood Free Campylobacter Medium Selective Supplement

An antimicrobial supplement recommended for the selective isolation of Campylobacter species

Composition

(per vial, sufficient for 500 ml medium):

Cefoperazone 16 mg

Directions

Rehydrate the contents of 1 vial aseptically with 2 ml of sterile distilled water. Mix well and aseptically add to 500 ml of sterile, molten Blood Free Campylobacter Selectivity Agar Base (B2426). Mix well and pour into sterile petri plates.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

