

Product Information

Nuclease micrococcal from *Staphylococcus aureus*

With reaction buffer and BSA solution

CS0004

Product Description

CAS Number: 9013-53-0

Enzyme Commission (EC) Number: 3.1.31.1

Synonyms: Endonuclease micrococcal, MNase, Micrococcal endonuclease

Micrococcal Nuclease (MNase) is an endo-exonuclease that preferentially digests single-stranded nucleic acids. MNase was historically isolated from *Micrococcus pyogenes* var. *aureus* (now known as *Staphylococcus aureus*).^{1,2} The end products from MNase activity are mononucleotides and oligonucleotides with terminal 3'-phosphates. MNase is also active against double-stranded DNA and RNA, and all nucleic acids will be ultimately cleaved. MNase can be used for the limited digestion of chromatin to identify the location of nucleosomes.³

Applications of MNase include:

- Chromatin Immuno-Precipitation (ChIP) analysis⁴
- Nucleosome preparation
- Removal of nucleic acids present in protein preparation
- Reduction of viscosity of cell lysates during non-mechanical cell lysis preparation
- RNA sequencing and transcriptomic profiling

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Components

- Nuclease Micrococcal (Component CS0004A): 50 units, in solution [300-500 units/mL in 10 mM Tris (pH 7.5), 50 mM NaCl, 1 mM EDTA, and 50% Glycerol]
- MNase Reaction Buffer, 10× (Component CS0004B): 1 mL [solution of 500 mM Tris (pH 8.0), with 50 mM CaCl₂]
- BSA (Component CS0004C): 1 mL (solution at 1 mg/mL BSA in water)

Storage/Stability

Store the kit components at -20 °C.

Preparation Instructions

For optimal activity, it is recommended to perform the reaction in 1× MNase reaction buffer with 0.1 mg/mL BSA. If dilution of enzyme is required:

1. Prepare fresh Diluent Solution (1× MNase reaction buffer with 0.1 mg/mL BSA solution). Store this Diluent Solution on ice.
2. Right before use, dilute the enzyme with the ice-cold Diluent Solution.

Procedure

- Nucleosome preparation can be performed as part of a ChIP procedure.
- Nucleosome fragmentation can be facilitated by a combination of sonication and digestion with MNase.

A generic procedure for nucleosome preparation is presented here. For specific materials and methodology, various publications in the literature may be consulted.⁵⁻⁸

Cell culture preparation and fixation


It is recommended to use this procedure with 4×10^6 to 5×10^6 cells. This assay was tested with HeLa cells and might require a calibration when other cells are used.

1. Crosslink cells:
 - 1.1. Add 1% formaldehyde for 10 minutes.
 - 1.2. Add 125 mM glycine for 2 minutes.
2. Remove the cell culture medium.
 - 2.1. Collect the cells.
 - 2.2. Centrifuge cells at $2000 \times g$ at 4°C for 5 minutes.
 - 2.3. Discard the supernatant.
 - 2.4. Continue to the next stage.

Nucleosome preparation

- Perform the MNase digest in a buffer composed of $1 \times$ MNase Reaction Buffer (Component CS0004B) with 0.1 mg/mL BSA (Component CS0004C).
 - If another buffer is used, make sure that it contains 1-5 mM CaCl_2 .
1. Wash the cells according to the ChIP protocol of choice. Transfer the cells to CaCl_2 -containing buffer.
 2. Add 0.5 μL of Nuclease Micrococcal (MNase, Component CS0004A). Mix well. Incubate for 20 minutes at 37°C .
 3. Add 10 μL of 0.5 M EDTA to stop MNase activity. Incubate on ice for 2 minutes.
 4. Centrifuge at $12,000 \times g$ for 1 minute at 4°C . Discard the supernatant.
 5. Resuspend the pellet in 100-500 μL of buffer as required in microcentrifuge tube.
 6. Sonicate the resuspended pellet with a probe sonicator, to break the nuclear membrane. This stage must be calibrated according to specific assay conditions. As a starting point, perform three sonication pulses, 10 seconds each, with 10-second intervals on ice. This was found to produce a homogenous preparation of monomeric nucleosomes (see Figure 1).

Figure 1. Chromatin preparation analysis

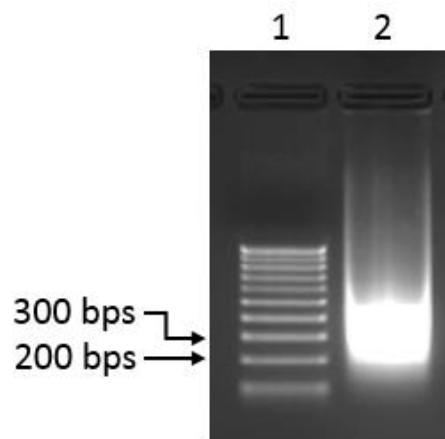

HeLa cells were formaldehyde-crosslinked.

Chromatin was prepared and digested as described:

- A 50 μL sample of DNA was incubated with RNase and Proteinase K.
- The sample was purified using a microspin column (such as Catalogue Number NA1020), according to the column instructions.
- A sample of 10 μL was separated by electrophoresis on a 2% agarose gel (Lane 2) and stained with ethidium bromide. Lane 2 shows that the chromatin was digested to 1 to 2 nucleosomes in length.

References

1. Cunningham, L. *et al.*, *J. Am. Chem. Soc.*, **78(18)**, 4642-4645 (1956).
2. Telford, D.J., and Stewart, B.W., *Int. J. Biochem.*, **21(2)**, 127-138 (1989).
3. Chung, H.-R. *et al.*, *PLoS ONE*, **5(12)**, e15754 (2010).
4. Long, L. *et al.*, *J. Immunother. Cancer*, **9(12)**, e003973 (2021).
5. O'Neill, L.P., and Turner, B.M., *Methods Enzymol.*, **274**, 189-197 (1996).
6. Zeng, P.-Y. *et al.*, *BioTechniques*, **41(6)**, 694-698 (2006).
7. Roca, H., and Franceschi, R.T., *Nucleic Acids Res.*, **36(5)**, 1723-1730 (2008).
8. Cosseau, C., and Grunau, C., *Methods Mol. Biol.*, **791**, 195-212 (2011).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

MilliporeSigma, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.
© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

CS0004pis Rev 07/22 NB,DT,LB,GCY,MAM

**Millipore
Sigma**