

Product Information

Concanavalin A-Agarose

From *Canavalia Ensiformis* (Jack Bean)

C7555

Storage temperature 2-8 °C

Synonym: ConA.

Product Description

Concanavalin A (Con A) is a lectin isolated from the jack bean. At pH > 7 it exists as a tetramer.¹ This is the species most likely attached to the resin. Con A is not a glycoprotein.³ It does not contain cysteine residues.⁴

Unlike most other lectins, Con A is a metalloprotein and requires divalent cations, such as manganese and calcium ions for binding.² Each subunit of Con A binds one calcium ion and one manganese ion. Removal of these cations under acidic conditions abolishes the carbohydrate-binding activity.⁵

Con A binds specifically to mannosyl and glucosyl residues of polysaccharides and glycoproteins.⁵ Unmodified hydroxyl groups at the C3, C4, and C6 positions of D-glucopyranosyl or D-mannopyranosyl rings may be essential for binding.⁶

General adsorption and desorption procedures and methods to prevent leakage of the lectin from the matrix have been described.^{7,8,9} This product has been used to isolate human fibroblast interferon¹⁰ and glycoprotein from sarcoplasmic reticulum.¹¹

Con A agarose has been used with SDS (0.05%)¹² and TRITON® X-100.¹³

This product is prepared with cross-linked 4% beaded agarose and is supplied in 0.1 M potassium phosphate, pH 6, containing 1 M NaCl, 1 mM CaCl₂, 1 mM MnCl₂, 1 mM MgCl₂, and preservative.

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Procedure

Since Con A requires Mn²⁺ and Ca²⁺ ions for carbohydrate binding, buffers should either include these metal ions or the gel should be equilibrated with these ions, immediately prior to use. This is particularly important if the sample may contain traces of chelating agent or if the resin had been treated under acidic conditions (< pH 5).

The procedure should be conducted at 2-8 °C.

1. Pre-wash column with 5 column volumes of wash solution (1 M NaCl, 5 mM MgCl₂, 5 mM MnCl₂, and 5 mM CaCl₂).
2. Equilibrate column in the buffer of choice (pH ranges generally between 6.5 to 7.5 although buffers as low as pH 4.1 and as high as 9.0 have been used successfully). A commonly used starting buffer is 20 mM Tris, pH 7.4, containing 0.5 M NaCl.
3. Load sample solution in equilibration buffer (protein concentrations 1-20 mg/mL, free of particulates).
4. Wash the resin with equilibration buffer until eluent solution is protein free.

Elution

5. Elute the target protein with gradient or stepwise elution with methyl α-D-glucopyranoside or methyl α-D-mannopyranoside, glucose, or mannose (5 mM - 500 mM).
6. Maximum recovery and cleaning of the resin may be achieved by using 1 M sucrose, glucose, mannose, or corresponding α-methyl glycoside. The addition of chaotropic agents (0.5 M to 6 M) may also be required for maximum recovery, but these denaturing conditions may severely damage the resin. Therefore, they should be used last.

Regeneration

7. After use, wash the column with 5-10 column volumes of 0.1 M Borate, pH 8.5, or 0.1 M Tris, pH 8.5, containing 0.5 M NaCl.
8. This should be followed by 0.1 M sodium acetate, pH 4.5, containing 1.0 M NaCl.
9. The acid and basic washes may need to be cycled 2-3 times for complete washing.
10. Finally, the gel should be stored at 2-8 °C in 1.0 M NaCl containing 5 mM CaCl₂, 5 mM MnCl₂, and 5 mM MgCl₂ or in 0.1 M acetate buffer, pH 6, containing 1 M NaCl, 1 mM CaCl₂, 1 mM MgCl₂, with preservative.

References

1. Senear, D.F. and Teller, D.C., Thermodynamics of concanavalin A dimer-tetramer self-association: sedimentation equilibrium studies. *Biochemistry*, 20(11), 3076-3083 (1981).
2. Sumner, J.B. and Howell, S.F., *J. Biol. Chem.*, 115, 583 (1936).
3. Olson, M.O. and Liener, I.E., Some physical and chemical properties of concanavalin A, the phytohemagglutinin of the jack bean. *Biochemistry*, 6(1), 105-111 (1967).
4. Agarwal, B. and Goldstein, I.J., Protein-carbohydrate interaction. VII. Physical and chemical studies on concanavalin A, the hemagglutinin of the jack bean. *Arch. Biochem. Biophys.*, 124(1), 218-229 (1968).
5. Goldstein, I.J., et al., *Biochemistry*, 4, 876 (1965).
6. Yahara, I. and Edelman, G.M., Restriction of the mobility of lymphocyte immunoglobulin receptors by concanavalin A. *Proc. Nat. Acad. Sci. USA*, 69(3), 608-612 (1972).
7. Bittiger, H. and Schnebli, H.P., ed. *Affinity Chromatography on Immobilized Concanavalin A. In Concanavalin A as a Tool*. Wiley and Sons, p. 323-329 (1976).
8. Kowal, R. and Parsons, R.G., Stabilization of proteins immobilized on Sepharose from leakage by glutaraldehyde crosslinking. *Anal. Biochem.*, 102, 72-76 (1980).
9. Marikar, Y., et al., Leaching of concanavalin A during affinity chromatographic isolation of cell surface glycoproteins from human fetal neurons and glial cells. *Anal. Biochem.*, 201(2), 306-310 (1992).
10. Davey, M.W., et al., Binding of human fibroblast interferon to concanavalin A-agarose. Involvement of carbohydrate recognition and hydrophobic interaction. *Biochemistry*, 15(3), 704-713 (1976).
11. Campbell, K.P. and MacLennan, D.H., Purification and characterization of the 53,000-dalton glycoprotein from the sarcoplasmic reticulum. *J. Biol. Chem.*, 256(9), 4626-4632 (1981).
12. Li, Y., et al., The p185neu-containing Glycoprotein Complex of a Microfilament-associated Signal Transduction Particle. *J. Biol. Chem.*, 274(36), 25651-25658 (1999).
13. Romero, P.A. et al., Glycoprotein biosynthesis in *Saccharomyces cerevisiae*. Partial purification of the alpha-1,6-mannosyltransferase that initiates outer chain synthesis. *Glycobiology*, 4(2), 135-140 (1994).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice).

Standard Warranty

The applicable warranty for the products listed in this publication may be found at [SigmaAldrich.com/terms](https://www.sigmaaldrich.com/terms).

Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://www.sigmaaldrich.com/offices).

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.