



---

## Product Information

### 5-HYDROXYEICOSATETRANOIC ACIDS and Derivatives

Product Nos. H 1017, H 6643, H 3894, H 4894,  
H 4144, H 2521, H 4269, H 7768, H 1515, H 1142,  
H 3023, H 7147, H 0270, H 2762, H 2645, and H 0153

#### Special Handling Instructions

All glassware and stainless steel syringes used for transferring HETE solutions should be cleaned, avoiding activating solvents such as strong acid, and then rinsed using absolute ethanol and air dried. Transfer and removal of samples should be performed in indirect light and under an inert atmosphere, with the bulk kept below 10 °C. Return bulk to freezer storage as soon as possible. Hydroxyeicosatetraenoic acid is unstable to oxygen and direct sunlight, therefore, extreme care should be taken in handling this product.

PCS/KMR 01/02

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.