

## Product Information

## 3CL, Mpro protease chromogenic substrate peptide

SARS-CoV2 Main protease chromogenic substrate peptide, pNA-peptide substrate, lyophilized

**SAE0181**

### Product Description

This hexapeptide, Thr-Ser-Ala-Val-Leu-Gln-pNA, is a substrate peptide for both the Main protease (Mpro) from the SARS coronavirus,<sup>1,2</sup> and also the Mpro of SARS-CoV2.<sup>3</sup> This substrate allows measurement of the activity of Mpro, using a standard plate reader or spectrophotometer. Proteolytic cleavage by Mpro will release the *p*-nitroaniline (pNA) group from the peptide (TSAVLQ), with a maximum absorption peak at 405 nm ( $A_{405}$  of 0.00916 at 1  $\mu$ M):



### Reagent

This product is lyophilized from 0.1% TFA in H<sub>2</sub>O.

### Storage/Stability

The product is stable in lyophilized form for at least 5 years when stored at -20 °C.

The product has a limited lifetime in solution. Long-term storage in solution should be avoided.

### Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

### Preparation Instructions

#### Solubilization:

The peptide is soluble in water at a concentration of 20 mg/mL. Alternatively, the peptide is soluble in DMSO at a concentration of 20-50 mg/mL.

### Procedure

#### Working reagent concentrations:

**Peptide:** suggested working concentration = 200  $\mu$ g/mL

**Mpro:** we recommend using Mpro, Cat. No. SAE0172, at a working concentration of 2-20  $\mu$ g/mL.

**Assay buffer:** Mpro protease is active under a wide variety of conditions and temperatures. However, we strongly advise against use of Tris-based buffers, because of their interactions with the given substrate.

Recommended assay buffer: 25 mM HEPES buffer, pH 7.0, 0.2% TWEEN® 20

**Reaction temperature:** 0-30 °C

The reaction can be stopped by addition of acetic acid to a final concentration of 2%.

### References

1. Liu, P. *et al.*, Potent inhibitors of SARS-CoV-2 3C-like protease derived from N-substituted isatin compounds. *Eur. J. Med. Chem.*, **206**, 112702 (2020).
2. Liu, Z. *et al.*, Virtual screening of novel noncovalent inhibitors for SARS-CoV 3C-like proteinase. *J. Chem. Inf. Model.*, **45**(1), 10-17 (2005).
3. Li, C. *et al.*, Maturation mechanism of severe acute respiratory syndrome (SARS) coronavirus 3C-like proteinase. *J. Biol. Chem.*, **285**(36), 28134-28140 (2010).

---

## Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

### Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://SigmaAldrich.com/techservice).

### Standard Warranty

The applicable warranty for the products listed in this publication may be found at [SigmaAldrich.com/terms](https://SigmaAldrich.com/terms).

### Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://SigmaAldrich.com/offices).

The life science business of Merck operates  
as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

SAE0181pis Rev 02/22 DK,DT,GCY

