

Product Information

5-Bromo-4-chloro-3-indolyl phosphate *p*-toluidine salt

≥99% (HPLC)

B8503

Product Description


CAS Registry Number: 6578-06-9

Molecular Formula: C₈H₆BrClNO₄P · C₇H₉N

Formula Weight: 433.62

Synonyms: BCIP® *p*-toluidine salt, X-phosphate *p*-toluidine salt

Storage temperature: -20 °C

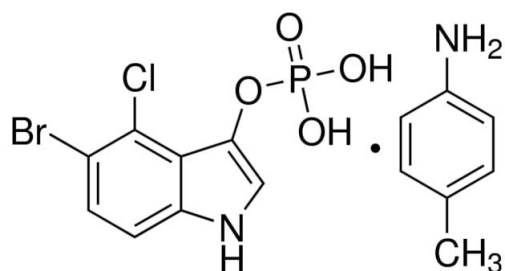

5-Bromo-4-chloro-3-indolyl phosphate (BCIP®) and nitro blue tetrazolium (NBT) are commonly used in tandem for the colorimetric detection of alkaline phosphatase-labeled molecules.¹⁻³ The BCIP®/NBT substrate system is versatile and functions in a variety of applications, including:

- Northern, Southern, and Western blotting^{4,5}
- *in situ* hybridization⁶
- immunohistochemistry⁷

BCIP® *p*-toluidine salt is soluble in dimethylformamide (DMF). DMF may be used to prepare a stock solution. A portion of the BCIP® stock solution is then combined with NBT in a reaction buffer to form a substrate solution for alkaline phosphatase. This substrate system, when incubated with alkaline phosphatase, produces an insoluble NBT diformazan product⁸ that is easily observable with its purple color. (See Figure 1 for a reaction scheme.)

BCIP® is prepared synthetically. Several theses⁹⁻¹³ and dissertations¹⁴⁻²⁸ have cited use of product B8503 in their protocols.

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Solubility

BCIP® *p*-toluidine salt is tested for solubility in dimethylformamide (DMF) at 20 mg/mL.

Storage/Stability

Store BCIP® lyophilized product at -20 °C, protected from light and moisture.


One publication has reported that 50 mg/mL stock solutions of BCIP in 100% DMF can be stored at 4 °C.²⁹ A BCIP® stock solution in DMF remains active for ~2 weeks kept in the dark at 2-8 °C. However, a working solution in aqueous buffer is only good for one day. Another publication has reported that 50 mg/mL stock solutions of BCIP in 100% DMF can be stored at -20 °C.³⁰

References

1. Blake, M.S. *et al.*, *Anal. Biochem.*, **136(1)**, 175-179 (1984).
2. Green, F.J. (ed.), *The Sigma-Aldrich Handbook of Stains, Dyes & Indicators*, Aldrich Chemical Co., (Milwaukee, WI), p. 523 (1990).
3. Horowitz, J.P. *et al.*, *J. Med. Chem.*, **9(3)**, 447 (1966).
4. Knecht, D.A., and Dimond, R.L., *Anal. Biochem.*, **136(1)**, 180-184 (1984).
5. Leary, J.J., *et al.*, *Proc. Natl. Acad. Sci. USA*, **80(13)**, 4045-4049 (1983).

6. Walters, C. *et al.*, *J. Clin. Pathol.*, **50(9)**, 749-754 (1997).
7. Meltzer, J.C. *et al.*, *J. Histochem. Cytochem.*, **45(4)**, 599-610 (1997).
8. McGadey, J., *Histochemie*, **23(2)**, 180-184 (1970).
9. Kubanek, Carol-Anne, "Antigenic analysis of cowpea severe mosaic virus and other selected viruses using monoclonal antibodies". Iowa State University, M.S. thesis, p. 66 (1987).
10. Colicos, Michael Adrian, "Construction of an adenovirus expression vector containing the T4 den V gene, which can complement the DNA repair deficiency of *Xeroderma pigmentosum* fibroblasts". McMaster University, M.Sc. thesis, p. 69 (1988).
11. Pursey, Jane Constance, "Characterization of Cytochrome P450 and a Putative Cytochrome P450 Gene in *Drosophila melanogaster*". McMaster University, M.Sc. thesis, p. 28 (1989).
12. Martin, Emily Ann Kean, "*Escherichia coli* in Tom Turkeys and Their Environment". University of Guelph, M.Sc. thesis, p. 129 (2000).
13. Nur Sami, Israk, "'O antigenic' Polysaccharide Specific Memory B Cell Responses in Bangladeshi Cholera Patients Infected with *Vibrio cholerae* O1". BRAC University (Bangladesh), M.Sc. thesis, p. 28 (2015).
14. Bos, Nicolaas Adrianus, "The Influence of Exogenous Antigenic Stimulation on the Murine 8 Cell Repertoire". Erasmus Universiteit Rotterdam, Ph.D. dissertation, p. 43 (1990).
15. Jonusys, Alison Michéle, "Immunological Characterisation of Erythrocyte Antigens". University of Wollongong, Ph.D. dissertation, p. 68 (1991).
16. Griffin, Jennifer K., "An analysis of tea domain proteins". University of Kansas, Ph.D. dissertation, p. 48 (1999).
17. Singleton, Richard Thomas, "B cell responses to respiratory viral infections". University College London, Ph.D. dissertation, p. 104 (2004).
18. Karreman, Robert Jan, "The Localization, Function and Applications of the Stress Response Protein Hsp12p in the yeast *Saccharomyces cerevisiae*". University of Cape Town, Ph.D. dissertation, p. 31 (2006).
19. Revilla-i-Domingo, Roger, "*cis*-Regulatory analysis of the sea urchin *delta* gene: validating the architecture of the gene regulatory network model for micromere lineage specification". California Institute of Technology, Ph.D. dissertation, p. 94 (2007).
20. Eckrich, Tobias, "Voltage gated Na⁺ channels and spontaneous action potential activity in cochlear hair cells during development". University of Sheffield, Ph.D. dissertation, p. 81 (2012).
21. Heimer-McGinn, Victoria, "Development and characterization of novel transgenic techniques for the study of neuronal circuitry". National University of Ireland (University College Cork), Ph.D. dissertation, p. 128 (2013).
22. Applebaum, Mordechai, "A Molecular Network Underlying Muscle Lineage Segregation from the Embryonic Dermomyotome". The Hebrew University of Jerusalem, Ph.D. dissertation, p. 27 (2014).
23. Fleming, Margaret Brigham, "The influence of extensin cross-linking on biomass recalcitrance". Colorado State University, Ph.D. dissertation, pp. 11, 93 (2015).
24. Filatova, Anna, "The role of *Fam20a* in the generation of Amelogenesis Imperfecta". Universität Zürich, Dr. sc. nat. dissertation, p. 32 (2016).
25. Zahabi, Azadeh, "Modeling of hematopoiesis using ES and patients-derived iPSCs". Eberhard-Karls-Universität Tübingen, Dr. rer. nat. dissertation, p. 41 (2016).
26. Mohamed, Adil, "Host and Viral Determinants of Reovirus Oncolysis". University of Alberta, Ph.D. dissertation, p. 56 (2018).
27. Jin, Ye, "The Role of Pitx Proteins in Early *Xenopus* Development". City University of New York, Ph.D. dissertation, p. 19 (2017).
28. Doddamani, Dadakhalandar, "The transcriptome analysis of Primordial Germ Cells of birds ". University of Edinburgh, Ph.D. dissertation, p. 46 (2019).
29. Hong, Y., and Scharl, M., *Methods Mol. Biol.*, **329**, 3-16 (2006).
30. Byrne, C., and Hardman, M.J., *Methods Mol. Biol.*, **289**, 127-136 (2005).

Figure 1. BCIP®/NBT Reactions

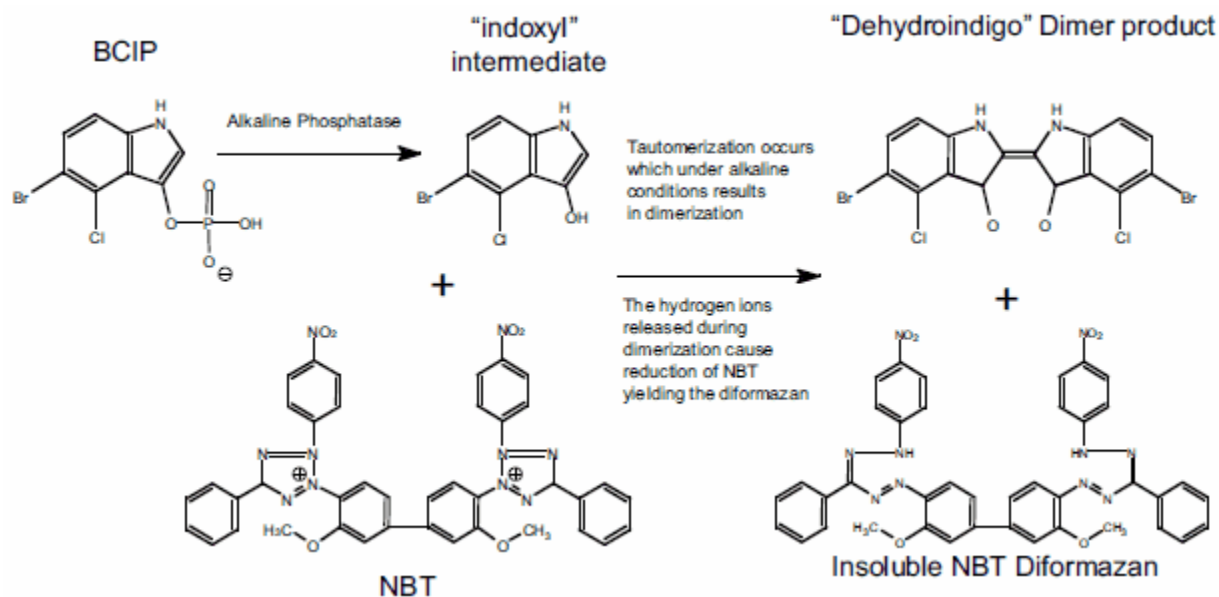

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice).

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at [SigmaAldrich.com/terms](https://www.sigmaaldrich.com/terms).

Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://www.sigmaaldrich.com/offices).

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

MilliporeSigma, BCIP, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2009-2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

B8503pis Rev 08/22 RBG,KTA,GCY,MAM

**MILLIPORE
SIGMA**