

59886 Anaerob Indicator Test

Anaerob Indicator Test is a sachet containing a test strip saturated with resazurin solution. It is for checking the anaerobic atmosphere in sealed containers used for incubation. The color change from pink to white indicates anaerob conditions.

Contents:

100 Test Strips per package

(Test Strips saturated with resazurin in a buffered aqueous solution containing preservatives)

Storage:

Store at 2°C to 8°C in the sachets provided. Allow to equilibrate to room temperature before opening.

Directions:

Cut as indicated. Using forceps, withdraw approximatly 10 mm of strip. Place in the anaerobic jar immediatly before applying the lid.

If after incubation, the indicatorstrip has not changed color, the integrity of the jar seals and catalyst function must be checked and the anaerobic culture repeated.

After use the sachet can be discarded with normal laboratory litter.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

