


Ascentis Express F5 HPLC Columns

Fused-Core® Technology Columns

Ascentis Express columns provide a breakthrough in HPLC column performance, as well as the benefits of sub 2 micron particles at much lower backpressure. Ascentis Express F5 brings together the highly efficient, robust 2.7 µm Fused-Core particle technology with F5 chemistry for the successful separation of polar and non-polar organic compounds.

The pentafluorophenylpropyl (PFP) stationary phase of Ascentis Express F5 provides a stable, reversed-phase packing with electron-deficient phenyl rings due to the presence of electronegative fluorines. In addition to forming π - π and mildly steric interactions, F5 phases also retain compounds by polar interactions. As a result of having both polar and non-polar character, F5 phases can show dual-mode retention behavior, sometimes producing a "U-shaped" retention as a function of acetonitrile content of the mobile phase, with retention increasing at both low and high concentrations of ACN (reversed-phase and HILIC retention modes). Ascentis Express F5 can be used for basic, acidic, or neutral compounds with alternate selectivity from C18.

Fused-Core Structure of Ascentis Express

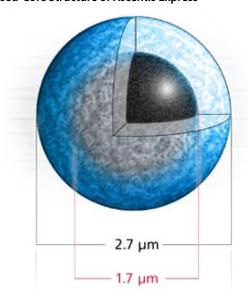

For more information on Ascentis Express HPLC Columns, call our technical experts at 800-359-3041/814-359-3041 or visit sigma-aldrich.com/express

Key Benefits of Ascentis Express

- Double the efficiency of 3 micron columns
- Same efficiency of sub 2 micron columns at half of the backpressure
- Rugged design capable of ultra high pressure operation

Key Benefits of F5

- Retains bases more and hydrophobic compounds less than C18
- Reversed phase, HILIC, and 100% aqueous applications
- Ideal for the separation of closely related compounds

Comparison of C18 and F5 for the Separation of Closely Related Steroids


columns: Ascentis Express F5 (53590-U) and Ascentis Express C18 (53827-U),

 $10 \text{ cm} \times 4.6 \text{ mm}$, I.D., $2.7 \mu \text{m}$ particle size

mobile phase A: water mobile phase B: methanol mobile phase ratio: A:B 50:50, v/v flow rate: 0.8 mL/min

flow rate: 0.8 mL/min temp.: 35 °C pressure: ~2400 psi det.: UV at 240 nm ini: 5 uL

sample: $10 \,\mu g/mL$ each in 90:10 water:methanol

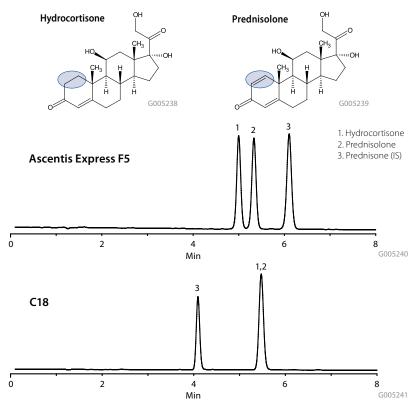

Ordering Information

Analytical Ascentis Express Columns

ID (mm)	Length (cm)	C18	C8	Phenyl-Hexyl	HILIC	RP-Amide	F5
2.1	2	53799-U	53795-U	53798-U	_	53797-U	53592-U
2.1	3	53802-U	53839-U	53332-U	53933-U	53910-U	53566-U
2.1	5	53822-U	53831-U	53334-U	53934-U	53911-U	53567-U
2.1	7.5	53804-U	53843-U	53335-U	53938-U	53912-U	53568-U
2.1	10	53823-U	53832-U	53336-U	53939-U	53913-U	53569-U
2.1	15	53825-U	53834-U	53338-U	53946-U	53914-U	53571-U
3.0	3	53805-U	53844-U	53341-U	53964-U	53915-U	53574-U
3.0	5	53811-U	53848-U	53342-U	53967-U	53916-U	53576-U
3.0	7.5	53812-U	53849-U	53343-U	53969-U	53917-U	53577-U
3.0	10	53814-U	53852-U	53345-U	53970-U	53918-U	53578-U
3.0	15	53816-U	53853-U	53346-U	53972-U	53919-U	53579-U
4.6	3	53818-U	53857-U	53347-U	53974-U	53921-U	53581-U
4.6	5	53826-U	53836-U	53348-U	53975-U	53922-U	53583-U
4.6	7.5	53819-U	53858-U	53351-U	53977-U	53923-U	53584-U
4.6	10	53827-U	53837-U	53352-U	53979-U	53929-U	53590-U
4.6	15	53829-U	53838-U	53353-U	53981-U	53931-U	53591-U

Capillary Ascentis Express Columns

	•	
Length (cm)	I.D. (μm)	Cat. No.
C18		
5	75	53982-U
15	75	54219-U
5	100	53985-U
15	100	54256-U
5	200	53989-U
15	200	54261-U
5	300	53992-U
15	300	54271-U
5	500	53998-U
15	500	54273-U
C8		
5	75	53983-U
15	75	54229-U
5	100	53987-U
15	100	54260-U
5	200	53991-U
15	200	54262-U
5	300	53997-U
15	300	54272-U
5	500	53999-U
15	500	54275-U

Trademarks

Ascentis – Sigma-Aldrich Biotechnology LP Fused-Core – Advanced Materials Technology Inc.

Ascentis Express Guard Columns



Universal Guard Holder

Description	Cat. No.
Holder w/EXP Titanium Hybrid Ferrule	53500-U
cartridge not included with holder	53500-0

Ascentis Express Guard Cartridges

Description	I.D. (mm)	Pkg. Size	Cat. No.
C18	2.1	3	53501-U
C18	3.0	3	53504-U
C18	4.6	3	53508-U
C8	2.1	3	53509-U
C8	3.0	3	53511-U
C8	4.6	3	53512-U
RP-Amide	2.1	3	53514-U
RP-Amide	3.0	3	53516-U
RP-Amide	4.6	3	53519-U
HILIC	2.1	3	53520-U
HILIC	3.0	3	53521-U
HILIC	4.6	3	53523-U
Phenyl-Hexyl	2.1	3	53524-U
Phenyl-Hexyl	3.0	3	53526-U
Phenyl-Hexyl	4.6	3	53531-U
F5	2.1	3	53594-U
F5	3.0	3	53597-U
F5	4.6	3	53599-U

Enabling Science to Improve the Quality of Life Order/Customer Service (800) 325-3010 • Fax (800) 325-5052 Technical Service (800) 325-5832 • sigma-aldrich.com/techservice Development/Custom Manufacturing Inquiries **SAFC*** (800) 244-1173 Safety-related Information sigma-aldrich.com/safetycenter World Headquarters 3050 Spruce St. St. Louis, MO 63103 (314) 771-5765 sigma-aldrich.com

©2010 Sigma-Aldrich Co. All rights reserved. SiGMA, SAFC, SiGMA-ALDRICH, ALDRICH, FLUKA, and SUPELCO are trademarks belonging to Sigma-Aldrich Co. and its affiliate Sigma-Aldrich Biotechnology, L.P. Sigma brand products are sold through Sigma-Aldrich, Inc. Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.

MSF T410095A