

Product Information

CM Sepharose® Fast Flow

CCF100

Product Description

CM Sepharose® Fast Flow is an ion exchange chromatography resin with a carboxymethyl (CM) functional group ($-\text{OCH}_2\text{COO}^-$) attached to a crosslinked 6% agarose matrix. The CM group serves as a weak cation exchanger, where the degree of dissociation, and thus the exchange capacity, varies markedly with pH. The terms "strong" and "weak" in ion exchange chromatography refer to the extent of ionization with pH, and **not** to the binding strength of the functional group to the target species.

The parent Sepharose® Fast Flow is a cross-linked derivative of Sepharose®. The particle size range is 45–165 μm . The average bead diameter is $\sim 90 \mu\text{m}$. The counterion in the product is sodium (Na^+).

Recommended anionic buffers to use with CM Sepharose® Fast Flow include acetate, barbiturate, citrate, glycine, or phosphate. In terms of pH, it is suggested to operate within 0.5 pH unit of the buffer's pK_a . In protein work, it is advised to operate at least 1 pH unit below the pI of the protein, to facilitate binding. Oxidizing agents, cationic detergents and cationic buffers should **not** be used with CM Sepharose® Fast Flow. Several publications¹⁻⁴ and theses^{5,6} cite use of CCF100 in their research.

Reagent

CM Sepharose® Fast Flow is offered as a suspension in 20% ethanol.

Approximate Exclusion Limit: average molecular mass of $\sim 4 \times 10^6$ Daltons

Ionic Capacity: 0.09–0.13 mmol/mL resin

Binding Capacity: 15 mg IgG (MW ~ 160 kDa) per mL gel; 50 mg RNase A (MW ~ 13.7 kDa) per mL gel

pH Stability: 4–13

Working temperature range: 4–40 °C

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

General Resin Preparation Procedure

1. Allow the ion exchange medium and ~ 10 column volumes (CV) of buffer to equilibrate to the temperature chosen for the chromatographic run.
2. Mix the pre-swollen suspension with starting buffer to form a moderately thick slurry, which consists of $\sim 75\%$ settled gel and 25% liquid.
3. Degas the gel under vacuum at the temperature of column operation.
4. Mount the column vertically on a suitable stand, out of the way of direct sunlight or drafts, which may cause temperature fluctuations.
5. Pour a small amount of buffer into the empty column. Allow the buffer to flow through spaces to eliminate air pockets.
6. Pour the suspension of ion exchange medium prepared in Step 3 into the column by allowing it to flow gently down the side of the tube, to avoid bubble formation.
7. For consistent flow rates and reproducible separations, connect a pump to the column.
8. Fill the remainder of the column to the top with buffer. Allow ~ 5 CV of buffer to drain through the bed at a flow rate at least 133% of the flow rate to be used in the procedure. The bed height should have settled to a constant height.
9. Using a syringe or similar instrument, apply the sample dissolved in starting buffer to the column. For isocratic separations, the sample volume should range between 1–5% of the column volume. If the chromatographic run involves elution with a gradient, the applied sample mass is of much greater importance than the sample volume, and the sample should be applied in a low ionic strength medium. Ion exchange is used both to concentrate and to fractionate the sample.
10. Elution:
 - If only unwanted substances in the sample are adsorbed, or if sample components are differentially retarded under isocratic conditions, the starting buffer can also be used as the eluent.

- Normally, however, separation and elution are achieved by selectively decreasing the affinity of the molecules for the charged groups on the resin by changing the pH and/or ionic strength of the eluent. This procedure is termed gradient elution.

11. Regeneration:

- Washing the column with a high ionic strength salt solution (such as 1 M NaCl), or changing the pH to the low and high extremes that are tolerable, is usually sufficient to remove all reversibly bound material.
- When necessary, lipids and precipitated proteins can be removed by washing with 1 CV of 1-2 M NaCl, followed by 1 CV of 0.1 M NaOH in 0.5 M NaCl.
- Rinse with several CV of water. Then re-equilibrate the resin with starting buffer.
- If base was used, adjust the pH of the resin to neutral before storing or using.

12. Storage: Sepharose® ion exchanger media may generally be stored at 2-8 °C in water or a dilute buffer of choice with 20% ethanol added as an antibacterial agent.

General Notes

Cation versus Anion Exchanger

- If sample components are most stable below their pI values, a cation exchanger should be used.
- If sample components are most stable above their pI values, an anion exchanger should be used.
- If stability is good over a wide pH range on both sides of the pI, either or both types of ion exchanger may be used.

Strong versus Weak Ion Exchanger

- Most proteins have pI values within the range 5.5-7.5, and can thus be separated on both strong and weak ion exchangers.
- When maximum resolution occurs at an extreme pH and the molecules of interest are stable at that pH, a strong ion exchanger should be used.

Choice of Buffer, pH, and Ionic Strength

- The highest ionic strength, which permits binding should normally be used.
- The required buffer concentration varies from substance to substance. Usually, an ionic strength of at least 10 mM is required to ensure adequate buffering capacity.
- As salts (such as buffers) help to stabilize proteins in solution, their concentration should be high enough to prevent denaturation and precipitation.

References

1. Wang, T. *et al.*, *Plant Physiol.*, **172**(4), 2107-2119 (2016).
2. Taglialaeta, A. *et al.*, *Mol. Cell*, **68**(2), 414-430.e8 (2017).
3. Nozawa, R.-S. *et al.*, *Cell*, **169**(7), 1214-1227.e18 (2017).
4. Krzyscik, M.A. *et al.*, *J. Vis. Exp.*, **(167)**, doi: 10.3791/61918 (2021).
5. Porkka, Kaija, "Solunulkoista ligniiniä tuottavaan kuusisolususpensioviljelmään (*Picea abies*) erittyvät kitinolyyttiset entsyymit ja niiden mahdolliset kasviperäiset substraattit" ("The chitinolytic enzymes secreted in the cell suspension culture of Norway spruce (*Picea abies*) producing apoplastic lignin, and their possible endogenous substrates"). University of Helsinki, M.Sc. thesis, p. 24 (2016).
6. Sims, Cynthe L., "Mechanism of Inhibition of Metallo-β-lactamase by Nucleic Acids". Texas Tech University, Ph.D. dissertation, pp. 24, 32 (2011).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Standard Warranty

The applicable warranty for the products listed in this publication may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

CCF100pis Rev 01/22 JJJ,MAM,GCY

