



3050 Spruce Street
Saint Louis, Missouri 63103 USA
Telephone 800-325-5832 • (314) 771-5765
Fax (314) 286-7828
email: techserv@sial.com
sigma-aldrich.com

Product Information

STREPTAVIDIN - 5 nm COLLOIDAL GOLD LABELED

Product Number **S 4188**

Storage Temperature 2 to 8 °C

Product Description

Streptavidin - 5 nm Colloidal Gold Labeled is prepared using purified Streptavidin from *Streptomyces avidinii* (Product No. S 4762) that is adsorbed to colloidal gold. It is suitable for detection of biotinylated compounds.

Colloidal gold is an electron-dense, non-fading marker useful as a probe in electron microscopy (TEM and SEM), light microscopy, and blotting.¹ It requires no additional processing for detection, but in some applications the signal can be enhanced by reaction with silver.^{2,3} It can be complexed with biomolecules by strong, non-covalent interactions.⁴

Generally, particles < 15 nm are most useful in transmission electron microscopy (TEM) and for applications where access to the probe may be hindered by larger particles.⁵ Particles >15 nm are more suitable for scanning electron microscopy (SEM), light microscopy, and blotting.

Reagent

Streptavidin - 5 nm Colloidal Gold Labeled is supplied as a clear, red, colloidal suspension in 0.01 M phosphate buffered saline, pH 7.4, containing 0.02% polyethylene glycol, 20 % glycerol and 15 mM sodium azide.

Particle size: 5 nm (Mean: 3 to 6 nm, the coefficient of variance of the particle size is <15% of the mean particle size).

Monodispersed with ≥ 50 % singlets

Gold conjugates are processed, after adsorption with proteins, to remove free protein and large aggregates. The concentration is expressed as absorbance at the absorption maximum (A_{520}).^{2,6}

Precautions and Disclaimer

Sodium Azide is considered highly toxic and highly reactive under certain conditions. Read the Material Safety Data Sheet carefully before use.

Preparation Instructions

Gold-conjugated proteins should be diluted for most applications. It is recommended that the diluent buffer contain 0.15 M saline, buffered at pH 6 to 8, and 0.05 % TWEEN[®] 20 to minimize background. Additional buffer supplements may be required for certain applications. Prior to application, allow the conjugate to equilibrate for at least 20 minutes in lower glycerol content. The optimum concentration of the conjugate should be determined empirically; a typical range is $A_{520}=1.0$ to 0.05 (1:2.5 to 1:50 dilution). Incubation times range from 0.5 to 12 hours.

Storage/Stability

Product may be stored for extended periods as packaged (undiluted) at 2 to 8 °C. Working dilution samples should be discarded if not used within 12 hours.

Procedure

Binding is evaluated by a "dot blot" assay modified from the method of Brada and Roth.² Serial dilutions are prepared from a 1 mg/ml positive control protein solution. One μ l of each solution is adsorbed onto a nitrocellulose membrane and allowed to dry. The gold conjugate is diluted to $A_{520} = 0.25$ with 0.01 M phosphate buffered saline, pH 7.4, containing 0.02% polyethylene glycol and 0.05 % TWEEN 20. The spotted membranes are incubated with the gold for 1 hour at 25 °C. The detection limit is the minimum amount of protein that can be detected as a pink-red spot on the membrane.

Product Profile

Streptavidin - 5 nm Colloidal Gold Labeled will detect 30 ng or less of biotinylated albumin (Product Number A 6043) by dot blot. References

References

1. Beesley, J.E., Proc. Royal Micr. Soc., **20**, 187 (1985).
2. Brada, D. and Roth, J., Anal. Biochem., **142**, 79 (1984).
3. Danscher, G., and Norgaard, J.D., Histochem. Cytochem., **31**, 1394 (1983).
4. Geoghegan, W.D., et al., Immunol. Comm., **7**, 1 (1978).
5. Slot, J.W., and Geuze, H.J., J. Cell. Biol., **90**, 533 (1981).
6. Tolson, N.D., et al., J. Microsc., **123**, 215 (1981).

TWEEN® is a registered trademark of the ICI Group.

LPG/MAM 3/02

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.