

Automation of Receptor-Ligand Binding Assays using the MultiScreen[®] HTS Filter Plate

Libby Kellard, Sonia Gil and Steven D. Sheridan, Ph.D.*
 Millipore Corporation, Life Sciences Division, Danvers, MA 01923

* Corresponding Author


Abstract

Perhaps the most critical screening parameter in the drug discovery process is the quantification of the specific affinity a drug has for a particular cellular receptor. Whether the process be carried out as a primary screening method for large compound libraries or as a secondary screening tool to rank compounds for binding affinity, high throughput, automation compatibility and accurate receptor-ligand binding analyses are required.

The design features of the new MultiScreen[®] HTS filter plate allows the plate to be easily gripped and moved around robotic decks and associated stackers. The plate can accommodate standard bar code labeling systems for ease of tracking when processing or storing numerous samples. The MultiScreen[®] HTS filter plate can be used for coincidence counting which provides the best signal to noise ratio.

The receptor ligand binding data presented here will demonstrate automation of the receptor ligand binding assay using the MultiScreen[®] HTS filter plate on the PerkinElmer Evolution[™] P³ Workstation. Scintillation counting and analysis were performed on a Wallac Microbeta[®] TriLux.

MultiScreen[®] HTS Filter Plate Assay System

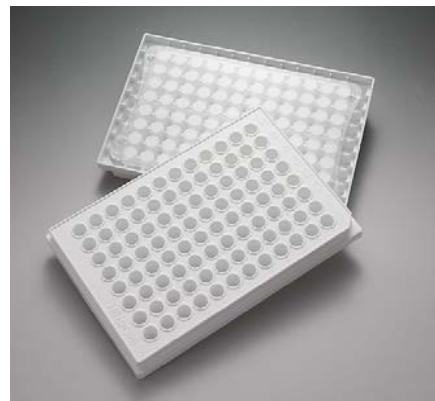

- 96 well format designed for highly sensitive and specific radiometric assay
- Removable underdrain for increased radiometric signal
- Completely isolated wells prevent radiometric signal crosstalk
- Completely ANSI-SBS compatible for full automation on robotic workstations
- Rigid side walls allow for easy barcode attachment
- Wide selection of filtration media including glass fiber


Methods and Materials

MultiScreen[®] (Millipore, FB glass fiber cat# MAFB N0B or HV membrane cat# MAHV N0B) and MultiScreen[®] HTS HV (Millipore, FB glass fiber cat# MSFB N0B or HV membrane cat# MSHV N0B) filter plates were pretreated with 0.1% polyethyleneimine (PEI) and washed with appropriate assay buffer.



Saturation Binding: 25 ng purified human estrogen β receptor (Sigma) or 8.75 μ g of a human Muscarinic M1 receptor expressing transgenic CHO cell membrane fragment bound receptor preparation (PerkinElmer) was incubated with serial dilutions of radiolabeled ligand [³H-estradiol (Sigma) and ³H-scopolamine (PerkinElmer NEN) respectively]. After 1 hour incubation in the filter plate, plates were washed 10 times by successive filling with binding buffer then vacuum. The plates were dried completely before the addition of Opti-Fluor Scintillation cocktail (PerkinElmer cat# 6013199). Non-specific binding was determined in a separate experiment with an excess of unlabeled competitor ligand. Specific binding was calculated as non-specific activity subtracted from total activity. Binding constants (K_d) were determined by fitting specific binding by free ligand concentration by non-linear regression and Scatchard analysis (shown) using Prism data software (www.Graphpad.com).

Displacement Binding: Receptor amounts used were the same as for above. Radioligand binding inhibition was determined with a constant radioligand concentration (0.2 nM ³H-estradiol or 0.6nM ³H-scopolamine) and serial dilutions of unlabelled competitor ligand (estradiol and pirenzepine, respectively) as compared to a control binding experiment without unlabelled ligand (% Control). Relative affinity values (IC₅₀) were determined by fitting binding inhibition values by non-linear regression using Prism data software.

In-Plate Receptor/Ligand Binding in MultiScreen[®] Filter Binding Plates

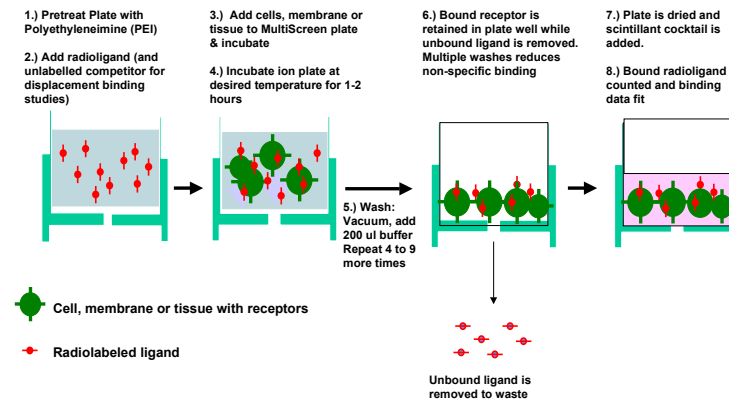

Automated Assays with MultiScreen[®] HTS on PerkinElmer Evolution[™] P³

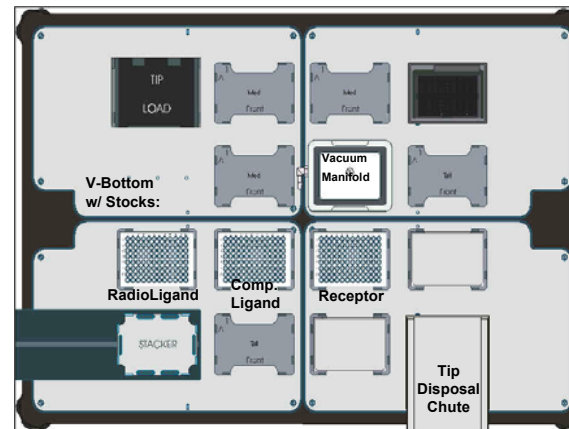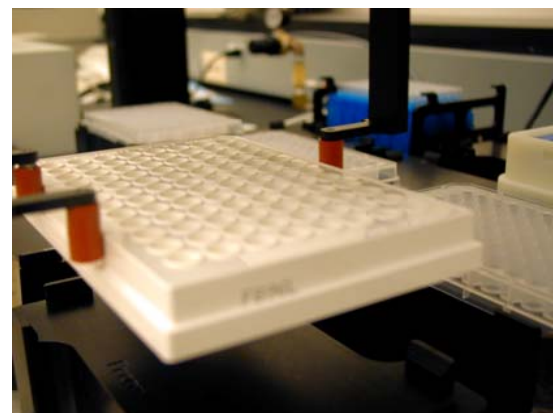


Radioligand, competitive non-radiolabeled ligand and receptor dilution, mixing, and filter plate washing are all performed on the PerkinElmer Evolution[™] P³. The program allows the user to input number of plates to run, number of filter plate washes, and whether the receptor-ligand binding reaction is performed directly in a filter plate or in a V-Bottom incubation plate followed by transfer to a filter plate

Solubilized Nuclear Hormone Receptor Binding on PEI Treated MultiScreen[®] with HV (0.45 μ M) Membrane Comparison of MultiScreen[®] and MultiScreen[®] HTS


Saturation Binding

Binding of Estradiol to Human Estrogen Receptor β

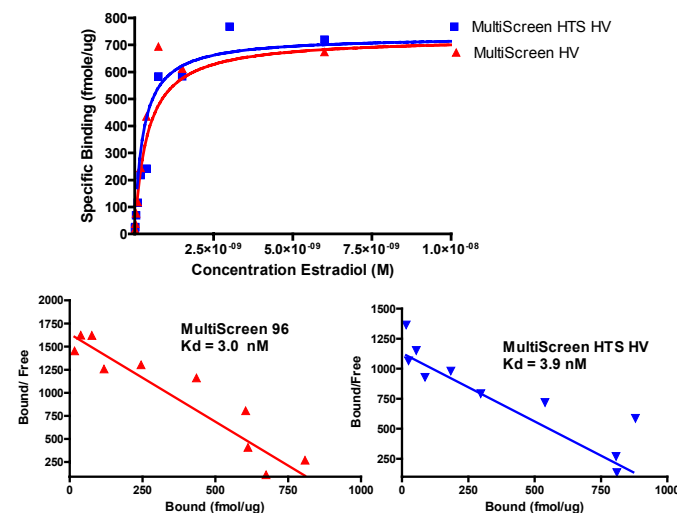

Displacement Binding

Unlabelled Estradiol Displacement Binding: Human Estrogen Receptor β

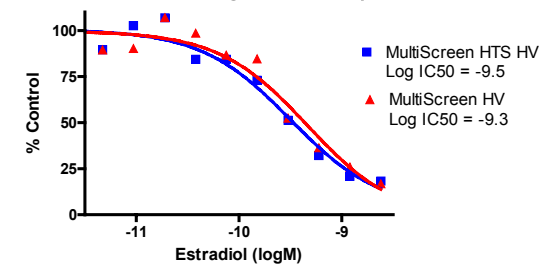


All Data Points in this Figure are the Average of Triplicate Experiments

G-Protein Coupled Receptor (GPCR) Binding on MultiScreen[®] with PEI Treated FB Glass Fiber Comparison of MultiScreen[®] and MultiScreen[®] HTS


Saturation Binding

Binding of Scopolamine to Human Muscarinic M1 Receptor

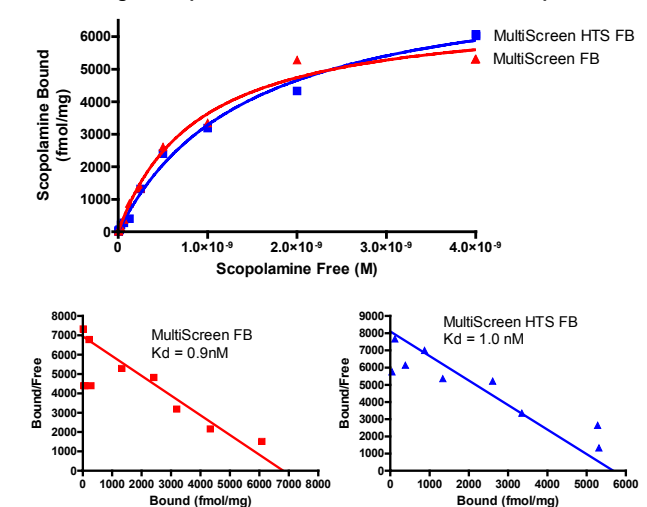

Displacement Binding

Unlabelled Pirenzepine Displacement Binding: Human Muscarinic M1 Receptor

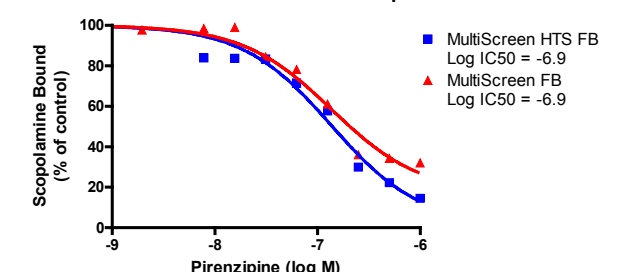

All Data Points in this Figure are the Average of Triplicate Experiments

Summary

- The Millipore MultiScreen[®] HTS 96 well filter plates with the PerkinElmer Evolution P³ Workstation allow for easy automation of receptor-ligand binding assays for both quantitative saturation binding and displacement screening experiments.
- Receptor ligand binding assays were performed with a soluble nuclear hormone receptor (the human estrogen β receptor) and a G-protein Coupled Receptor [(GPCR) the human Muscarinic M1 receptor]. Accurate and reproducible determinations of receptor specific activity (B_{max}), binding affinity (K_d) and IC₅₀ values of competitor ligands were measured.
- The MultiScreen[®] HTS filter plate is ANSI/SBS 2004 1-4 standard compliant thus allowing for the use of robotic gripper arms and can accommodate bar code labels.
- The new MultiScreen[®] HTS filter plate provides the same quantitative, robust and reproducible data acquisition as the original MultiScreen[®] filter plate platform.
- Receptor-ligand binding reactions can be performed entirely in the filter plate, saving the expense of an extra incubation plate and limiting the number of steps in the protocol for time savings.
- The availability of various filter media allows for the characterization of soluble receptors (e.g. Nuclear Hormone Receptors) and cell membrane embedded GPCRs.