

3050 Spruce Street, St. Louis, MO 63103 USA
Tel: (800) 521-8956 (314) 771-5765 Fax: (800) 325-5052 (314) 771-5757
email: techservice@sial.com sigma-aldrich.com

## **Product Information**

# Fibrinogen from human plasma

Catalog Number **F3879** Storage Temperature –20 °C

CAS RN 9001-32-5

#### **Product Description**

Molecular mass: ~340 kDa  $\lambda_{max}$  = 280 nm Extinction coefficient:  $E^{1\%}$  = 15.1 (solvent not given)

Fibrinogen, or Factor I, is a blood protein that is involved in clotting and is converted to fibrin by thrombin. Fibrinogen has an approximate molecular mass of 340 kDa. It consists of three non-identical pairs of disulfide-bonded chains. The  $\alpha$ -chain has an approximate molecular mass of 63.5 kDa, the  $\beta$ -chain 56 kDa, and the  $\gamma$ -chain 47 kDa. At the amino termini, the three chains are connected in a dimeric disulfide knot (DSK). A second DSK occurs later in the molecule. Fibrinogen has ~4% carbohydrate content. The Stokes radius for human fibrinogen is 10.7 nm.  $^6$ 

In general, fibrinogen from any mammalian source will cross-react with thrombin from any mammalian source. When any mammalian thrombin is injected into a different mammal, clotting will occur.

#### **Precautions and Disclaimer**

This product is for R&D use only, not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

### **Preparation Instructions**

The optimal way to solubilize fibrinogen is to layer it on top of warm (37 °C) saline; fibrinogen will not dissolve in water. The saline concentration can be in the range of 0.85–0.9%. The fibrinogen-saline solution can be gently agitated, but it must not be vortexed. The fibrinogen will slowly dissolve to give a hazy solution. Fibrinogen may be sterile-filtered, but may not go through a 0.1  $\mu m$  filter; a 0.2  $\mu m$  filter is suggested, with positive pressure using a syringe and "button" filter. Vacuum filtration should not be used, since this will lead to breakdown of the molecule during the filtration.

#### Storage/Stability

Store the product at -20 °C.

After sterile filtration, stock solutions of fibrinogen at 2.5 mg/ml may be stored refrigerated for ~1 week.

#### References

- The Plasma Proteins, I, 2nd ed., Putnam, F.W., ed., Academic Press (New York, NY: 1975), pp. 66 and
- Doolittle, R.F., The Structure and Evolution of Vertebrate Fibrinogen. Ann. N. Y. Acad. Sci., 408, 13-27 (1983).
- Henschen, A. et al., Covalent Structure of Fibrinogen. Ann. N. Y. Acad. Sci., 408, 28-43 (1983).
- 4. Blomback, M. et al., Fibrinogen Detroit--a Molecular Defect in the N-terminal Disulphide Knot of Human Fibrinogen? Nature, **218(137)**, 134-137 (1968).
- Takagi, T., and Doolittle, R.F., Amino Acid Sequence Studies on Plasmin-derived Fragments of Human Fibrinogen: Amino-terminal Sequences of Intermediate and Terminal Fragments. Biochemistry, 14(5), 940-946 (1975).
- Axelsson, I., Characterization of Proteins and other Macromolecules by Agarose Gel Chromatography. J. Chromatogr., 152, 21-32, (1978).

TD,GCY,RXR,MAM 12/15-1