

94029 Legionella Selective Supplement IV

An antibiotic supplement recommended for the selective isolation of Legionella species.

Composition
(per vial, sufficient for 500 ml medium)

Polymyxin B sulphate	25'000 units
Glycine	1.5 g
Anisomycin	40 mg
Vancomycin	0.5 mg
Bromo thymol blue	5 mg
Bromo cresol purple	5 mg

Directions

Rehydrate the contents of 1 vial aseptically with 5 ml of 50% ethanol. Mix well and aseptically add to 500 ml of sterile molten Buffered Charcoal Yeast Extract Agar, Base (Cat. No. 86558). Mix well and pour into sterile petri plates.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

