

## 67515 STEC Millichrome™ plus Supplement

Selective supplement for detection of Shiga-Toxin producing *E. coli* (STEC)

Package information:

5 vials (1 vial is sufficient for 1 L of final media)

Directions

- Aseptically rehydrate 1 vial with 10 mL of sterile water.
- Swirl well until complete dissolution.
- Add 1 vial rehydrated solution to 1L STEC Millichrome™ plus Agar base cooled at 45-50 °C.
- Swirl gently to homogenize.

### Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

