

Product No. I-7139
HUMAN INTERLEUKIN-2
From a Human T-Leukemic Cell Line

Lot 071H40231

Description

Interleukin-2 (IL-2, also known as T-cell growth factor) is an immunomodulatory factor produced by certain subsets of T-lymphocytes.¹ This lymphokine has proved useful in promoting long term growth of activated T-cells and related cell types. Interleukin-2 also has been shown to effect the activation and proliferation of NK cells, induce γ -interferon and B-cell growth factor secretion,^{2,3,4,5} and modulate the expression of the IL-2 receptor.⁶

Interleukin-2 has been isolated from a number of cell types^{7,8,9} and has also been produced by recombinant DNA technology.¹⁰ Human IL-2 (Natural, 15-17 kD glycoprotein) is secreted by a human leukemia T-cell line, after stimulation with T-cell mitogens.^{11,12}

Reagent

Interleukin-2 is supplied as a frozen solution containing 0.1% bovine serum albumin in 1ml of phosphate buffered saline containing 500 or 2,000 units of IL-2. The contents of the vial can be diluted further using a buffered solution or tissue culture containing 0.1-1% BSA or 1-10% serum. If aseptic technique is used, additional filtration should not be necessary and should be avoided due to possible adsorption to the filter membrane.

Storage

Store at -70°C. Prolonged storage of product or repeated freezing and thawing is **not** recommended and will result in decreased biological activity. After the initial thaw of the vial contents, store working aliquots at -70°C.

Performance Characteristics

Human Interleukin-2 has been tested in culture using a variation of the biological assay of Gillis, et al.¹³ IL-2 dependent cells are plated at low density in medium containing various dilutions of IL-2. After a specific incubation period, the cultures are pulsed with thymidine (³H), incubated further and then counted. One unit of biological activity is defined as the amount of the IL-2 required to induce half-maximum incorporation of thymidine.

IL-2 can be used at an approximate concentration of 20U/ml of medium for the culture of murine T-cells and approximately 100-200U/ml in LAK/NK cell activation or in proliferation studies.

This product has been tested in culture to ensure the absence of bacteria, yeasts, mold and mycoplasma.

References

1. Smith, K., Science, **240**, 1169 (1988).
2. Morgan, D., et al., Science, **193**, 1007 (1976).
3. Ortaldo, J., et al., J. Immunol., **133**, 779 (1984).
4. Farrar, J., et al., Immunol. Rev, **63**, 129 (1982).
5. Inaba, K., et al., J. Exp. Med., **158**, 2040 (1983).
6. ~~Smith, K., et al., Proc. Natl. Acad. Sci. USA, **82**, 1864 (1985).~~
7. Welte, K., et al., J. Exp. Med., **156**, 454 (1982).
8. Kiehl, E., et al., Eur. J. Biochem., **143**, 199 (1984).
9. ~~Smith, K., et al., Proc. Natl. Acad. Sci., USA, **80**, 5990 (1983).~~
10. Taniguchi, T., et al., Nature (London), **302**, 305 (1983).
11. Sieder, U., et al., Int. J. Cancer, **19**, 621 (1977).
12. Robb, R., Methods Enzymol., **116**, 493 (1985).
13. Gillis, S., et al., J. Immunol., **120**, 2027 (1978).

Potential Biohazard

Handle as if capable of transmitting infectious agents. Starting material tested and found negative for HBsAg and antibody to HIV.

Refer to material safety data sheet (MSDS).

Sigma warrants that its products conform to the information contained in this and other Sigma products. Purchaser must determine the suitability of the products for its particular use. See reverse side of invoice or packing slip for additional terms and conditions of sale.

Reissued 8/91