



Murine RNase Inhibitor

Product Number: RNAINH

Product Description

Murine RNase Inhibitor is mouse-source recombinant protein expressed in *E.coli*. It can inhibit the activity of RNase A, B or C in a non-competitive way, protecting RNA from degradation. Murine RNase Inhibitor is a thermo-stable RNase inhibitor. It still has enzyme activity when using thermo-stable reverse transcriptase. It is compatible with various commercialized reverse transcriptase and DNA polymerase. Compared with human-source RNase inhibitor, mouse-source RNase inhibitor doesn't contain 2 oxidation-sensitive Cys. And therefore, it has a higher antioxidant activity and is more suitable for high-DTT-sensitive experiments (i.e. qPCR).

Components

Components	RNAINH-MC2	RNAINH-MC10
Murine RNase Inhibitor (40 U/μl)	50 μl	50 μl

Storage

Store at -30 ~ -15°C and transport at ≤ 0°C

Application

It can be used to protect RNA from degradation.

- 1st strand cDNA synthesis in RT-PCR, PCR and qPCR.
- In vitro reverse transcription/translation.
- RNA isolation and purification.
- RNase protection assay.

This product won't interfere with other enzymes during application.

Source

It is purified from *E.coli*.

Unit Definition

One unit (U) is defined as the enzyme needed for inhibiting 50% activity of 5 ng RNase A.

The activity of RNase A is detected by quantifying the hydrolysis of Cyclic 2', 3'-CMP to 3'-CMP.

Notes

- The reaction temperature is 25 ~ 55°C. The RNase Inhibitor will be inactivated at temperature ≥ 65°C.
- Murine RNase Inhibitor can inhibit RNase activity under a broad spectrum of pH (pH 5.0 - 9.0). The highest inhibitory activity is obtained at pH 7.0 - 8.0.
- It can be inactivated by bubbling or stirring intensely (i.e. Vortexing).
- No inhibitory activity for RNase H and RNase T1.

Experiment Process

- Mix the following components in a RNase-free centrifuge tube and mix gently:

RNase-free ddH ₂ O	to 20 μl
5 × HiScript Buffer	1 μl
Oligo (dT) ₁₈ (50 μM)	1 μl
dNTP Mix (10 mM each)	1 μl
Murine RNase Inhibitor (40 U/μl)	1 μl
HiScript Reverse Transcriptase (200 U/μl)	2 μl
Template RNA	10 pg - 2.5 μg

- Incubate at 50°C for 45 min, then at 70°C for 15 min.
- The products can be stored at -20°C.

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

To place an order or receive technical assistance

In Europe, please call Customer Service:

France: 0825 045 645

Germany: 069 86798021

Italy: 848 845 645

Spain: 901 516 645 Option 1

Switzerland: 0848 645 645

United Kingdom: 0870 900 4645

For other countries across Europe, please call: +44 (0) 115 943 0840

Merck KGaA

Frankfurter Strasse 250

64293 Darmstadt, Germany

For other countries across Europe and the world, please visit:

[SigmaAldrich.com/offices](https://www.SigmaAldrich.com/offices)

For Technical Service, please visit:

[SigmaAldrich.com/techservice](https://www.SigmaAldrich.com/techservice)

[SigmaAldrich.com](https://www.SigmaAldrich.com)