

Technical Bulletin

ExtrAvidin®-Peroxidase

Buffered aqueous solution

E2886

Product Description

ExtrAvidin® is prepared from egg white avidin. It is a tetrameric protein containing four high affinity binding sites for biotin. ExtrAvidin® is a modified form of affinity purified avidin. ExtrAvidin® combines the high specific activity of avidin with the low background staining of streptavidin, a biotin binding protein produced by the bacteria *Streptomyces avidinii*. ExtrAvidin® binds biotin with the high affinity of egg white avidin, however, it does not exhibit the unwanted non-specific binding reported for egg white avidin at physiological pH, such as the staining of mast cells.

Uses

ExtrAvidin®-Peroxidase may be used in conjunction with biotin in techniques such as ELISA, immunohistochemistry, protein blotting or with biotinylated-DNA probes. It provides a convenient yet highly sensitive and very specific detection system.

Reagent

Supplied as a solution in 0.01 M phosphate buffered saline, pH 7.4, with 0.05% MIT as a preservative.

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Storage

Store at 2-8 °C. **Do Not Freeze.**

If slight turbidity occurs upon prolonged storage, clarify the solution by centrifugation before use.

Produce Profile

Immunohistology

A dilution of 1:100 of the conjugate was determined.

Dot blot

A minimum dilution of 1:4,000 of the conjugate was determined.

ELISA

A minimum dilution of 1:2,000 of the conjugate was determined.

Note: In order to obtain best results, it is recommended that each individual user determine their optimum working dilution by titration assay.

Specificity

The product is found to be reactive with Anti-Avidin, Cat. No. A5170, and non-reactive with Anti-Streptavidin, Cat. No. S6390, by immunoelectrophoresis.

Molar Ratio (ExtrAvidin®:Peroxidase)

0.5-0.9

Protein Concentration:

2.0-2.5 mg/mL by $E_{280}^{1\%} = 12.4$

Enzyme Activity

Enzyme activity is determined using 5% pyrogallol, Cat. No. P0381, in deionized water, pH 6.0, at 20 °C. One purpurogallin unit will form 1 mg of purpurogallin from pyrogallol in 20 seconds at pH 6.0, 20 °C.

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Standard Warranty

The applicable warranty for the products listed in this publication may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.