

Product Information

Pyruvate Kinase M2 human

Recombinant, expressed in *E. coli*, specific activity ≥ 100 units/ μ g protein

SAE0021

Product Description

CAS Registry Number: 9001-59-6

Enzyme Commission (EC) Number: 2.7.1.40

Synonyms: PKM2, Tumor M2-PK, OPA-interacting protein 3 (OIP-3), Pyruvate kinase 3 (PK3), Pyruvate kinase muscle isozyme, Cytosolic Thyroid hormone-binding protein (CTHBP)

Pyruvate kinase (PK) is an important glycolytic enzyme that catalyzes the last step of glycolysis, transferring a phosphate group from phospho(enol)pyruvate to ADP to yield one molecule of ATP and one molecule of pyruvate. Human Pyruvate kinase appears in four major forms:

- L (found in liver)
- R (found in erythrocytes)
- M1 (found in muscle and brain tissues)
- M2 (the dominant isoform, found in proliferating tissues)

The M2 isoform (PKM2) is involved in the regulation of cell proliferation, cell size, and apoptosis.¹ PKM2 is predominantly found in many human cancers and in embryonic tissues.²

While all PK enzymes require potassium ions for their activity, PKM2 also requires fructose-1,6-bisphosphate (FBP) to stabilize its active tetrameric form.³ PKM2 functions as a cytosolic thyroid hormone binding protein (CTHBP)⁴ and as a phosphotyrosine binding protein.⁵ The binding of phosphotyrosine peptides to PKM2 results in the release of the allosteric activator FBP leading to inhibition of PKM2 activity. This regulation of PKM2 diverts glucose metabolites from energy production to anabolic processes when cells are stimulated by certain growth factors such as IGF.⁵

This product has been used as a standard for an activity assay for serum tumor PKM2.⁶ Other research publications⁷⁻⁹ and dissertations¹⁰ have cited use of product SAE0021 in their research protocols.

Precautions and Disclaimer

This product is for R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Product

This recombinant human PKM2 (P14618) product is expressed in *E. coli*, with a predicted molecular mass of 58 kDa. The product is lyophilized from a buffered solution containing phosphate buffer, pH 7.5, NaCl, DTT, and a carbohydrate stabilizer.

Unit definition: One unit will convert 1.0 μ mole of phospho(enol)pyruvate to pyruvate per minute at pH 7.6 at 37 °C, in the presence of 1 mM fructose-1,6-bisphosphate.

Storage/Stability

Store the product at -20 °C.

Preparation Instructions

Reconstitute in 30% glycerol or in 30 mM fructose-1,6-bisphosphate to a final concentration of 1-2 mg protein/mL.

After reconstitution, the product can be stored at 2-8 °C for at least 24 hours with minimal loss of activity or frozen in working aliquots at -20 °C. Frozen aliquots can be kept for at least one month. Repeated freezing and thawing is not recommended. Do not store in a frost-free freezer.

References

1. Spoden, G. A. *et al.*, *Exp. Cell Res.*, **315(16)**, 2765-2774 (2009).
2. Iqbal, M. A. *et al.*, *FEBS Lett.*, **588(16)**, 2685-2692 (2014).
3. Ashizawa, K. *et al.*, *Biochemistry*, **30(29)**, 7105-7111 (1991).
4. Kato, H. *et al.*, *Proc. Nat. Acad. Sci. USA*, **86(20)**, 7861-7865 (1989). (Correction: *Proc. Nat. Acad. Sci. USA*, **87(4)**, 1625 (1990).)
5. Christofk, H. R. *et al.*, *Nature*, **452(7184)**, 181-186 (2008).
6. Guan, M. *et al.*, *Ann. Clin. Lab. Sci.*, **47(6)**, 676-686 (2017).
7. Wubben, T. J. *et al.*, *Sci. Rep.*, **10**, 2990 (2020).
8. Patel, S. *et al.*, *Eur. J. Pharm. Sci.*, **170**, 106112 (2022).
9. Ko, Y. *et al.*, *Proc. Nat. Acad. Sci. USA*, **120(20)**, e2300763120 (2023).
10. Biyik-Sit, Rumeysa, "Investigating a novel function for phosphoserine aminotransferase 1 (PSAT1) in epidermal growth factor receptor (EGFR)-mediated lung tumorigenesis". University of Louisville, Ph.D. dissertation, p. 33 (2021).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Terms and Conditions of Use

Warranty, use restrictions, and other conditions of sale may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2018-2023 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.
SAE0021dat Rev 07/23 DT,GCY,MAM

