

Instruction Manual For

phospho-Src (Tyr418) STAR ELISA Kit Catalog # 17-468

Sufficient reagents for 96 assays per kit

Contents		Page
Ι.	TEST PRINCIPLE	2
II.	BACKGROUND	3
III. Re	ASSAY SENSITIVITY, DETECTION LIMITS and SPECIES	4
IV.	STORAGE OF KIT COMPONENTS	4
V.	KIT COMPONENTS, PRECAUTIONS AND TECHNICAL NOTES	4
VI.	PREPARATION OF SAMPLE	5
VII.	REAGENT PREPARATION	6
VIII.	ASSAY PROTOCOL	8
IX.	CALCULATION OF RESULTS	9

upstate

The expertise of Upstate® is now a part of Millipore


sales orders: 800 233 3991 outside the U.S. 434 975 4300 tech support: 800 548 7853 fax orders: 866 831 3991 www.millipore.com FOR RESEARCH USE ONLY.

NOT RECOMMENDED OR INTENDED FOR DIAGNOSIS OF DISEASE IN HUMANS. DO NOT USE IN HUMANS.


I. TEST PRINCIPLE

The UPSTATE[®] colorimetric STAR (Signal Transduction Assay Reaction) ELISA kit is a solid phase sandwich enzyme linked immunosorbent assay that provides a fast, sensitive method to detect specific levels of signaling targets in whole cell extracts. The Src plate is coated with a specific mouse monoclonal Src capture antibody on the microwells of the 96-well clear plate. Sample lysate or the standard included in the kit are incubated in the microwells allowing Src antigen to be captured in the plate wells. The plate is then washed to remove any non-bound unspecific material. The wells are then incubated with a specific rabbit anti-Src antibody to detect the captured phospho-Src (Tyr418) on the plate well. The unbound detection antibody is washed away followed by incubation with an HRP-conjugated anti-rabbit antibody. This allows for a sensitive enzymatic detection of the sample. After the addition of TMB substrate and stop solution the absorbance is measured at 450 nm using a plate reader.


The entire assay takes less than 5 hours to complete with minimal hands-on time. Many of the reagents are supplied in ready-to use formulations for ease of use. The kit also includes a standard that is run as both a positive control and to develop a standard line of detection.

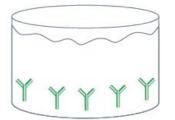


Step 1. 96 well clear plates coated with a specific mouse monoclonal capture antibody.

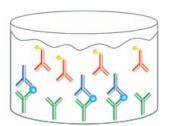


Step 4. Wash unbound detection antibody and add HRP-conjugated anti-rabbit antibody.

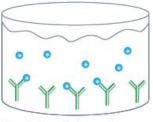


Step 2. Sample lysate (or standard) is incubated in the microwells binding to the capture antibody.

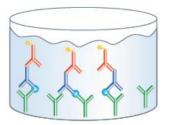

Step 5. Wash and incubate with TMB enzymatically activated detection reagent.

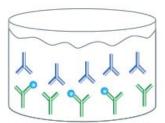

Step 3. Wash away unbound, non-specific material. Add specific rabbit total or phosphospecific antibody in each sample well.

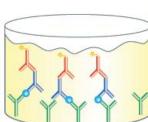


Vell. Capture Antibody Analyte Capture Antibody Detection Antibody HRP-conjugated secondary antibody

Step 6. Add stop solution and measure absorbance.

II. Src BACKGROUND

Discovery of the Rous Sarcoma Virus (RSV) and the virus encoded v-Src kinase lead to discovery of the cellular homologue c-Src, a non-receptor tyrosine kinase associated with the development of numerous human cancers. pp60c-Src is the prototype non-receptor tyrosine kinase. Homology to Src in three regions is shared by a number of proteins, SH1 being the catalytic domain, SH2 being a phospho-tyrosine binding domain, and SH3 being a proline-binding domain. Highly activated forms of c-Src kinase are associated with colon cancers and metastatic progression. Tyrosine residue 416 (equivalent to tyrosine residue 418 in human c-Src) is phosphorylated by autocatalysis and is highly correlated with the kinases activation state and transforming ability. The assay kit is designed to measure the activation state, and thereby transforming potential, of the Src kinase by assessing the phosphorylation levels at tyrosine 418.

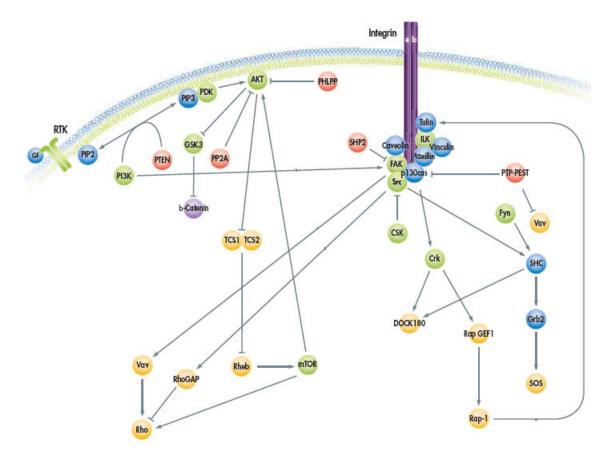

III. ASSAY SENSITIVITY, DETECTION LIMITS and SPECIES REACTIVITY

Sensitivity: 1 unit/mL.

Range of Detection: 1.6 to 100 units/mL

Species Reactivity: Human, mouse and rat

NOTE: This data is presented for reference use only and should not be used to interpret actual assay results. A standard curve must be generated for each assay.

IV. STORAGE OF KIT COMPONENTS

Maintain the unopened kit at 2-8°C until expiration date.

V. KIT COMPONENTS

- 1. <u>Capture Plate pre-coated with anti-Src antibody</u>: (Part No. 17-468A) One pre-coated 96-stripwell immunoplate sealed in a foil pouch.
- <u>Anti-phospho-Src (Tyr418) detection antibody</u>: (Part No. 17-468B) One bottle (11 mL) of anti-_phospho-Src (Tyr418) detection antibody containing sodium azide, ready to use.
- 3. <u>ELISA Diluent</u>: (Part No. 17-468C) One bottle (25 mL) of ELISA Diluent containing sodium azide, ready to use.
- 4. <u>25X ELISA Wash Buffer</u>: (Part No. 17-468D) One bottle (50 mL) of 25X ELISA Wash Buffer.
- 5. <u>Anti-Rabbit IgG HRP conjugate</u>: (Part No. 17-468E) One vial (125 μL) of 100X anti-rabbit HRP conjugate containing thimerosol.
- 6. <u>HRP Diluent</u>: (Part No. 17-468F) One bottle (25 mL) of HRP Diluent containing thimerosol.
- 7. <u>TMB Solution</u>: (Part No. 17-468G) One bottle (25 mL) of stabilized tetramethylbenzidine (TMB), ready to use.
- 8. <u>Stop Solution</u>: (Part No. 17-468H) One bottle (25 mL) of stop solution, ready to use.
- 9. <u>Phospho-Src (Tyr418) Standard</u>: (Part No. 17-468I) Two vials of phospho-Src (Tyr418) standard, lyophilized.
- 10. <u>Plate Covers</u>: Two plate covers.

Materials Not Supplied

- 1. Multi-channel or repeating pipettes
- 2. Plate shaker (optional)
- 3. Pipettors & tips capable of accurately measuring 1-1000 L
- 4. Graduated serological pipettes
- 5. 96-well microtiter Plate Reader with 450 nm filter
- 6. Graphing software for plotting data or graph paper for manual plotting of data
- 7. Microfuge tubes for standard and sample dilutions
- 8. Mechanical vortex
- 9. 1 liter container
- 10. Distilled or deionized water

Precautions

- The instructions provided have been designed to optimize the kit's performance. Deviation from the instructions may result in suboptimal performance of the kit and the failure to produce accurate data.
- Caustic Material: Stop Solution. Caution: Eye, hand, face, and clothing protection should be worn when handling this material.
- Safety Warnings and Precautions: This kit is designed for research use only and not recommended for internal use in humans or animals. All chemicals should be considered potentially hazardous and principles of good laboratory practice should be followed.
- The Detection Antibody and Elisa Diluent contain sodium azide. Sodium azide may react with copper and lead plumbing to form highly explosive metal azides. Upon disposal, flush with large amounts of water to prevent azide build-up. Avoid contact with skin.
- The Anti-Rabbit IgG HRP Conjugate and HRP Diluent contain thimerosal. Thimerosal is highly toxic by inhalation, contact with skin or if swallowed. Thimerosal is a possible mutagen and should be handled accordingly.

Technical Notes

- All kit reagents should be at room temperature (20°C to 25°C) prior to use.
- Do not use reagents beyond the expiration date of the kit.
- Do not mix or interchange reagent from various kit lots.
- Manual Plate Washing: Vigorous washing and complete removal of all liquid by

aspiration at the end of each washing step is very important to obtain low background values. Gentle agitation during the wash steps or a 2-3 minute soak may reduce background values.

• The desiccant enclosed in the 96-well capture plate pouch will keep the plate stable when stored at 2° to 8°C should the plate loose its seal during shipping.

VI. PREPARATION OF SAMPLE

- 1. Culture cells stimulating Src activation as desired.
- Remove culture media and wash cells twice ice-cold with 1X TBS (Tris Buffered Saline) or PBS (Phosphate Buffered Saline). Discard supernatant
- 3. Add 5-10 mL of cold 1X RIPA containing protease inhibitors per 150 mm tissue culture plate.

Note: 10 mL of 1X RIPA containing protease inhibitors can be prepared by adding 10 μ L of 1 μ g/ μ L Leupeptin, 10 μ L of 1 μ g/ μ L Aprotinin, 10 μ L of 1 μ g/ μ L Pepstatin, 100 μ L of 100mM PMSF, 1 mL 10X RIPA (Cat. No. 20-188) to 8.87 mL of distilled or deionized water.

- 4. Scrape cells from plate with a rubber policeman.
- Transfer cells in RIPA buffer to a microcentrifuge tube and incubate on ice for 15 minutes. Vortex tube for 10 seconds or sonicate briefly for 10 seconds.
- 6. Clarify lysate by centrifugation at 12,000 rpm for 10 minutes at 4°C in a microcentrifuge prior to use.
- The assay tolerates a maximum of 20 μL per well of lysate prepared in 1X RIPA (20-188) diluted in ELISA Diluent.
- 8. Cell extract containing SDS must be diluted to 0.01% SDS using ELISA Diluent prior to use.
- 9. Collect the supernatant and calculate protein concentration using a Bradford Assay or by densitometry.
- 10. Keep samples cold and use immediately if not freezing away for later use. Samples can be frozen and stored at -80° C. It is suggested to use the cell lysate be immediately used or use within 6 months if storing at -80° C. Avoid repeated freeze thaws.

Further information of lysate preparation protocols can be obtained at <u>http://www.millipore.com</u>

VII. REAGENT PREPARATION

1. 1X Wash Buffer

Warm the 25X ELISA Wash Buffer to room temperature and mix to ensure that any precipitated salts have re-dissolved. For 500 mL of Wash Buffer, combine 20 mL of 25X ELISA Wash Buffer and 480 mL distilled or deionized water. Stir to homogeneity. Wash Buffer can be stored for up to 4 weeks at 2-8°C. Discard the Wash Buffer if it becomes turbid or if a precipitate develops.

2. Anti-Rabbit IgG HRP Conjugate

Dilute the anti-Rabbit IgG HRP Conjugate 100-fold with HRP Diluent immediately before use. Prepare 1 mL for each strip used.

3. Standard

<u>Note</u>: When opening lyophilized Standard, remove rubber stopper gently as the lyophilizate may have become dislodged during shipping.

Reconstitute the standard with the volume of ELISA Diluent specified on the vial label to give a concentration of 100 units/mL. Gently swirl the vial and allow the vial to sit for 10 minutes to ensure the material is completely reconstituted. The standard should be reconstituted immediately before the assay. This stock material (tube #1) is then used to generate a standard curve. A suggested 2-fold dilution scheme is as follows:

- a) Label 7 test tubes #2-7 and "0 dose". Add 250 L of the ELISA Diluent to tubes #2-7 and "0 dose".
- b) Add 250 L of the stock Standard solution [100 units/mL] to tube #2 and vortex. This is Standard tube #2 with a concentration of 50 units/mL.
- c) Standards #2-7 are then prepared by performing a 2-fold serial dilution of the preceding standard. Refer to Fig. 1. For example, to make Standard #3, remove 250 L of Standard #2 and add it to tube #3 and vortex and so on. Do not add any Standard to the "0 Dose" Standard tube.

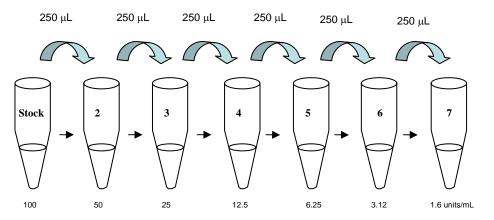


Figure 1: Recommended 2-fold Serial Dilution of Standard

Note: The Standard curve can set up with a different serial dilution scheme by making appropriate adjustments to the dilution pattern.

VIII. ASSAY PROTOCOL

- 1. Prepare the reagents as described in the Reagent Preparation section.
- 2. Place the desired number of strips in the strip well plate holder. (Re-bag the extra strips and return unused strips to refrigerator for future use.)
- 3. Add 100 μ L of either the Standards 1 through 7 or the samples to wells. Add 100 μ L of the zero dose to the control wells. It is recommended that standards and samples be run in duplicate. Wells reserved for TMB blanks should be left empty.

Note: A standard curve must be run at each setting.

4. Seal the plate with a plate sealer. Incubate the plate for 2 hours at room temperature or at 2-8°C overnight (on a shaker if possible).

5. **IMPORTANT WASH STEP:**

Gently remove the plate sealer and wash the plate at least 4 times. A thorough washing of the plate is extremely important to reduce background. We recommend using a multi-channel pipette to fill each well with 250 μ L of diluted Wash Buffer. Fluid removal from the wells is best accomplished by inverting the plate over a sink and flicking the fluid out of the wells and then blotting the plate on clean paper towels. Using the multichannel pipet add 250 μ L of Wash Buffer to each well; flick and blot the plate. Repeat this procedure for a total of 4 times.

For users of automatic plate washers: It is important to ensure that the wash apparatus is properly maintained and operating correctly. Tubing and tips can easily become clogged, leading to incomplete washing and inadequate aspiration of wells. The result may be poor precision and an unsuitable standard curve. For best results, we recommend at least 4 wash cycles.

- 6. Add 100 μ L of the detection antibody to each well. Cover the plate and incubate at room temperature for 1 hour (on shaker if possible). Wash as described in Step 5.
- 7. Add 100 μ L of a 1:100 dilution of the anti-Rabbit IgG HRP Conjugate to each well. Cover the plate and incubate at room temperature for 45 minutes (on shaker with mild agitation if possible). Wash as described in Step 5. Remove all fluid from the wells and blot the wells dry.
- 8. Add 100 μ L of the TMB Solution to each well. Incubate at room temperature in the dark for 10 to 45 minutes, monitor the color development. Stop the reaction by adding 100 μ L of Stop Solution to each well. Immediately read the plate at 450 nm. Plate should be read within 1 hour of adding the stop solution.

9. The plate reader may be blanked against a TMB blank prepared by adding 100 μ L of stop solution to 100 μ L of the TMB solution.

CAUTION: Bubbles in the wells will cause inaccurate readings. Ensure that all bubbles are removed prior to taking the absorbance reading.

NOTE: For very low starting protein levels, samples can be placed at 37°C during the final incubation to obtain greater sensitivity.

VII. CALCULATION OF RESULTS

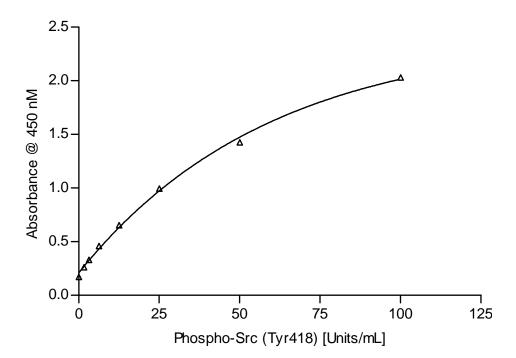


Figure 2. Typical phospho-Src (Tyr418) Standard Curve 100 μ L of progressive 2 fold dilutions of the phospho-Src (Tyr418) standard included in the kit and run as described in the assay instructions.

NOTE: This data is presented for reference use only and should not be used to interpret actual assay results. A standard curve must be generated for each assay.

Warranty

Millipore Corporation ("Millipore") warrants its products will meet their applicable published specifications when used in accordance with their applicable instructions for a period of one year from shipment of the products. MILLIPORE MAKES NO OTHER WARRANTY, EXPRESSED OR IMPLIED. THERE IS NO WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE. The warranty provided herein and the data, specifications and descriptions of Millipore products appearing in Millipore's published catalogues and product literature may not be altered except by express written agreement signed by an officer of Millipore. Representations, oral or written, which are inconsistent with this warranty or such publications are not authorized and if given, should not be relied upon.

In the event of a breach of the foregoing warranty, Millipore's sole obligation shall be to repair or replace, at its option, the applicable product or part thereof, provided the customer notifies Millipore promptly of any such breach. If after exercising reasonable efforts, Millipore is unable to repair or replace the product or part, then Millipore shall refund to the Company all monies paid for such applicable Product. **MILLIPORE SHALL NOT BE LIABLE FOR CONSEQUENTIAL, INCIDENTAL, SPECIAL OR ANY OTHER DAMAGES RESULTING FROM ECONOMIC LOSS OR PROPERTY DAMAGE SUSTAINED BY ANY COMPANY CUSTOMER FROM THE USE OF ITS PRODUCTS.**

©2008 - 2009: Millipore Corporation. - By Millipore Corporation. All rights reserved. No part of these works may be reproduced in any form without permissions in writing.

Unless otherwise stated in our catalog or other company documentation accompanying the product(s), our products are intended for research use only and are not to be used for any other purpose, which includes but is not limited to, unauthorized commercial uses, in vitro diagnostic uses, ex vivo or in vivo therapeutic uses or any type of consumption or application to humans or animals.