

3050 Spruce Street
Saint Louis, Missouri 63103 USA
Telephone 800-325-5832 • (314) 771-5765
Fax (314) 286-7828
email: techserv@sial.com
sigma-aldrich.com

# **ProductInformation**

# Phleomycin from Streptomyces verticillus

Product Number **P 9564** Storage Temperature 2-8 °C

## **Product Description**

Molecular Formula:  $C_{55}H_{85}O_{21}N_{20}S_2Cu \bullet HCI$ 

Molecular Weight: 1526.5 CAS Number: 11006-33-0

Phleomycin is a glycopeptide antibiotic, structurally related to the antibiotic, bleomycin. This purified phleomycin is mainly composed of phelomycin D1 (terminal amine is agmatine) in a copper chelate and hydrochloride salt form.

Phleomycin blocks S-phase entry in the cell cycle. While phleomycin can damage DNA similar to bleomycin, it is not used as an anticancer agent. Although the mechanism of action of phleomycin has not been clearly defined, it is reportedly unable to intercalate DNA due to a modified bithiazole moiety. The RAD6 DNA repair gene is essential for phleomycin resistance in mutant yeast.

Phleomycin is used as a selective agent in molecular genetics experiments. Phleomycin has a broad spectrum of toxicity, exhibiting *in vivo* activity against bacteria, eukaryotic organisms, and plant and animal cells. Phleomycin, can thus be useful for identification and selection of a variety of cell types carrying a phleomycin resistant gene, like the *Sh ble* gene.

Phleomycin has toxicity against bacteria ( $E.\ coli$ ), and is used at 5 µg/ml in media for selecting resistant transformants. It is also used at 5-50 µg/ml for eukaryotic microorganisms, at 10 µg/ml for yeasts, at 10-50 µg/ml for fungi, and at 5-25 µg/ml for plants. Since the sensitivity of cultured cells to phleomycin is increased at higher pH, a lower concentration of phleomycin can be used in a higher pH medium, for complete cell growth inhibition. The activity of phleomycin is reduced by a factor of 2-3 in hypertonic media.

#### **Precautions and Disclaimer**

For Laboratory Use Only. Not for drug, household or other uses.

# **Preparation Instructions**

Phleomycin is freely soluble in water (20 mg/ml), yielding a light, blue solution. It is also slightly soluble in methanol.

### Storage/Stability

The powder is very hygroscopic and should be stored desiccated. Phleomycin is sensitive to high concentrations of acids, but solutions can be exposed for a short-term to dilute acids without adverse affects.

Solutions can be sterile filtered and stored as singleuse aliquots at -20 °C.

#### References

- He, C.H., et al., A Saccharomyces cerevisiae phleomycin-sensitive mutant, ph140, is defective in the RAD6 DNA repair gene. Can. J. Microbiol., 42, 1263-1266 (1996).
- Povirk, L.F., et al., Copper(II) bleomycin, iron(III) bleomycin, and copper(II) phleomycin: comparative study of deoxyribonucleic acid binding. Biochemistry, 20, 665-671 (1981).
- Robson, C.N., et al., Bleomycin and X-rayhypersensitive Chinese hamster ovary cell mutants: genetic analysis and cross-resistance to neocarzinostatin. Mutat. Res., 193, 157-165 (1988).
- Sleigh, M.J., and Grigg, G.W., Sulphydrylmediated DNA breakage by phleomycin in *Escherichia coli*. Mutat. Res., 42, 181-190 (1977).

ARO/JRC 1/03