

3050 Spruce Street, St. Louis, MO 63103 USA
Tel: (800) 521-8956 (314) 771-5765 Fax: (800) 325-5052 (314) 771-5757
email: techservice@sial.com sigma-aldrich.com

Product Information

IgG from porcine serum

reagent grade

Catalog Number I4381

Product Description

Purified pig IgG is isolated from pooled normal pig serum by fractionation and ion-exchange chromatography.

Reagent

Supplied as an essentially salt-free lyophilized powder.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

Storage

Prior to reconstitution, store at 2-8 °C. After reconstitution, store aliquots at -20 °C. Repeated freezing and thawing is not recommended.

Reconstitution

The product is to be reconstituted to a minimum of 1 mg/ml with a 150 mM NaCl solution. Protein concentration should be determined after reconstitution by absorbance at 280nm (E280 1% = 14).

Uses

Purified pig IgG may be used as a reference antigen, standard, blocking agent, or coating protein in a variety of immunoassays including ELISA, dot immunobinding, Western immunoblotting, immunodiffusion, and immunoelectrophoresis. Other applications include starting materials for the preparation of immunogens and solid-phase immunoadsorbents.

Performance Characteristics

Purity: ≥ 95% (SDS-PAGE)

Identity (Immunoelectrophoresis): Single arc of precipitation in the gamma region versus anti-porcine whole serum and anti-porcine IgG. Minor gamma arcs may be present versus anti-porcine whole serum.

Sodium (ICP): ≤1.0%

IDC,PHC 06/12-1