

HYCON® Agar Strips YM

Technical Data Sheet

Ordering number: 1442420050

HYCON® Agar Strips YM are ready-to-use culture media for assessment of airborne yeasts and molds with HYCON® Microbial Air Samplers, i.e. RCS® High Flow Touch, RCS® High Flow, RCS® Plus, RCS® Plus Ex, and RCS® Standard.

Each agar strip is individually sealed in the transparent, primary package. The agar is filled in a flexible backing film. The formulation of the medium is based on a Rose Bengal Agar for selective detection of yeasts and molds.

Mode of Action

Rose-Bengal Chloramphenicol Agar is a selective medium for the enumeration of yeasts and molds. The medium has a neutral pH. Chloramphenicol, streptomycin and rose bengal are included as selective agents to inhibit bacterial growth. Rose bengal also restricts the size and height of colonies of the more rapidly growing molds, which facilitates the detection of slow-growing fungi by preventing overgrowth by more rapidly growing species.

Typical Composition

Soytone	6 g/l
D(+) Dextrose	15 g/l
Malt extract	5 g/l
Magnesium sulfate	0.5 g/l
Yeast extract	1 g/l
Agar	15 g/l
Selective Mix (Rose Bengal, Streptomycin, Chloramphenicol)	
Supplements and buffers	

The appearance of the medium is clear and pink. The pH value is in the range of 7.1 to 7.5. The medium can be adjusted and/or supplemented according to the performance criteria required.

Application and Interpretation

Prior to use the agar strip should be equilibrated to room temperature. Please check each agar strip before use to verify sterility and take care on aseptic handling in order to avoid false positive results. Contaminated or dehydrated agar strips should not be used for sampling.

Open the wrapper approximately at 1/3 by peeling back the plastic seal at the rounded side of the wrapper. Remove the agar strip with the coated side facing downwards. Insert the agar strip into the opening of the rotor, or the impeller drum according to the directions outlined in the user manual of the respective microbial air sampler. Place the instrument into required position, choose the appropriate sample volume and start the air sampling procedure.

When sampling is finished, remove the agar strip and place it back into the original wrapper. Seal the wrapper with an adhesive tape or Cover Slides (Order. No. 1.44111.0100). Label the wrapper e.g. with a waterproof pen for identification. The closed agar strips are transferred to an incubator.

Several recommendations are given by different guidelines for incubation: according to USP <1116> the agar media used for environmental monitoring should be incubated between 20 and 35 °C for not less than 72 hours. According to the FDA Aseptic Guide the agar strips for determination of the total yeast and mold count should be incubated at 20 to 25 °C for 5 to 7 days. Individual incubation conditions can be chosen and should be validated at the application site.

Finally, the number of CFU per slide is examined.

Grown colonies may be identified using suitable methods related to root cause analysis programs or to support sanitizing management.

Important Notes

- Practice aseptic technique when handling agar strips.
- The coated surface of the agar strips should face down during incubation in order to avoid the formation of satellites by condensing water.

Storage

The product can be used until the expiry date if stored in the original box, protected from light and properly sealed at the temperature range indicated on the box label. The total shelf from the date of production is 6 months.

Condensation can be prevented by avoiding quick temperature shifts and mechanical stress. Upon storage agar strips should not be placed near heat sources such as refrigerators with heat-emitting condensers. Boxes should be stored with the coated side of the agar strip facing downwards.

Disposal

Please mind the respective regulations for the disposal of used culture medium (e.g. autoclave for 20 min at 121°C, disinfect, incinerate etc.).

Quality Control

Control Strains	ATCC #	Inoculum	Incubation	Recovery
<i>Candida albicans</i>	10231	10-100 CFU	≤ 72h at 20-25°C	50-200%
<i>Aspergillus brasiliensis</i>	16404	10-100 CFU	≤ 72h at 20-25°C	50-200%
<i>Staphylococcus aureus</i>	6538	> 100 CFU	24-48h at 30-35°C	≤ 30%
<i>Bacillus subtilis</i>	6633	> 100 CFU	24-48h at 30-35°C	≤ 30%

Please refer to the actual batch related Certificate of Analysis.

Quality

This product is manufactured in a Millipore SAS facility whose quality management system is approved by an accredited registration body to ISO 9001 quality standard.

This product is manufactured in a Millipore SAS facility whose environmental management is approved by an accredited registration body to the appropriate ISO 14001 systems standard.

Literature

EUROPEAN COMMISSION (2008) Volume 4 EU Guidelines to Good Manufacturing Practice Medicinal Products for Human and Veterinary Use Annex 1 Manufacture of Sterile Medicinal Products (corrected version)

FDA (2004) Guidance for industry: Sterile Drug Products Produced by Aseptic Processing — Current Good Manufacturing Practice

ISO 14698-1:2003: Cleanrooms and associated controlled environments - Biocontamination control - Part 1: General principles and methods

United States Pharmacopoeia 39 NF 34 (2016): <1116> Microbiological Control and Monitoring of Aseptic Processing Environments

Merck KGaA
64271 Darmstadt, Germany Fax:
+49 (0) 61 51 / 72-60 80
mibio@merckgroup.com

Find contact information for your country at:
www.merckmillipore.com/offices

For Technical Service, please visit:
www.merckmillipore.com/techservice

www.merckmillipore.com/biomonitoring

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

HYRON, Merck, Millipore, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. Detailed information on trademarks is available via publicly accessible resources.

© 2019 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

The life science business of Merck operates as
MilliporeSigma in the U.S. and Canada.