



3050 Spruce Street
Saint Louis, Missouri 63103 USA
Telephone (800) 325-5832 (314) 771-5765
Fax (314) 286-7828
email: techserv@sial.com
sigma-aldrich.com

Product Information

14,15-EPOXY-(5Z,8Z,11Z)- EICOSATRIENOIC ACID (14,15-EET)

Product No. **E 5891**

SPECIAL HANDLING INSTRUCTIONS

All glassware and stainless steel syringe needles to be used in the transfer of 14,15-epoxyeicosatrienoic acid solution should be prerinsed with absolute ethanol and air dried. Transfer and removal of samples should be carried out under an inert atmosphere, keeping the bulk at 0 °C or below whenever possible.

14,15-Epoxyeicosatrienoic acid is unstable to oxygen and direct sunlight, therefore extreme care must be taken in handling this product.

Pcs3/00

Sigma brand products are sold through Sigma-Aldrich, Inc.
Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications.
Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply.
Please see reverse side of the invoice or packing slip.