

Product Information

SIGMAFAST™ BCIP®/NBT Tablet

B5655

Product Description

5-bromo-4-chloro-3-indolyl phosphate (BCIP®) and nitro blue tetrazolium (NBT) are reagents that are widely used in tandem for detection of alkaline phosphatase conjugates.¹⁻³ SIGMAFAST™ BCIP®/NBT tablets are useful in immunochemistry to detect alkaline phosphatase. Common uses are in Western blotting or dot blotting, and to a lesser extent in immunohistology.

SIGMAFAST™ BCIP®/NBT tablets require no additional buffers or steps to prepare an active substrate solution. One tablet, dissolved in 10 mL of water, provides 10 mL of ready-to-use buffered substrate solution. The substrate solution contains:

- BCIP (0.15 mg/mL)
- NBT (0.30 mg/mL)
- Tris buffer (100 mM)
- MgCl₂ (5 mM)

The solution pH is in the range of 9.25 – 9.75.

Several theses⁴⁻⁶ and dissertations⁷⁻¹⁶ have cited use of product B5655 in their protocols.

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Storage/Stability

Store the tablets at -20 °C.

Preparation Instructions

- Remove the tablet package from the freezer. Allow it to warm to room temperature.
- Open the foil pack. Drop the SIGMAFAST™ BCIP®/NBT tablet into an appropriate container. **Do not touch the tablet with your fingers.**
- Add 10 mL of water. Vortex until dissolved (2-5 minutes).

The SIGMAFAST™ BCIP®/NBT Substrate Solution is now ready for use. For best results, the Substrate Solution should be used within one hour.

Procedure

1. Pour the SIGMAFAST™ BCIP®/NBT Substrate Solution into a suitable container.
2. Lay the nitrocellulose paper in the solution. Make sure the paper is completely covered with the Substrate Solution.
3. Remove the nitrocellulose paper when sufficient color has developed (5-10 minutes).
4. Rinse in water. Although several blots can be developed in this manner, the potential for carryover from blot to blot does exist and should be evaluated carefully.
5. Blots stained with SIGMAFAST™ BCIP®/NBT Substrate Solution may be dried and stored away from light for future reference.

Troubleshooting

Background is too high

- Use a blocking step prior to application of the primary antibody. Normal serum (10% v/v) from the same species as the second antibody generally produces the best results.
- Additional blocking agents for immunoblotting include:
 - 10% BSA
 - 0.05% TWEEN® 20
 - 3% nonfat-dried milk
 - Do not use milk as a blocking agent when using avidin-biotin systems.
- Decrease staining time.
- Titer the conjugate to optimize the working dilution.

No color develops or color is too faint

- Adjust the concentration of the primary antibody.
- Adjust the concentration of the secondary antibody.

- Determine if the enzyme conjugate is active.
- Consider using an amplifying system such as avidin-biotin.
- Increase the staining time.
- Adjust the transfer time of the samples to the nitrocellulose membrane.
- Increase the amount of sample.

References

1. Blake, M.S. *et al.*, *Anal. Biochem.*, **136(1)**, 175-179 (1984).
2. Horowitz, J.P. *et al.*, *J. Med. Chem.*, **9(3)**, 447 (1966).
3. Smejkal, G.B., and Kaul, C.A., *J. Histochem. Cytochem.*, **49(9)**, 1189-1190 (2001).
4. Duench, Kellie, "Unravelling the Induction and Development of Intramembranous Bones in the Chicken Eye". Saint Mary's University, M.Sc. thesis, p. 90 (2010).
5. Muller, Manuel, "Enzymatic and enzyme inhibitory activity on a paper-based lateral flow device". University of Rhode Island, M.S. thesis, p. 40 (2015).
6. Fields, Christopher J., "Functional and Expression Analysis of a Novel Basement Membrane Degrader in *Drosophila melanogaster*". Western Kentucky University, M.S. thesis, p. 26 (2016).
7. Harvey, Lucy, "Exploring the *in vitro* behaviour of endothelial cells in different cell culture models". University of Bristol, M.Res thesis, p. 49 (2018).
8. Moragoda, Lathika Pinto, "Characterization of the expression of two maize root-preferential genes at the protein level". Iowa State University, Ph.D. dissertation, p. 69 (1997).
9. Williams, Bryony A.P., "Molecular evolution and the fate of the mitochondrion in microsporidia parasites". University College London, Ph.D. dissertation, p. 77 (2003).
10. Chang, Li-Wen, "Development of molecular diagnostic tests for inclusion body disease in boid snakes". University of Florida, Ph.D. dissertation, p. 57 (2012).
11. Arulmoli, Janahan, "Scaffolds for Neural Stem Cell Tissue Engineering". University of California Irvine, Ph.D. dissertation, p. 97 (2016).
12. Brod, Florian, "Identification and assessment of antigens involved in membrane traversal by malaria parasites as vaccine candidates". University of Oxford, Ph.D. dissertation, p. 59 (2018).
13. Philipson, Alice Rae, "Towards Rapid Label-free Enrichment of Specific Stem Cell Populations for Autologous Cell Therapies". University of Leeds, Ph.D. dissertation, p. 125 (2019).
14. Rieger, Jan Hendrik, "Targets of tumor-infiltrating lymphocyte reactivity in pancreatic cancer". Ruperto Carola University Heidelberg, Dr. rer. nat. dissertation, p. 26 (2019).
15. Bullen, Louise Frances, "An interrogation into the role of p62 in the pluripotent state". Manchester Metropolitan University, Ph.D. dissertation, p. 70 (2020).
16. Visconti, Richard, "A Functional Three-Dimensional Microphysiological Model of Myeloma Bone Disease". Seton Hall University, Ph.D. dissertation, p. 17 (2020).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Standard Warranty

The applicable warranty for the products listed in this publication may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck KGaA, Darmstadt, Germany operates as MilliporeSigma in the U.S. and Canada.

MilliporeSigma, SIGMAFAST, BCIP, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2009-2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

B5655dat Rev 08/22 RBG,GCY,MAM

**MILLIPORE
SIGMA**