

3050 Spruce Street Saint Louis, Missouri 63103 USA Telephone (800) 325-5832 (314) 771-5765 Fax (314) 286-7828 email: techserv@sial.com sigma-aldrich.com

ProductInformation

L-Proline Cell culture tested, meets EP & USP testing specifications

Product Number P 5607

Product Description

Molecular Formula: $C_5H_9NO_2$ Molecular Weight: 115.1 CAS Number: 147-85-3 pK_a : 1.95 (COOH), 10.64 $(NH_2)^1$ pl: 6.3¹ Synonyms: (S)-Pyrrolidine-2-carboxylic acid, Pro

This product is cell culture tested (0.03 mg/ml) and is tested for endotoxin levels.

The amino acid L-Proline is unique among the 20 amino acids normally found in proteins because it is the only imino acid among them.² The side chain is bonded to both the α -carbon and nitrogen atoms, giving a secondary amino group in the structure, and thus its characterization as an imino acid. *In vivo*, proline is both synthesized from and degraded to glutamate, with glutamate γ -semialdehyde and Δ^1 -pyrroline-5-carboxylate as intermediates, by separate pathways.^{3,4}

In protein structure, collagen and elastin are notably rich in proline content. In collagen in particular, the proline ring structure helps to stabilize the collagen α -helix conformation.⁵ The cis-trans isomerization of proline residues in ribonuclease A has been studied by fluorescence methods.⁶

L-Proline is used in cell culture media and is a component of MEM non-essential amino acids solution (Product No. M 7145). In cell culture research, L-proline has been used in studies of MOLT-4 cells in starvation-induced apoptosis, maintenance of hematopoietic cell cultures derived from the prawn *Nephrops norvegicus*, growth behavior of various *Escherichia coli* strains at various L-proline concentrations, and rat primary hepatocytes in a collagen gel sandwich culture.^{7,8,9,10}

A procedure for the asymmetric self-aldolization of acetaldehyde that uses L-proline as a catalyst has been reported.¹¹ Ground- and transition-state analogue inhibitors of cyclophilin have been synthesized using L-proline as the starting compound.¹² The use of L-proline to synthesize an optically active 1-oxo-2-oxa-5-azaspiro[3.4]octane compound, related to the antibiotic oxazolomycin, has been described.¹³

Precautions and Disclaimer

For Laboratory Use Only. Not for drug, household or other uses.

Preparation Instructions

This product is soluble in water (50 mg/ml), yielding a clear, colorless solution.

References

- Molecular Biology LabFax, Brown, T. A., ed., BIOS Scientific Publishers Ltd. (Oxford, UK: 1991), p. 29.
- 2. The Merck Index, 12th ed., Entry# 7963.
- 3. Biochemistry, 3rd ed., Stryer, L., W. H. Freeman (New York, NY: 1988), pp. 18, 505, 579.
- Textbook of Biochemistry with Clinical Correlations, 5th ed., Devlin, T. M., ed., Wiley-Liss (New York, NY: 2002), pp. 97, 792.
- 5. Molecular Biology of the Cell, 3rd ed., Alberts, B., et al., ed., Garland Publishing (New York, NY: 1994), pp. 115, 979, 984.
- Bhat, R., et al., Proline isomerization in bovine pancreatic ribonuclease A. 2. Folding conditions. Biochemistry, **42(19)**, 5722-5728 (2003).
- Franek, F., et al., Antiapoptotic and proapoptotic action of various amino acids and analogs in starving MOLT-4 cells. Arch. Biochem. Biophys., 398(1), 141-146 (2002).
- Mulford, A. L., et al., Development and characterization of primary cell cultures from the hematopoietic tissues of the Dublin Bay prawn, *Nephrops norvegicus*. Methods Cell Sci., 22(4), 265-275 (2000).

- Rothen, S. A., et al., Growth characteristics of Escherichia coli HB101[pGEc47] on defined medium. Biotechnol. Bioeng., 58(1), 92-100 (1998).
- Lee, J., et al., Proline-mediated enhancement of hepatocyte function in a collagen gel sandwich culture configuration. FASEB J., 7(6), 586-591 (1993).
- Cordova, A., et al., Proline-catalyzed one-step asymmetric synthesis of 5-hydroxy-(2E)-hexenal from acetaldehyde. J. Org. Chem., 67(1), 301-303 (2002).
- Wang, H. C., et al., Structure-activity studies of ground- and transition-state analogue inhibitors of cyclophilin. J. Med. Chem., 44(16), 2593-2600 (2001).
- Papillon, J. P., and Taylor, R. J., The first syntheses of the 1-oxo-2-oxa-5azaspiro[3.4]octane ring system found in oxazolomycin. Org. Lett., 2(14), 1987-1990 (2000).

GCY/RXR 11/05

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.