

P493-6 Human Lymphoblastoid Cell Line

Cancer Cell Line
Cat. # SCC279

Pack size: $\geq 1 \times 10^6$
viable cells/vial

FOR RESEARCH USE ONLY.
NOT FOR USE IN DIAGNOSTIC PROCEDURES.
NOT FOR HUMAN OR ANIMAL CONSUMPTION.
THIS PRODUCT CONTAINS GENETICALLY MODIFIED ORGANISMS.

Store in liquid nitrogen



Data Sheet

page 1 of 3

Background

The proto-oncogene *c-myc* is an activator of cell proliferation, promoting cell cycle progression, and is thus significant factor in many cancers. Burkitt's lymphoma is a rare cancer of human B cells associated with Epstein-Barr virus (EBV) that is the most rapidly growing human tumor, due to transcriptional activation of *c-myc* by chromosomal translocation.¹ Because of its highly aggressive nature, Burkitt's lymphoma has a high mortality rate if left untreated. Cellular models of *c-myc*-activated lymphomas are essential for understanding mechanisms of proliferation and for developing treatments for this disease.

The P493-6 Human Lymphoblastoid Cell Line expresses *c-myc* under control of a tetracycline-inducible promoter.² P493-6 cells constitutively express *c-myc* when cultured without tetracycline. Addition of tetracycline switches off *c-myc* expression and arrests cells in G0/G1 phase. P493-6 cells demonstrate the morphology and surface antigen expression typical of Burkitt lymphoma cells, including CD10 and CD38.³ The P493-6 B cell line permits conditional *myc* expression allowing for detailed studies of proliferative mechanisms in B cell lymphoma.

Source

The P493-6 cell line was established from a clone of the EREB2-5 EBV-positive lymphoblastoid cell line stably transfected with a *c-myc-tet* plasmid construct.²

Short tandem repeat (STR) Profile

D3S1358: 15, 18	D16S539: 12
TH01: 8, 9.3	CSF1PO: 9, 12
D21S11: 28, 32.2	Penta D: 11, 13
D18S51: 13	vWA: 14, 18
Penta E: 12, 14	D8S1179: 9, 13
D5S818: 11, 12	TPOX: 8
D13S317: 11, 12	FGA: 22
D7S820: 8	Amelogenin: X

Cancer cell lines are inherently genetically unstable. Genetic instability may arise in the form of loss of heterozygosity of alleles at one or more genetic sites with increased passages.

Storage & Handling

P493-6 Human Lymphoblastoid Cell Line should be stored in liquid nitrogen. The cells can be cultured for at least 10 passages after initial thawing without significantly affecting the cell marker expression and functionality.

Quality Control Testing

- Each vial contains $\geq 1 \times 10^6$ viable cells.
- Cells are tested negative for infectious diseases by a Human Essential CLEAR panel by Charles River Animal Diagnostic Services.
- Cells are verified to be of human origin and negative for inter-species contamination from rat, mouse, chinese hamster, Golden Syrian hamster, and non-human primate (NHP) as assessed by a Contamination CLEAR panel by Charles River Animal Diagnostic Services.
- Cells are negative for mycoplasma contamination.
- Each lot of cells is genotyped by STR analysis to verify the unique identity of the cell line.

Representative Images

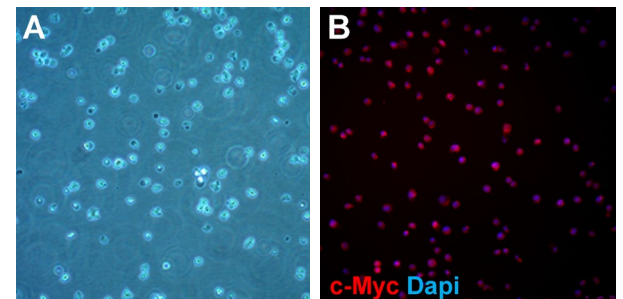


Figure 1. Brightfield image of cells one day after thaw (A). Cells express *c-myc* (B, MAB8864) under control of a tetracycline-inducible promoter.

References

1. *Lancet*. 2012; 379(9822): 1234-1244.
2. *Int J Cancer*. 2000; 87(6): 787-793.
3. *Int J Cancer*. 2001; 93(6): 810-816

Please visit www.milliporesigma.com for additional product information and references.

Submit your published journal article and earn credit toward future purchases. Visit www.milliporesigma.com/publicationrewards to learn more!

P493-6 Human Lymphoblastoid Cell Line

Cat # SCC279

Protocols

P493-6 Human Lymphoblastoid Cell Line grows as suspension cells and thus do not require enzymatic detachment or dissociation. Passage when the cell density reaches 1–1.5 million cells/mL. Optimal plating density should be ~200,000 - 250,000 cells/mL. The cells should not be grown at excessively high densities.

1. Do not thaw the cells until the recommended medium is on hand.

P493-6 Expansion Medium: Cells are thawed and expanded in RPMI-1640 (Sigma Cat. No. R0883) supplemented with 2 mM Glutamine (Cat. No. TMS-002-C) and 10% FBS (Cat. No. ES-009-B).

2. Remove the vial of frozen P493-6 cells from liquid nitrogen and incubate in a 37°C water bath. Closely monitor until the cells are completely thawed. Maximum cell viability is dependent on the rapid and complete thawing of frozen cells.

IMPORTANT: Do not vortex the cells.

3. As soon as the cells are completely thawed, disinfect the outside of the vial with 70% ethanol. Proceed immediately to the next step.

4. In a laminar flow hood, use a 1 or 2 mL pipette to transfer the cells to a sterile 15 mL conical tube. Be careful not to introduce any bubbles during the transfer process.

5. Using a 10 mL pipette, slowly add dropwise 9 mL of P493-6 Expansion Medium (Step 1 above) to the 15 mL conical tube.

IMPORTANT: Do not add the entire volume of media all at once to the cells. This may result in decreased cell viability due to osmotic shock.

6. Gently mix the cell suspension by slowly pipetting up and down twice. Be careful not to introduce any bubbles.

IMPORTANT: Do not vortex the cells.

7. Centrifuge the tube at 300 x g for 2-3 minutes to pellet the cells.

8. Decant as much of the supernatant as possible. Steps 5-8 are necessary to remove residual cryopreservative (DMSO).

9. Resuspend the cells in 15 – 20 mL of P493-6 Expansion Medium.

10. Transfer the cell suspension to a T75 flask.

11. Incubate the cells at 37°C in a humidified incubator with 5% CO₂. P493-6 suspension cells require media replenishment every 2-3 days. Passage cells when the cell density is at 1 -1.5 million cells/mL.

12. Cells are typically plated at a density of 200,000 - 250,000 cells/mL

Cryopreservation of Cells

P493-6 Human Lymphoblastoid Cell Line may be frozen in the expansion medium plus 10% DMSO using a Nalgene slow freeze Mr. Frosty container.

■ antibodies ■ Multiplex products ■ biotools ■ cell culture ■ enzymes ■ kits ■ proteins/peptides ■ siRNA/cDNA products

Please visit www.milliporesigma.com for additional product information, test data and references

MilliporeSigma, 28820 Single Oak Drive, Temecula, CA 92590, USA 1-800-437-7500

Technical Support: T: 1-800-MILLIPORE (1-800-645-5476) • F: 1-800-437-7502

FOR RESEARCH USE ONLY. Not for use in diagnostic procedures. Not for human or animal consumption. Purchase of this Product does not include any right to resell or transfer, either as a stand-alone product or as a component of another product. Any use of this Product for purposes other than research is strictly prohibited.

EMD Millipore®, the M mark, Upstate®, Chemicon®, Linco® and all other registered trademarks, unless specifically identified above in the text as belonging to a third party, are owned by Merck KGaA, Darmstadt, Germany. Copyright ©2008-2020 Merck KGaA, Darmstadt, Germany. All rights reserved.



We Buy 100% Certified Renewable Energy

ACADEMIC USE AGREEMENT
(subject to local law)

THIS PRODUCT MAY ONLY BE USED BY INDIVIDUALS EMPLOYED BY AN ACADEMIC INSTITUTION AND IS INTENDED SOLELY TO BE USED FOR ACADEMIC RESEARCH, WHICH IS FURTHER DEFINED BELOW. BY OPENING THIS PRODUCT, YOU (“PURCHASER”) HEREBY REPRESENT THAT YOU HAVE THE RIGHT AND AUTHORITY TO LEGALLY BIND YOURSELF AND/OR YOUR EMPLOYER INSTITUTION, AS APPLICABLE, AND CONSENT TO BE LEGALLY BOUND BY THE TERMS OF THIS ACADEMIC USE AGREEMENT. IF YOU DO NOT AGREE TO COMPLY WITH THESE TERMS, YOU MAY NOT OPEN OR USE THE PRODUCT AND YOU MUST CALL MILLIPORESIGMA (“SELLER”) CUSTOMER SERVICE (1-800-645-5476) TO ARRANGE TO RETURN THE PRODUCT FOR A REFUND.

“Product” means P493-6 Human Lymphoblastoid Cell Line (SCC279)

“Academic Research” means any internal *in vitro* research use by individuals employed by an academic institution. Academic Research specifically excludes the following uses of whatever kind or nature:

- Re-engineering or copying the Product
- Making derivatives, modifications, or functional equivalents of the Product
- Obtaining patents or other intellectual property rights claiming use of the Product
- Using the Product in the development, testing, or manufacture of a Commercial Product
- Using the Product as a component of a Commercial Product
- Reselling or licensing the Product
- Using the Product in clinical or therapeutic applications including producing materials for clinical trials
- Administering the Product to humans
- Using the Product in collaboration with a commercial or non-academic entity

“Commercial Product” means any product intended for: (i) current or future sale; (ii) use in a fee-for-service; or (iii) any diagnostic, clinical, or therapeutic use.

Access to the Product is limited solely to those officers, employees, and students of PURCHASER’s academic institution who need access to the Product to perform Academic Research. PURCHASER shall comply with all applicable laws in its use and handling of the Product and shall keep it under reasonably safe and secure conditions to prevent unauthorized use or access.

These use restrictions will remain in effect for as long as PURCHASER possesses the Product.

COMMERCIAL OR NON-ACADEMIC ENTITIES INTERESTED IN PURCHASING OR USING THE PRODUCT MUST CONTACT licensing@emdmillipore.com AND AGREE TO SEPARATE TERMS OF USE PRIOR TO USE OR PURCHASE.

GMO

This product contains genetically modified organisms.
Este producto contiene organismos genéticamente modificados.
Questo prodotto contiene degli organismi geneticamente modificati.
Dieses Produkt enthält genetisch modifizierte Organismen.
Ce produit contient des organismes génétiquement modifiés.
Dit product bevat genetisch gewijzigde organismen.
Tämä tuote sisältää geneettisesti muutettuja organismeja.
Denna produkt innehåller genetiskt ändrade organismer.

■ antibodies ■ Multiplex products ■ biotools ■ cell culture ■ enzymes ■ kits ■ proteins/peptides ■ siRNA/cDNA products

Please visit www.milliporesigma.com for additional product information, test data and references

MilliporeSigma, 28820 Single Oak Drive, Temecula, CA 92590, USA 1-800-437-7500

Technical Support: T: 1-800-MILLIPORE (1-800-645-5476) • F: 1-800-437-7502

FOR RESEARCH USE ONLY. Not for use in diagnostic procedures. Not for human or animal consumption. Purchase of this Product does not include any right to resell or transfer, either as a stand-alone product or as a component of another product. Any use of this Product for purposes other than research is strictly prohibited.

EMD Millipore®, the M mark, Upstate®, Chemicon®, Linco® and all other registered trademarks, unless specifically identified above in the text as belonging to a third party, are owned by Merck KGaA, Darmstadt, Germany. Copyright ©2008-2020 Merck KGaA, Darmstadt, Germany. All rights reserved.



We Buy 100% Certified Renewable Energy