

Product Information

STRIP-FUGE™



OPERATING INSTRUCTIONS

Catalog #C5980 – Mini-centrifuge – 220V
Catalog #C5855 – Mini-centrifuge – 110V

- Place the Strip-Fuge on a flat, clear surface.
- To open, press firmly down on release tab while lifting dome.
- Close dome and activate side switch. Device will not operate with dome open.
- To stop, deactivate side switch. Wait until completely stopped to release dome. Remove samples.



Loading *The BUTTERFLY Rotor*™ - 8 x 200 µl Tube Strips

- Load all tubes on both strips with equal amounts of sample. Close the dome securely.
- Insert strips into rotor. If necessary, add a strip of dummy-tubes to provide balance.

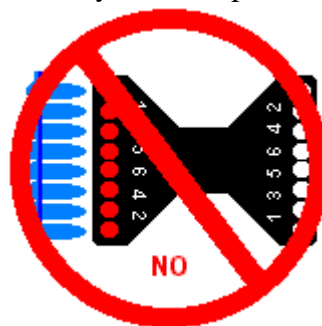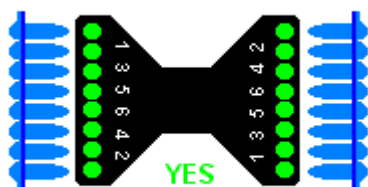


Loading *The BUTTERFLY Rotor*™ - 200 µl Single Tubes

- Load tubes with equal amounts of sample in opposing positions as indicated on the rotor.

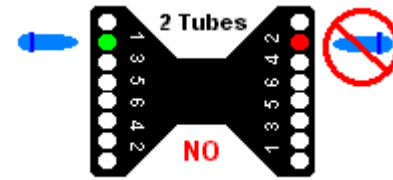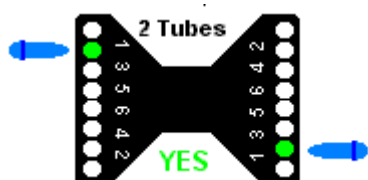

- Never place single tubes in corner positions. Your unit has been designed to minimize footprint. Single tubes may make contact with the dome and cause damage if placed in corner positions.

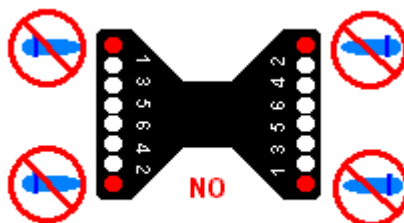

Safeguard

- Do not immerse device in water. Injury due to shock or fire may result.
- Do not operate device with dome open, or open dome before rotor has ceased spinning.
- Do not insert non-standard tubes or other foreign articles into rotor.
- Do not operate under uneven loads. Destabilized operation may result in damage or injury.
- Store unit at room temperature in a dry area. Avoid sun, moisture or extreme temperatures.

Sigma-Aldrich • P.O. Box 14508 • St. Louis, Missouri 63178

Phone: (800) 325-3010 • Fax: (800) 325-5052 • www.sigma-aldrich.com

Our one-year warranty assures you reliable quality at an unbeatable value!