

Technical Data Sheet

Test Agar pH 6.0 for the Inhibitor Test

Ordering number: 1.10663.0500

For the detection of antimicrobial inhibitors in meat and organ samples together with *Bacillus subtilis* (BGA) spore suspension and *Micrococcus luteus* ATCC 9341 as test organisms.

The nutrient media are suitable both, for the inhibitor test (LEVETZOW, 1971) according to the German Meat Inspection Law as well as for the EEC Four-Plate-Test (BOGAERTS and WOLF, 1980) suggested by the Scientific Veterinary Commission of the European Economic Community.

Test agar pH 7.2 with addition of trimethoprim is used particularly for determination of sulfonamide residues.

Mode of Action

The test is carried out according to the agar diffusion procedure. Small slices of the meat sample are placed on the inoculated Test Agar plates and incubated. Antimicrobial inhibitors contained in the samples diffuse into the nutrient media and cause growth-free inhibition zones to develop on the otherwise thickly covered plates. Repeated tests with pH 6.0, pH 8.0 and pH 7.2 are necessary, as penicillin and streptomycin are optimally active at pH 6.0 and 8.0 respectively (PICHNARCIK et al., 1969) and the activity optimum of sulfonamide is found at pH 7.2. Addition of trimethoprim to Test Agar pH 7.2 considerably increases the sensitivity of the test system to sulfonamides (GUDDING, 1976; EBRECHT, 1982).

Typical Composition (g/L)

Test Agar pH 6.0 for the Inhibitor Test	
Peptone from casein, tryptic	3.45
Peptone from meat, tryptic	3.45
Sodium chloride	5.1
Agar-agar	13.0

**Agar-agar is equivalent to other different terms of agar.

Preparation

Suspend 25 g/litre Test Agar pH 6.0, autoclave (15 min at 121°C), test the pH and if necessary adjust. Cool to 50-45 °C, mix in 1ml/litre *Bacillus subtilis* (BGA) spore suspension.

After mixing the spore suspension, immediately pour the plates and place in the refrigerator.

The plates are clear and yellowish-brown.

Experimental Procedure and Evaluation

Inhibitor Test

Test Agar pH 6.0 with *Bacillus subtilis* (BGA).

EEC Four-Plate-Test

Test Agar pH 6.0 with *Bacillus subtilis* (BGA).

For details regarding the collection of samples, transportation as well as the execution of the test, see the Fleischbeschauengesetz (Meat Inspection Law) or BOGAERTS and WOLF (1980).

According to these specifications the cylinder-shaped tissue sections (8 mm in diameter and 2 mm thick) are stamped out under conditions, which are as aseptic as possible and laid on the plates; according to BOGAERTS and WOLF two sections are required per plate. As a control, one test disc with 10 IU of penicillin G-sodium is placed on a plate with pH 6.0, one test disc with 10 µg of streptomycin on each of the two plates with pH 8.0 and one test disc with 0.5 µg sulfadimidine on a plate with pH 7.2. The test discs can also be made by the user using filter paper discs, 6 mm in diameter.

Incubation: 18-24 hours at 30 °C (*Bacillus subtilis* BGA).

The inhibition zones between the edge of the tissue section or the test disc and the growth limit of the test organism are measured. Complete inhibition of growth with an inhibitory zone of at least 2 mm is regarded as positive result, an inhibitory zone of 1-2 mm should be considered as doubtful. This only applies, however, if the controls, prepared at the same time, display zones of inhibition measuring about 6 mm.

For possible methodological improvements see FORSCHNER and SEIDLER (1976).

Storage

Store at +15 °C to +25 °C, dry and tightly closed. Do not use clumped or discolored medium. Protect from UV light (including sun light). For *in vitro* use only.

The ready-to-use plates can be sealed with air-tight adhesive tape and, when cooled (+4 to +6 °C) may be kept for up to 2 weeks. Additional packing into plastic bags is recommended. If stored for longer periods, the temperature should not be more than +3 °C; however, freezing of the culture medium must be avoided



We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

Merck Millipore, the M logo, are registered trademarks of Merck KGaA, Darmstadt, Germany. ATCC is a registered trademark of ATCC, Manassas, VA, USA. Lit. No. TN6093EN00

Quality Control

Parameter	Diameter
Inhibition zone diameter Gentamicin 10 µg Bacillus subtilis BGA	20-28 mm
Inhibition zone diameter Gentamicin 30 µg; Bacillus subtilis BGA	22-30 mm
Inhibition zone diameter Penicillin G 10 U/IE Bacillus subtilis BGA	36-48 mm
Inhibition zone diameter Streptomycin 10 µg Bacillus subtilis BGA	19-27 mm

Please refer to the actual batch related Certificate of Analysis.



Gentamicin 30µg



Penicillin 10IE



We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

Merck Millipore, the M logo, are registered trademarks of Merck KGaA, Darmstadt, Germany. ATCC is a registered trademark of ATCC, Manassas, VA, USA. Lit. No. TN6093EN00

Literature

Arbeitsgruppe des Instituts für Lebensmitteltechnologie und Verpackung der Technischen Universität München: Merkblätter für die Prüfung von Packmitteln, Merkblatt 18 "Prüfung auf antimikrobielle Bestandteile in Packstoffen". - **Verpackungs.-Rundschau 25**; Techn.-wiss. Beilagen; 5-8 (1974).

BAUR, E.: Untersuchungen von Fleisch- und Wurstwaren mit dem Hemmstofftest im Rahmen der tierärztlichen Lebensmittelüberwachung. - **Fleischwirtschaft 55**; 843-845 (1975)

BOGAERTS, R., u. WOLF, F.: Eine standardisierte Methode zum Nachweis von Rückständen antibakteriell wirksamer Substanzen in frischem Fleisch. - **Fleischwirtschaft 60**; 667-675 (1980).

Deutsches Fleischbeschaugesetz: Ausführungsbestimmungen A über die Untersuchung und gesundheitspolizeiliche Behandlung der Schlachttiere und des Fleisches bei Schlachtungen im Inland; Anlage 4 zu § 20 Abs. 4: Rückstandsuntersuchung.

EBRECHT, A.: Verbesserung des Hemmstofftestes durch Zusatz von Trimethoprim zum Nachweis von Sulfonamiden. - **Arch. Lebensmittelhyg. 33**; 109-115 (1982).

FORSCHNER, E., u. SEIDLER, M.: Alternativvorschläge zum Hemmstofftest. Rationalisierung und Absicherung. - **Fleischwirtschaft 56**; 1008-1013 (1976).

GUDDING, R.: An improved bacteriological method for the detection of sulfonamide residues in food. - **Acta Vet. Scand. 17**; 458-464 (1976).

LEVETZOW, R.: Untersuchungen auf Hemmstoffe im Rahmen der Bakteriologischen Fleischuntersuchung (BU). - **Bundesgesundheitsblatt 14**; 15/16, 211-213 (1971).

PICHNARCIK, J., WENZEL, S., u. GISSKE, W.: Beitrag zur Methodik des Hemmstoffnachweises in Organen und Muskulatur von Schlachttieren. - **Arch. Lebensmittelhyg. 20**; 272-279 (1969).

Ordering Information

Product	Cat. No.	Pack size
Test Agar pH 6.0 for the Inhibitor Test	1.10663.0500	500 g
Bacillus subtilis (BGA)-spore suspension	1.10649.0001	15 x 2 ml
Penicillin G, Potassium Salt	Calbiochem 5161-25MU	25 MU
Gentamycin sulfate	Calbiochem 345814-1GM	1 g
Streptomycin sulfate	Calbiochem 5711-100GM	100 g

Merck KGaA, 64271 Darmstadt, Germany
Fax: +49 (0) 61 51 / 72-60 80
mibio@merckgroup.com
www.merckmillipore.com/biomonitoring

Find contact information for your country at:
www.merckmillipore.com/offices
For Technical Service, please visit:
www.merckmillipore.com/techservice



We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

Merck Millipore, the M logo, are registered trademarks of Merck KGaA, Darmstadt, Germany. ATCC is a registered trademark of ATCC, Manassas, VA, USA. Lit. No. TN6093EN00