


MC-Media Pad® Rapid Aerobic Count Test


Interpretation Guide

The MC-Media Pad® is a convenient method for indicator organism testing. Compliant with international food and beverages standards such as AOAC-PTM, AOAC-OMA and ISO 16140 (MicroVal), MC-Media Pad® tests streamline your microbial quality control workflow.


Accurate results in 4 steps

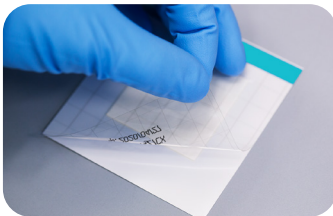


1. Open the cover film diagonally

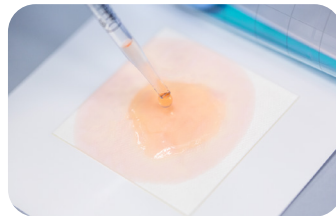

2. Inoculate your sample, it will diffuse automatically

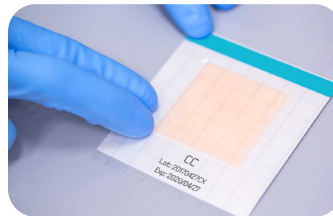

3. Close the cover film and incubate

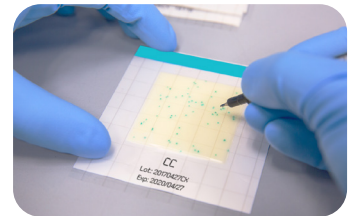

4. Count your results



Staphylococcus aureus
ATCC® 6538™ 50 CFUs



Staphylococcus aureus
ATCC® 6538™ 150 CFUs




Staphylococcus aureus
ATCC® 6538™ 300 CFUs



Escherichia coli ATCC®
6538™ 80 CFUs

With MC-Media Pad[®], it is easy to see when a dilution is needed.

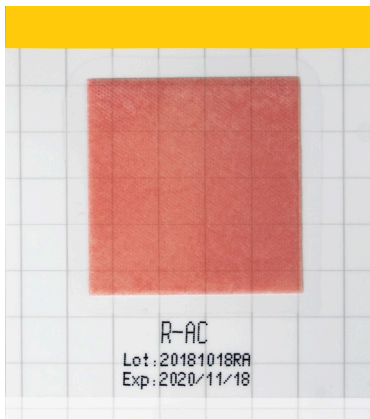

Dilution for easier read-out




Dilution factor 1:10



Dilution for easier read-out



Dilution factor 1:10

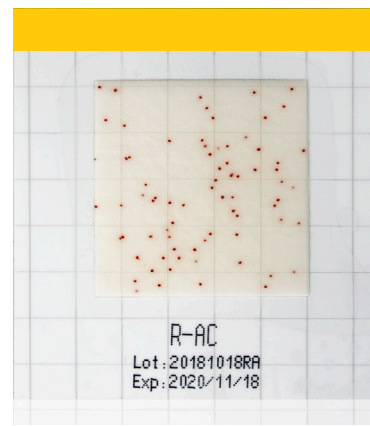

Ordering information	Product name	Packaging
1.32359.001	MC-Media Pad [®] RAC (Rapid Aerobic Count)	100 pcs

Please visit: SigmaAldrich.com/MC-MediaPad

