Product Information

sigma-aldrich.com

3050 Spruce Street, Saint Louis, MO 63103 USA Tel: (800) 521-8956 (314) 771-5765 Fax: (800) 325-5052 (314) 771-5757 email: techservice@sial.com sigma-aldrich.com

Arabinogalactan from larch wood

for density gradient separations

Catalog Number **A6606** Store at Room Temperature

CAS RN 9036-66-2 Synonym: Polyarabinogalactan

Product Description

In the early seventies a crude form of arabinogalactan was purified for use in density gradient separation of red blood cells.¹ Many researchers have since found arabinogalactan to be an appropriate medium for use in cell separations.

This product is a highly purified arabinogalactan prepared specifically for use in density gradient separations of cells, viruses, and organelles. Arabinogalactan is nontoxic, completely compatible with biological materials, readily soluble in water, and can be sterilized by filtration or autoclaving. Arabinogalactan solutions have low viscosities. They are not sensitive to changes in pH or dissolved salts, and are readily adjusted to the desired osmolality.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

Preparation Instructions

General Density Gradient Solution [28.5% (w/v) Arabinogalactan (Catalog Number A6606) solution in water] - Heat as needed to dissolve and cool to room temperature. Bring to final volume with water.

Density: 1.120±0.002 g/ml

General Density Gradient Diluent - May also be used as diluent for solution for platelet or lymphocyte separation.

KCI (Catalog Number P4504)	28 g/L
Na ₂ HPO ₄ (Catalog Number S0876)	1.22 g/L
KH ₂ PO ₄ (Catalog Number P5379)	0.22 g/L
D(+)-Glucose (Catalog Number G8270)	2.00 g/L

Heat as needed to dissolve and cool to room temperature. If necessary, adjust the pH to 7.3-7.5 prior to bringing to final volume with water.

Density: 1.040±0.002 g/ml Osmolality: 295±5 mOsm/kg

Isotonic Arabinogalactan Solution -	
NaCI (Catalog Number S9625)	5.50 g/L
Na ₂ HPO ₄ (Catalog Number S0876)	1.22 g/L
KH ₂ PO ₄ (Catalog Number P5379)	0.22 g/L
Arabinogalactan (Catalog Number A6606)	29.5%

Heat as needed to dissolve and cool to room temperature. Bring to final volume with water. Density: 1.120±0.002 g/ml Osmolality: 295±5 mOsm/kg

Heat as needed to dissolve and cool to room temperature. If necessary, adjust the pH to 7.3-7.5 prior to bringing to final volume with water. Density: 1.040±0.002 g/ml

Solution for Platelet Separation -	
NaCI (Catalog Number S9625)	4.9 g/L
Na ₂ HPO ₄ (Catalog Number S0876)	1.20 g/L
KH ₂ PO ₄ (Catalog Number P5379)	0.20 g/L
Sodium citrate tribasic dihydrate	4.00 g/L
(Catalog Number C7254)	-
D(+)-Glucose (Catalog Number G8270)	2.00 g/L
Arabinogalactan (Catalog Number A6606)	20.0%

Heat as needed to dissolve and cool to room temperature. If necessary, adjust the pH to 7.3-7.5 prior to bringing to final volume with water.

Density: 1.082±0.001 g/ml Osmolality: 290±5 mOsm/kg



Solution for Lymphocyte Separation -	
NaCl (Catalog Number S9625)	6.8 g/L
KCI (Catalog Number P4504)	0.40 g/L
Sodium bicarbonate	0.35 g/L
(Catalog Number S8875)	
Na ₂ HPO ₄ (Catalog Number S0876)	0.048 g/L
KH ₂ PO ₄ (Catalog Number P5379)	0.060 g/L
D(+)-Glucose (Catalog Number G8270)	1.0 g/L
Arabinogalactan (Catalog Number A6606)	18.5%

Heat as needed to dissolve and cool to room

temperature. If necessary, adjust the pH to 7.3-7.5 prior to bringing to final volume with water.

Density: 1.077±0.001 g/ml Osmolality: 290±5 mOsm/kg

Storage/Stability

Store the product at room temperature. When stored at room temperature, the arabinogalactan powder is stable indefinitely.

Stock solutions are stable for at least one year at -20 °C. Filter the solutions through a 0.45 μ m filter and store in 50 ml aliquots in capped containers at -20 °C.

References

The general references cite uses of arabinogalactan in cell research.

- Corash, L.M., et al., Separation of erythrocytes according to age on a simplified density gradient. *J. Lab. Clin. Med.*, 84, 147-51 (1974).
- Corash, L.M., et al., Heterogeneity of human whole blood platelet subpopulations. I. Relationship between buoyant density, cell volume, and ultrastructure. *Blood*, **49**, 71-87 (1977).
- Corash, L.M., et al., Heterogeneity of human whole blood platelet subpopulations. II. Use of a subhuman primate model to analyze the relationship between density and platelet age. *Blood*, **52**, 726-84 (1978).

- 4. Bunn, H.F., et al., Molecular and cellular pathogenesis of hemoglobin SC disease. *Proc. Natl. Acad. Sci. USA*, **79**, 7527-31 (1982).
- Clark, M.R., et al., Study on the dehydrating effect of the red cell Na⁺/K⁺-pump in nystatin-treated cells with varying Na⁺ and water contents. *Biochim. Biophys. Acta*, 646, 422-32 (1981).
- Clark, M.R., et al., Monovalent cation transport in irreversibly sickled cells. *J. Clin. Invest.*, **62**, 329-37 (1978).
- Clark, M.R., and Shohet, S.B., Hybrid erythrocytes for membrane studies in sickle cell disease. *Blood*, 47, 121-131 (1976).
- Embury S.H., et al., Concurrent sickle cell anemia and α-thalassemia. Effect on pathological properties of sickle erythrocytes. *J. Clin. Invest.*, **73**, 116-23 (1984).
- 9. Friedman, S., and Roll, F.J., Isolation and culture of hepatic lipocytes, Kupffer cells, and sinusoidal endothelial cells by density gradient centrifugation with Stractan. *Anal. Biochem.*, **161**, 207-18 (1987).
- Galili, U., et al., Excessive binding of natural anti-α-galactosyl immunoglobin G to sickle erythrocytes may contribute to extravascular cell destruction. *J. Clin. Invest.*, **77**, 27-33 (1986).
- 11. Lane, P.A., et al., Cellular dehydration and immunoglobulin binding in senescent neonatal erythrocytes. *Pediatr. Res.*, **23**, 288-92 (1988).
- Matovik, L.M., et al., Drug-induced endocytosis of neonatal erythrocytes. *Blood*, 65, 1056-1063 (1985).
- Mohandas, N., et al., A technique to detect reduced mechanical stability of red cell membranes: relevance to elliptocytic disorders. *Blood*, **59**, 768-774 (1982).
- Corash, L.M., et al., High resolution separation of blood cell types by density gradient centrifugation, American Society of Hematology Sixteenth Annual Meeting abstracts. *Blood*, **42**, 1006 (1973).

VNC,MAM 10/07-1

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.