

(Catalog Number: CYL4039)

CATALOG NUMBER: CYL4039 QUANTITY: 200 units (assayed with [³H]-astemizole)

Note: this quantity is equivalent to 400 units

assayed with [125]-BeKm.

LOT NUMBER: RI08030014 VOLUME/CONCENTRATION: 2 mL, 1 mg/mL

BACKGROUND:

The human ether-a-go-go related gene (hERG) is a potassium ion channel which is essential for normal cardiac repolarization. In drug screening models, the hERG K⁺ channel has been indicated to inhibit a wide variety of compounds, and its blockage can lead to cardiac QT interval prolongation and life threatening arrhythmias (Murphy *et al.* 2006). Cardiac safety relating to I_{Kr} K⁺ channels has become a major concern of regulatory agencies, as hERG channel inhibition has been identified as the firmest link to QT prolongation (Chiu *et al.* 2004). Millipore's hERG membrane preparations are crude membrane preparations made from HEK293 stable recombinant cell lines (Millipore cat. # CYL3039), which are ideal HTS tools for screening antagonists against the hERG channel. The membrane preparations exhibit a $K_{\rm d}$ of 5.2 nM for [3 H]-Astemizole. With 10 $\mu \rm g/well$ hERG Membrane Prep and 3.0 nM [3 H]-Astemizole, a greater than 4-fold signal-to-background ratio was obtained.

APPLICATIONS: Radioligand binding assay

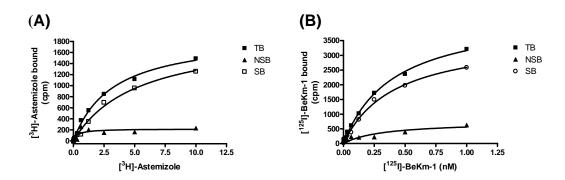


Figure 1. Saturation binding for hERG with [³H]-astemizole and [¹²⁵I]-BeKm. (A), one unit of hERG membrane preparation was incubated with increasing amount of [³H]-astemizole in the absence (total binding, TB) or presence (nonspecific binding, NSB) of 500-fold excess unlabeled astemizole. (B), half a unit of hERG Membrane Preparation was incubated with increasing amount of [¹²⁵I]-BeKm-1 in the absence or presence of 500-fold excess unlabeled BeKm-1. In both curves, specific binding (SB) was determined by subtracting NSB from TR



(Catalog Number: CYL4039)

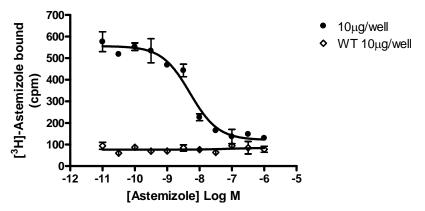


Figure 2. Competition binding for hERG. hERG Membrane Preparation (10 μ g/well) or Wild-Type HEK293 membrane preparation was incubated with 3.0 nM [3 H]-Astemizole and increasing concentrations of unlabeled Astemizole. More than 4- fold signal: background was obtained.

Table 1. Signal: background and specific binding values obtained in a competition binding assay for the hERG membrane prep.

	10μg/well
Signal: background	4.6
Specific binding (cpm)	435

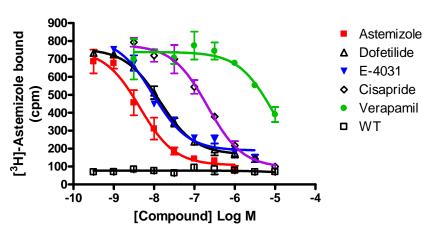


Figure 3. Rank ordering small molecule inhibitors for hERG. hERG Membrane Preparation ($10\mu g$ /well) was characterized by evaluating the activity for known hERG small molecule inhibitors in a competition binding assay. The membranes were incubated with 3.0 nM [3 H]-Astemizole and increasing concentrations of unlabeled compounds to determine sample activity and rank order.



(Catalog Number: CYL4039)

Table 2. Rank order comparison of various small molecule hERG inhibitors using [³H]-astemizole binding assay and E-Phys assays:

	Binding Ki (nM) with hERG membrane preps		Electrophysiology IC50 (nM) with hERG cell lines		
	Millipore CYL4039M value	Literature value*	Millipore CYL4039 by IonWorks PPC	Millipore CYL4039 by PatchExpress	Literature value by Patch Clamp
Astemizole	3.4 ± 0.7	3.3 ± 0.7	4.6 ± 0.003	ND	0.5
Dofetilide	5.8 ± 2.4	28 ± 6	ND	ND	15.3 ± 2.5
E-4031	6.1 ± 0.8	58 ± 5	115 ± 0.02	144 ± 44	14
Pimozide	18.8 ± 2.3	14 ± 4	ND	ND	18
Cisapride	104.9 ± 30.9	123 ± 25	75.6 ± 0.02	5 ± 1	44.5 ± 10.6
Haloperidol	241.4 ± 21.4	234 ± 20	ND	ND	63
Risperidone	> 4000	4302 ± 319	ND	ND	167
Verapimil	> 4000	3902 ± 529	1332 ± 120	3100 ± 400	830

^{*} Data obtained from Chiu et al., 2004

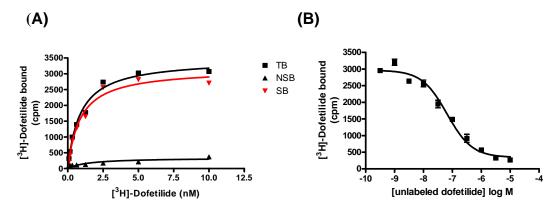


Figure 4. Binding of [³H]-dofetilide to hERG membrane preparation. (A), five units (50 μg) of hERG membrane preparation was incubated with increasing amount of [³H]-Dofetilide (American Radiolabeled Chemicals) in the absence (total binding, TB) or presence (nonspecific binding, NSB) of 1000-fold excess unlabeled Dofetilide. A Kd of 0.9 nM and Bmax of 0.84 pmol/mg were obtained. (B), five units (50 μg) of hERG membrane preparation was incubated with 5 nM [3H]-dofetilide and increasing concentrations of unlabeled dofetilide.

SPECIFICATIONS: 1 unit = $10 \mu g$ membrane preparation with [3 H]-astemizole

B_{max} with [³H]-astemizole: 6.7 pmol/mg

K_d for [³H]-astemizole: 5.2 nM

Species: Human ERG (Accession number U04270)

HOST CELLS: HEK293

RECOMMENDED ASSAY CONDITIONS: Membranes are mixed with radioactive ligand and unlabeled competitor (see Figures 1 and 2 for concentrations tested) in binding buffer in a



(Catalog Number: CYL4039)

nonbinding 96-well plate, and incubated for 1-2 h. Prior to filtration, an FB 96-well harvest plate (Millipore cat. # MAHF B1H) is coated with 0.33% polyethyleneimine for 30 min, then washed with 25 mM Tris, pH 7.4, 130 mM NaCl, 5 mM KCl, 0.8 mM CaCl₂, 0.1% BSA. Binding reaction is transferred to the filter plate, and washed 6 times (250 μ L per well per wash) with Wash Buffer. The plate is dried and counted.

Binding buffer: 10 mM Hepes, pH 7.4, 130 mM NaCl, 5 mM KCl, 0.8 mM MgCl₂, 1 mM NaEGTA, 10 mM glucose, 0.1% BSA, filtered and stored at 4°C

Radioligand: [3H]-Astemizole (Perkin Elmer # NET-1140)

Wash Buffer: 25 mM Tris, pH 7.4, 130 mM NaCl, 5 mM KCl, 0.8 mM CaCl₂, 0.1% BSA, filtered and stored at 4°C.

One package contains enough membranes for at least 200 assays (units), where a unit is the amount of membrane that will yield greater than 4-fold signal: background with ³H-labeled Astemizole.

PRESENTATION:

Liquid in packaging buffer: 50 mM Tris pH 7.4, 10% glycerol and 1% BSA with no

preservatives.

Packaging method: Membrane protein was adjusted to the indicated concentration in packaging buffer, rapidly frozen, and stored at -80°C.

STORAGE/HANDLING:

Maintain frozen at -70°C up to expiration date indicated on the label. Do not freeze and

thaw.

REFERENCES:

Ackerman MJ *et al.* (1998) The long QT syndrome: ion channel diseases of the heart. *Mayo Clin. Proc.* 73: 250-269.

Chiu PJS *et al.* (2004) Validation of a [³H]-Astemizole Binding Assay in HEK293 Cells Expressing hERG K⁺ Channels. *J. Pharmacol. Sci.* 95: 311-319.

Chouabe C *et al.* (1998) HERG and KvLQT1/IsK, the cardiac K⁺ channels involved in long QT syndromes, are targets for calcium channel blockers. *Mol. Pharmacol.* 54: 695-703.

Murphy SM *et al.* (2006) Evaluation of functional and binding assays in cells expressing either recombinant or endogenous hERG channel. *J. Pharmacol. Toxicol. Methods* 54: 42-55

Rampe D *et al.* (1997) A mechanism for the proarrhythmic effects of cisapride (Propulsid): high affinity blockade of the human cardiac potassium channel HERG. *FEBS Lett.* 417: 28-32.

Yao JA *et al.* (2005) Estimation of potency of HERG channel blockers: impact of voltage protocol and temperature. *J. Pharmacol. Toxicol. Methods.* 52: 146-153.

For research use only; not for use as a diagnostic.

Unless otherwise stated in our catalog or other company documentation accompanying the product(s), our products are intended for research use only and are not to be used for any other purpose, which includes but is not limited to, unauthorized commercial uses, in vitro diagnostic uses, ex vivo or in vivo therapeutic uses or any type of consumption or application to humans or animals.

©2008: Millipore Corporation. All rights reserved. No part of these works may be reproduced in any form without permission in writing.