

MultiScreen® Solvinert and MultiScreen Deep Well Solvinert Filter Plates

96-well chemically compatible sample prep platforms

- ▶ Automation-compatible plate designs
- ▶ Low binding and low extractables
- ▶ Membrane offerings accommodate solvents and aqueous samples
- ▶ Designed for extended incubations

High Performance Plate Designs

Automation-compatible MultiScreen Solvinert filter plates are available in 96-well deep well (2 mL) and standard (500 μ L) formats. The one-piece plate designs incorporate rigid sidewalls to facilitate use with robotic gripper arms and provide space for barcode labels.


All MultiScreen Solvinert plate materials are selected for broad chemical compatibility and enable extended in-plate incubations with no leaking.

Optimized for Drug Discovery and HPLC Sample Prep

MultiScreen Solvinert filter plates – both deep well and standard – are optimized for drug discovery applications including total drug analysis, NCE cleavage from solid phase libraries, and sample preparation prior to HPLC or LC/MS/MS. The plates and membranes demonstrate low binding, low extractables and high recoveries. For chromatography separations, the MultiScreen Column Loader accessory can be used to create 96 mini-columns per plate.


Maximum Assay Versatility

MultiScreen Solvinert Filter Plates

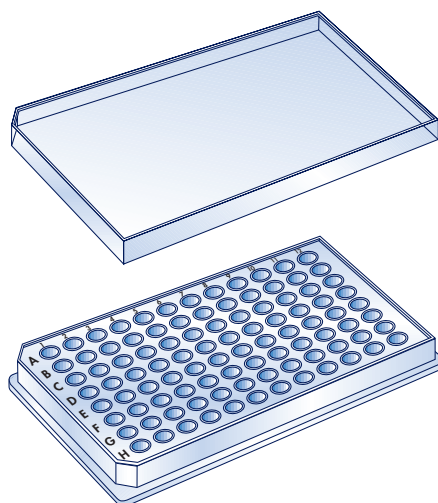

MultiScreen Solvinert Filter Plate holds up to 500 μ L/well.

- Solvent-resistant plate and membrane materials
- Rigid sidewalls designed for robotic gripper arms
- Plastic skirt protects contact with filtration zone and enables easy stacking
- Plates accommodate bar code labels
- One piece design for reliable incubations
- Deep well plate features clear top wells for visual inspection

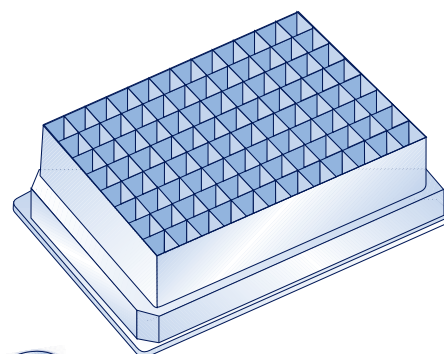

MultiScreen Deep Well Solvinert Filter Plate holds up to 2 mL/well.

MultiScreen Deep Well with hydrophobic membrane and prefilter is also well suited for total drug analysis in plasma by protein precipitation. An optimized protocol is available from Millipore for this application (Millipore Protocol Note AN1218EN00).

Hydrophilic and Hydrophobic PTFE Membranes Available

MultiScreen Solvinert filter plates are available with chemically-resistant hydrophilic and hydrophobic membranes to accommodate aqueous and non-aqueous samples. MultiScreen Deep Well Solvinert filter plates are also available with an optional

pre-filter to use with highly particulate media. Both membranes are highly retentive (>99% retention of acid precipitated BSA).

For the broadest chemical resistance, **hydrophobic PTFE** membrane is recommended. Designed for extended sample incubation times and the lowest extractables, hydrophobic PTFE membrane is ideal for NCE cleavage and cleanup. It is also suitable for in-plate protein precipitation and sample recovery following in-plate compound cleavage from solids or beads. The plates are also used for peptide synthesis.

Hydrophilic PTFE membrane is optimized for low drug and protein binding with excellent throughput in typical aqueous and solvent sample preparation. High sample recoveries and low extractables provide for optimum analysis by HPLC and LC/MS/MS. The membrane is suitable for applications including natural product screening, aqueous solubility testing and total drug analysis.

Filter Selection by Application Needs

MultiScreen Solvinert filter plates are available with hydrophobic and hydrophilic 0.45 µm PTFE membrane. Deep well plates are available with an optional pre-filter with either membrane type. This quick selection table details optimum performance recommendations. For additional selection help, contact your local Millipore office.

Application Need	Hydrophilic 0.45 µm	Hydrophobic 0.45 µm
Highest protein recovery	Y	—
Low drug binding	Y	Y
Low solvent extractables (HPLC/UV)	Y	Y
Aqueous vacuum filtration without prewetting	Y	—
Chemical compatibility	Y	Y
Centrifugal Filtration	Y	Y
Precipitate and particle retention	Y	Y
In-plate protein precipitation with ACN	—	Y
Extended solvent incubations	—	Y
High particulate load*	Y	Y
	(Deep Well with pre-filter)	(Deep Well with pre-filter)

*MultiScreen Deep Well Solvinert filter plates are available with an optional pre-filter.

Performance

Broad Chemical Compatibility

MultiScreen Solvinert and MultiScreen Deep Well Solvinert Filter Plates

Solvent	Filter Plate	Solvent	Filter Plate
Acids		Glycols	
Acetic acid, glacial	R	Glycerine (Glycerol)	R
Hydrochloric acid (37%)	R	Aromatic Hydrocarbons	
Hydrochloric acid (1 N)	R	Benzene	R/TST*
Hydrochloric acid (6 N)	R	Ketones	
Hydrochloric acid (conc.)	R	Acetone	R
Nitric acid (30%)	R	Cyclohexanone	R
Nitric acid (6 N)	R	Oils	
Nitric acid (conc.)	R	Cottonseed oil	R
Sulfuric acid (conc.)	R	Miscellaneous	
Trifluoroacetic acid	R	Acetonitrile (ACN)	R
Alcohols		Dimethylformamide	R
Ethyl Alcohol	R	Dimethyl Sulfoxide (DMSO)	R
Isopropyl Alcohol	R	Dioxane	R/TST*
Methyl Alcohol	R	Formaldehyde (35%)	R
Bases		Hexane	R/TST*
Sodium Hydroxide (70%)	R	Methylene Chloride	L
Sodium Hydroxide (3 N)	R	Pyridine	R
Sodium Hydroxide (conc.)	R		
Ammonium Hydroxide (conc.)	R		

R = Recommended: No significant change was observed in flow rate, nor was there any visible indication of chemical attack.

L = Limited Recommendation: Do not expose for long periods of time. The plate did not show significant physical change after a one-hour incubation.

TST = Testing Recommended

Note: The storage lid does not have the same solvent resistance properties as the filter plate and can be damaged by exposure to some solvent vapors.

Table 1. The data presented in this chart are a compilation of testing by Millipore with certain chemicals and manufacturers' compatibility recommendations. These data are intended to provide expected results when filtration devices are exposed to chemicals under static conditions for 48 hours at 25 °C (77 °F), unless otherwise noted.

Extended Incubation Times With No Leaking

Solution	24-hour Incubation Performance	
	Hydrophilic PTFE Membrane	Hydrophobic PTFE Membrane
Milli-Q 18 MΩ Water	●	●
EtOH, 100%	●	●
MeOH, 100%	●	●
MeCl ₂	○	●
DMSO/PBS (5%/95% v/v)	●	●
ACN, 100%	●	●
ACN/H ₂ O (75%/25% v/v)	●	●
TFA/ACN (80%/20% v/v)	○	●
DMSO, 100%	●	●
DMF, 100%	●	●
NaOH, 1.75N	●	●

● Recommended
○ Not Recommended. Some wells exhibited partial or complete drip out in 24 hour testing. May be compatible for shorter incubation times.

Table 2. MultiScreen Solvinert Filter Plates are designed to support extended incubations. This solvent compatibility table shows results for 200 μL of liquid incubated for 24 hours at room temperature.

Reliable Particle Capture for In-plate Plasma Protein Precipitation and Removal

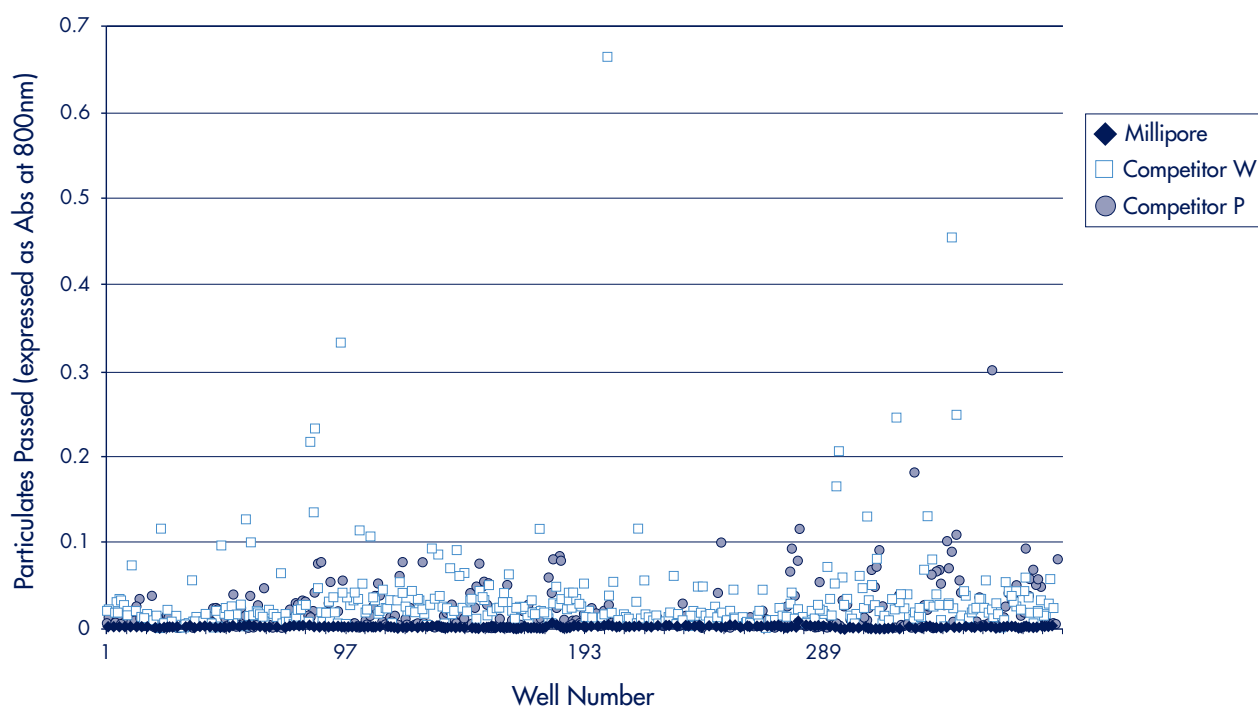


Figure 1. A 1 mL 4:1 mixture of acetonitrile and plasma was incubated by shaking in the plate for 10 minutes followed by vacuum filtration to remove precipitated protein and other insoluble materials. The clarified filtrate was then analyzed for particulates by scattering of visible light at 800 nm. The results are graphed as the $-\log(\%T)$. Millipore has reliable membrane retention for all wells where competitive plates show inconsistent results including passage of measurable amounts of precipitated protein that would harm a HPLC column.

High Recovery of Drug in Solvent-based Preparations

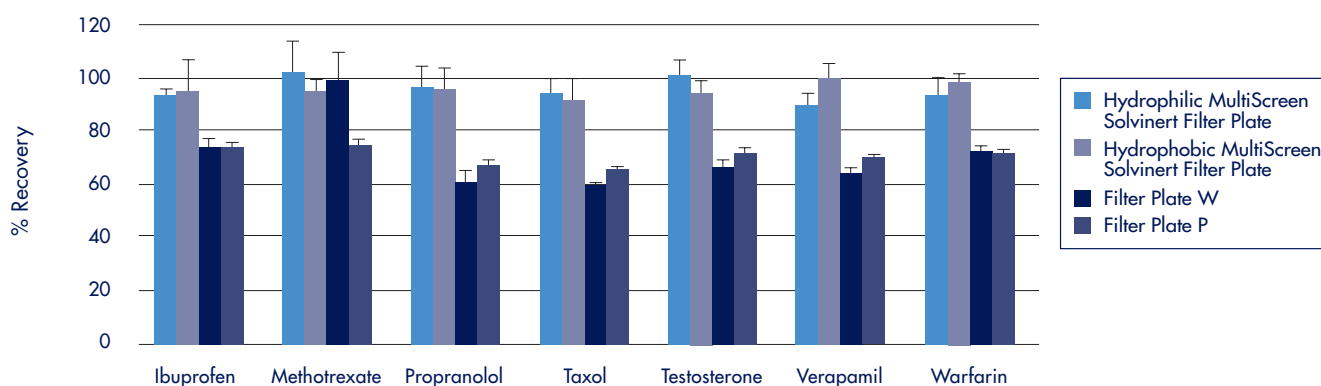


Figure 2. Seven drugs were tested for percent recovery by acetonitrile plasma precipitation. Plasma stock samples (5 mL) were spiked with drug (100 μM stock and tritiated) to a final 5 μM drug concentration followed by 1-hour incubation. Protein was precipitated by the addition of acetonitrile (15 mL) and the solution was vortexed vigorously. For each drug a 300 μL aliquot of the supernatant was added to 8 wells per plate. The samples were filtered by vacuum filtration (12" Hg) and filtrates were collected. Percent recovery was determined by comparing an aliquot (100 μL) of filtrate to an aliquot of the precipitated stock ($\text{cpm filtrate}/\text{cpm precipitated stock}$). Results show superior drug recovery (>90%) for both hydrophilic and hydrophobic MultiScreen Solvinert filter plates. (The sample was pre-precipitated by the addition of ACN followed by vigorous mixing.) Equivalent results are seen with MultiScreen Deep Well Solvinert filter plates.

Specifications

MultiScreen Deep Well Solvinert Filter Plate MultiScreen Solvinert Filter Plate

Materials of Construction

Base plate:	Polyolefin	Polyolefin
Membrane:	PTFE	PTFE
Pre-filter	PP (optional)	N/A

Dimensions

Plate length:	127.8 mm	127.8 mm
Plate width:	85.5 mm	85.5 mm
Plate depth:	40.7 mm	14.6 mm

Sample Volume Per Well

Recommended:	50 µL – 1.8 mL	50 µL – .45 mL
Maximum:	1.9 mL	0.5 mL


Filtration Parameters

Maximum relative centrifugal force (RCF):	3000 x g	3000 x g
Maximum vacuum:	24" Hg	24" Hg


Membrane Area

0.28 cm ²	0.28 cm ²
----------------------	----------------------

Ordering Information

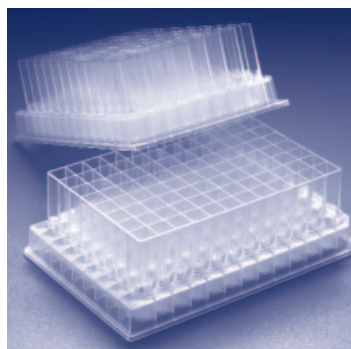

MultiScreen Deep Well Solvinert Filter Plates

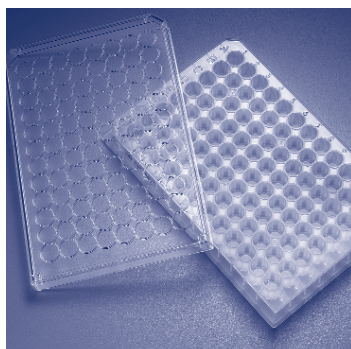

MultiScreen Solvinert Filter Plates

Description	Qty/Pk	Catalogue No.
MultiScreen Deep Well Solvinert Filter Plates (2 mL wells)		
Hydrophobic membrane	5	MDRP NO4 05
	10	MDRP NO4 10
With prefilter	5	MDRP NP4 05
	10	MDRP NP4 10
Hydrophilic membrane	5	MDRL NO4 05
	10	MDRL NO4 10
With prefilter	5	MDRL NP4 05
	10	MDRL NP4 10
MultiScreen Solvinert Filter Plates (500 µL wells)		
Hydrophobic PTFE membrane	10	MSRP NO4 10
	50	MSRP NO4 50
Hydrophilic PTFE membrane	10	MSRL NO4 10
	50	MSRL NO4 50
Accessories		
MultiScreen Vacuum Manifold		MAVM 096 0R
MultiScreen Vacuum Manifold Deep Well Collar		MAVM 096 0T
MultiScreen Vacuum/Pressure Pump	115 Volts, 60 Hz	WP61 115 60
	220 Volts, 50 Hz	WP61 220 50
MultiScreen Column Loader, 25 µL		WP61 220 50
MultiScreen Column Loader, 45 µL		MACL 096 45
MultiScreen Column Loader, 80 µL		MACL 096 80
MultiScreen Column Loader, 100 µL		MACL 096 00
MultiScreen Deep Well Plate Lid (5 pk)		MALI DPP 05
Plate Sealing Tape		MATA HCL 00
Vacuum Flask, 1L		XX10 047 05

To Place an Order or Receive Technical Assistance

For additional information call your nearest Millipore office:

In the U.S. and Canada, call toll-free **1-800-MILLIPORE (1-800-645-5476)**

In the U.S., Canada and Puerto Rico, fax orders to **1-800-MILLIFX (1-800-645-5439)**

Internet: www.millipore.com

Tech Service:

www.millipore.com/techservice

pure e-commerce

Now you can buy Millipore products online @

www.millipore.com/purecommerce



Millipore Worldwide

AUSTRALIA

Tel. 1 800 222 111
or (02) 9888 8999
Fax (02) 9878 0788

AUSTRIA

Tel. 0820 87 44 64
Fax 0820 87 44 44

BALTIC COUNTRIES

Tel. +358 203 05 645
Fax +358 203 05 644

BELGIUM

Tel. 070 225 645
Fax 070 225 644

BRAZIL

Tel. (011) 5548-7011
Fax (011) 5548-7923

CANADA

Tel. (800) 645-5476
Fax (800) 645-5439

CHINA, PEOPLE'S REPUBLIC OF

Beijing:
Tel. (86-10) 8519 1250
(86-10) 6518 1058
Fax (86-10) 8519 1255

Guangzhou:
Tel. (86-20) 8752 0187
(86-20) 8752 0173
Fax (86-20) 8752 0172

Hong Kong:
Tel. (852) 2803 9111
Fax (852) 2513 0313
Shanghai:
Tel. (86-21) 5306 9100
Fax (86-21) 5306 0838

CZECH REPUBLIC

Tel. 02 2051 3841
Fax 02 2051 4298

DENMARK

Tel. 7010 5645
Fax 7010 5644

EASTERN EUROPE, C.I.S., AFRICA, MIDDLE EAST AND GULF

Tel. +33 3 88 38 9536
Fax +33 3 88 38 9539

FINLAND

Tel. 0203 05 645
Fax 0203 05 644

FRANCE

Tel. 0825 045 645
Fax 0825 045 644

GERMANY

Tel. 01 805 045 645
Fax 01 805 045 644

HUNGARY

Tel. 01 209 04 33
Fax 01 209 02 95

INDIA

Tel. 080 839 46 57
Fax 080 839 63 45

IRELAND

Tel. +44 1923 816375
Fax +44 1923 818297

ITALY

Tel. 848 845 645
Fax 848 845 644

JAPAN

Tel. (03) 5442-9711
Fax (03) 5442-9736

KOREA

Tel. (822) 551-0230
Fax (822) 551-0228

LUXEMBOURG

Tel. +32 2726 88 40
Fax +32 2726 98 84

MALAYSIA

Tel. 03-7957-1322
Fax 03-7957-1711

MEXICO

Tel. (55) 5576 9688
Fax (55) 5576 8706

THE NETHERLANDS

Tel. 0900 7 645 645
Fax 0900 7 645 644

NORWAY

Tel. 810 62 645
Fax 810 62 644

POLAND

Tel. 22 669 12 25
Fax 22 663 70 33

PORTUGAL

Tel. +34 91 728 39 60
Fax +34 91 729 29 09

PUERTO RICO

Tel. (787) 273-8495
Fax (787) 747-6553

SINGAPORE

Tel. 6842-1822
Fax 6842-4988

SPAIN

Tel. 901 516 645
Fax 902 011 644

SWEDEN

Tel. 0771 200 645
Fax 0771 200 644

SWITZERLAND

Tel. 0848 645 645
Fax 0848 645 644

TAIWAN

Tel. 886-2-2792-9333
Fax 886-2-2792-6555

U.K.

Tel. 0870 900 46 45
Fax 0870 900 46 44

U.S.A.

Tel. (800) 645-5476
Fax (800) 645-5439

IN ALL OTHER COUNTRIES

Millipore International
Tel. +1 781 533 8622
Fax +1 781 533 8630

MILLIPORE

Millipore and MultiScreen are registered trademark of Millipore Corporation.

Lit. No. PF2813EN00 Printed in U.S.A. 01/04 04-003

Copyright 2004 Millipore Corporation, Billerica, MA 01821 U.S.A. All rights reserved.