

92650 Strep B *ChromoSelect* Selective Supplement

An antibiotic supplement recommended for the selective isolation of Group B Streptococci from clinical samples.

Composition (per vial sufficient for 1000 ml medium)

Colistin	10mg
Nalidixic acid	10mg
Gentamicin	2mg

Directions

Rehydrate the contents of 1 vial aseptically with 5 ml of sterile distilled water. Mix well and aseptically add to 1000 ml of sterile, molten cooled (45-50°C) Strep B *ChromoSelect* Selective Agar Base (Cat. No. 56921). Mix well and pour into sterile Petri plates.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

