

Product Information

IceTEV Protease

TEV protease optimized for low temperatures, Recombinant, expressed in *E. coli*

SAE0195

Product Description

Synonyms: Tobacco Etch Virus protease, TEVp

TEV protease is a highly sequence-specific serine protease from Tobacco Etch Virus.¹ Due to its high specificity, TEV protease is popular for cleavage of recombinant fusion proteins. The optimal sequence for TEV protease cleavage is ENLYFQ\S. However, TEV protease is active on a range of substrates with a consensus sequence of EXLYΦQ\φ where:

- X is any residue
- Φ is any large or medium hydrophobic residue
- φ is any small hydrophobic or polar residue (such as Glycine, Serine, Alanine, Valine, Cysteine)²

Fusion tag removal *in vitro* is the most popular application of TEV protease. One publication has reported that standard TEV protease diminishes in activity by around 10-fold at 4 °C.³

IceTEV protease is specifically designed for cleavage at low temperatures (2-10 °C). After cleavage, this histidine-tagged IceTEV protease can be removed with Ni-agarose beads (such as Cat. No. P6611).

This IceTEV Protease product is supplied at a concentration of ≥ 3 U/ μ L in an aqueous buffer which contains 20 mM Tris-HCl (pH 7.5), 50 mM Sodium Chloride, 1 mM TCEP, 1 mM EDTA and 50% (V/V) glycerol.

Unit Definition

One unit of TEV protease cleaves > 85% of 3 μ g of control substrate in one hour at pH 8.0 at 0-5 °C.

Storage/Stability

The product retains activity for at least 2 years when stored at -20 °C.

Preparation Instructions

TEV Protease is active over a wide range of pH values, ionic strengths, and temperatures. IceTEV protease is an excellent choice for temperature-sensitive proteins. However, the activity toward substrate proteins depends on the substrate identity and the reaction conditions. Since TEV protease is a cysteine protease, the use of low concentration of reducing agent, such as 0.1-2 mM DTT, in the reaction buffer is suggested, to keep the enzyme active during prolonged incubations.

A good starting point for optimization is to use 20 units of TEV Protease per mg of target protein at 2-10 °C for 12-24 hours. Depending on the results, increase or decrease the concentration of TEV protease in subsequent experiments.

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

References

1. Dougherty, W.G. *et al.*, *Virology*, **171**(2), 356-364 (1989).
2. Kapust, R.B. *et al.*, *Biochem. Biophys. Res. Comm.*, **294**(5), 949-955 (2002).
3. Raran-Kurussi, S. *et al.*, *Anal. Biochem.*, **436**(2), 142-144 (2013).

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Standard Warranty

The applicable warranty for the products listed in this publication may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck operates
as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2021 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

SAE0195pis Rev 10/21 DK,DT,SS,GCY

For research use only. Not for use in diagnostic procedures.

The Merck logo, consisting of the word "MERCK" in a bold, red, sans-serif font.