

## Technical Data Sheet

### Colour reference solution Y Certipur®

1.00267.0001

for testing the colour intensity related to Ph Eur.

The preparation of the solutions required is complex and time-consuming. Certipur® reference solutions save time and money.

The Certipur® colour reference solutions are developed for 'Method 1' as described in chapter 2.2.2 of the European Pharmacopeia.



The colour reference solution set 1.00267.0001 contains 7 cuvettes of reference solutions Y1- Y7 and 3 empty sample cells.

#### Benefits:

- Stable solutions with a minimum shelf life of 3 years.
- Ready-to-use solutions
- Certificate of analysis is included in the set.

#### Handling:

To perform the analysis, the sample is filled in the empty cell. This cell is an identical colourness neutral glass grade as the reference tubes. Hold the sample cell alongside the colour comparison solutions in the rack provided. Check the colour in diffused daylight, viewing horizontally against a white background.

#### Ordering Information:

Ordering number: 1.00267.0001

Empty cells with screw caps (25 units) are separately available.

Ordering number: 1.14724.0001

This Certipur® product is produced and analyzed by Merck KGaA, Darmstadt, Germany.