

### SIGMA QUALITY CONTROL TEST

# **ProductInformation**

# Enzymatic Assay of DEOXYRIBONUCLEASE I (EC 3.1.21.1)

#### PRINCIPLE:

**CONDITIONS:** T = 25°C, pH = 5.0,  $A_{260nm}$ , Light path = 1 cm

METHOD: Continuous Spectrophotometric Rate Determination

### **REAGENTS:**

- A. 1 M Sodium Acetate Buffer, pH 5.0 at 25°C
   (Prepare 100 ml in deionized water using Sodium Acetate, Trihydrate, Sigma Prod. No S-8625. Adjust to pH 5.0 at 25°C with 5 M HCl.)
- B. 100 mM Magnesium Sulfate Solution (MgSO<sub>4</sub>)
   (Prepare 5 ml in deionized water using Magnesium Sulfate, Heptahydrate, Sigma Prod. No. M-1880.)
- C. 0.033% (w/v) Deoxyribonucleic Acid Solution (DNA)
   (Dissolve a 1 mg vial of Deoxyribonucleic Acid, Sodium, Sigma Prod. No. D-3664, in 3.1 ml of deionized water.)
- D. 150 mM Sodium Chloride (NaCl)
   (Prepare 50 ml in deionized water using Sodium Chloride, Sigma Prod. No. S-9625.)
- E. Deoxyribonuclease I Enzyme Solution (DNase I) (Immediately before use, prepare a solution containing 400 - 500 Kunitz units/ml in cold Reagent D.)
- F. Standard DNase Solution (Std Soln)<sup>1</sup> (Reconstitute a vial of Deoxyribonuclease I, Standardized vial, containing 2000 Kunitz units, Sigma Prod. No. D-4263, with 1 ml of cold Reagent D. Immediately before use, dilute to 400-500 Kunitz units/ml with cold Reagent D.)

SP-DNA02 Page 1 of 4 Revised: 01/03/97

# Enzymatic Assay of DEOXYRIBONUCLEASE I (EC 3.1.21.1)

### PROCEDURE:

Prepare a reaction cocktail by pipetting (in milliliters) the following reagent into a suitable container:

Reagent A (Buffer)	2.50
Reagent B (MgSO <sub>4</sub> )	1.25
Reagent C (DNA)	3.00
Deionized Water	18.25

Mix by gentle stirring and adjust to pH 5.0 at 25°C with 1 M HCl or 1 M NaOH, if necessary. Then pipette (in milliliters) the following reagents into suitable quartz cuvettes:

	<u>Test</u>	<u>Std</u>	<u>Blank</u>
Reagent D (NaCl)			0.50
Reaction Cocktail	2.50	2.50	2.50

Mix by inversion and equilibrate to  $25^{\circ}$ C. Monitor the A<sub>260nm</sub> until constant, using a suitably thermostatted spectrophotometer. Then add:

Immediately mix by inversion and record the increase in  $A_{260nm}$  for approximately 10 minutes. Obtain the  $\Delta A_{260nm}$ /minute using the maximum linear rate for the Test, Standard, and Blank.

## **CALCULATIONS:**

Units/mI Test Soln = 
$$\frac{(\Delta A_{260nm}/min \text{ Test - } \Delta A_{260nm}/min \text{ Blank}) \text{ (3) (df)}}{(0.001) \text{ (0.5)}}$$
Units/mI Std Soln = 
$$\frac{(\Delta A_{260nm}/min \text{ Std - } \Delta A_{260nm}/min \text{ Blank}) \text{ (3) (df)}}{(0.001) \text{ (0.5)}}$$

3 = Volume (in milliliters) of assay

df = Dilution factor (includes dilution of the standard vial)

 $0.001 = \Delta A_{260nm}$  as per the Unit Definition

0.5 = Volume (in milliliter) of enzyme used

SP-DNA02 Page 2 of 4

Revised: 01/03/97

# Enzymatic Assay of DEOXYRIBONUCLEASE I (EC 3.1.21.1)

CALCULATIONS: (continued)				
CF = -	(2000)			
	Experimental uni	ts/Std vial		
	CF = Correction 2000 = Theoretic			
Corrected Units/ml enzyme = (CF) (Units/ml Test Solution)				
Units/mg solid =	units/ml enzyme			
	mg solid/ml enzyme			
		units/ml enzyme		

## **UNIT DEFINITION:**

One Kunitz unit will produce a  $\Delta A_{260}$  of 0.001 per minute per ml at pH 5.0 at 25°C, using DNA, Types I or III as substrate. [Mg<sup>++</sup>] = 4.2 mM.

## **FINAL ASSAY CONCENTRATION:**

Units/mg protein = -

In a 3.00 ml reaction mix, the final concentrations are 83 mM sodium acetate, 4.2 mM magnesium sulfate, 25 mM sodium chloride, 0.003% (w/v) deoxyribonucleic acid, and 200 - 250 units deoxyribonuclease I.

## **REFERENCE:**

Kunitz, M. (1950) Journal of General Physiology 33, 349-362

mg protein/ml enzyme

Kunitz, M. (1950) Journal of General Physiology 33, 363-377

Lindberg, U. (1964) Biochimica et Biophysica Acta 82, 237-248

SP-DNA02 Revised: 01/03/97

# Enzymatic Assay of DEOXYRIBONUCLEASE I (EC 3.1.21.1)

### **NOTES:**

- 1. There is no absolute standard for the assay of DNase. When the procedure of Kunitz is used, the result is affected by the particular lot of substrate used. For DNase studies, we offer a standard DNase I vial, Sigma Prod. No. D-4263, which has been standardized to contain 2,000 Kunitz units using our DNA, Sigma Prod. No. D-1501 as substrate.
- 2. This assay is based on the cited references.
- Where Sigma Product or Stock numbers are specified, equivalent reagents may be substituted.

Sigma warrants that the above procedure information is currently utilized at Sigma and that Sigma products conform to the information in Sigma publications. Purchaser must determine the suitability of the information and products for its particular use. Upon purchase of Sigma products, see reverse side of invoice or packing slip for additional terms and conditions of sale.

SP-DNA02 Page 4 of 4

Revised: 01/03/97