## Safety Sheet

# Temperature Module

Select language:

|                                    |    |                                   |    |
|------------------------------------|----|-----------------------------------|----|
| English. . . . .                   | 2  | Italiano (Italian) . . . . .      | 30 |
| العربية (Arabic) . . . . .         | 4  | 日本語 (Japanese) . . . . .          | 32 |
| Български (Bulgarian) . . . . .    | 6  | 한국어 (Korean) . . . . .            | 34 |
| 中文 (Chinese, Simplified) . . . . . | 8  | Latviski (Latvian) . . . . .      | 36 |
| Hrvatski (Croatian) . . . . .      | 10 | Lietuvių (Lithuanian) . . . . .   | 38 |
| Čeština (Czech) . . . . .          | 12 | Malti (Maltese) . . . . .         | 40 |
| Dansk (Danish) . . . . .           | 14 | Polski (Polish) . . . . .         | 42 |
| Nederlands (Dutch) . . . . .       | 16 | Português (Portuguese) . . . . .  | 44 |
| Eesti keel (Estonian) . . . . .    | 18 | Română (Romanian) . . . . .       | 46 |
| Suomi (Finnish) . . . . .          | 20 | Srpski (Serbian) . . . . .        | 48 |
| Français (French) . . . . .        | 22 | Slovenčina (Slovak) . . . . .     | 50 |
| Deutsch (German) . . . . .         | 24 | Slovenščina (Slovenian) . . . . . | 52 |
| Ελληνικά (Greek) . . . . .         | 26 | Español (Spanish) . . . . .       | 54 |
| Magyar (Hungarian) . . . . .       | 28 | Svenska (Swedish) . . . . .       | 56 |

## Intended Use

The Opentrons Temperature Module GEN2 is a hot and cold plate module. The module can reach and maintain temperatures ranging from 4 °C to 95 °C. It is often used in protocols that require heating, cooling, temperature changes, or for long duration storage of samples and reagents at a specific, constant temperature. The module is compatible with various thermal block adapters and the Opentrons Flex and OT-2 liquid handling robots.

## Specifications

### Power Requirements

#### Power Adapter

|                       |                            |
|-----------------------|----------------------------|
| Input voltage         | 100-240 VAC (+/- 10%)      |
| Input frequency       | 50/60 Hz                   |
| Input current         | 4 A/115 VAC<br>2 A/230 VAC |
| Output voltage        | 36 VDC                     |
| Output current        | 6.1 A                      |
| Output wattage        | 219.6 W maximum            |
| Power source Required | C13                        |

#### Instrument

|                  |        |                      |
|------------------|--------|----------------------|
| Outer dimensions | Length | 194 mm (7.64 inches) |
|                  | Width  | 90 mm (3.54 inches)  |
|                  | Height | 84 mm (3.31 inches)  |
| Weight           |        | 1.5 kg (3.31 lbs)    |

## Environmental Conditions

For indoor use only. The module is not rated for wet locations or any other outdoor parameter. See Manual for recommended parameters for system operation.

|                     | Acceptable for Operation            | Storage/Transport                   |
|---------------------|-------------------------------------|-------------------------------------|
| Ambient Temperature | 20 °C to 25 °C<br>68 °F to 77 °F    | -10 °C to 60 °C<br>14 °F to 140 °F  |
| Relative Humidity   | 30-80% non-condensing (below 30 °C) | 10-85% non-condensing (below 30 °C) |
| Altitude            | Up to 2000 m above sea level        | Up to 2000 m above sea level        |

IEC61010-1 Pollution Degree II

Overvoltage Category II

Ingress Protection Rating IP2X

## Precautions

Please read these safety instructions and manual carefully before using Flex and/or any module. Keep this Safety Sheet near the instrument for reference.

- This module generates enough heat to cause serious burns. Allow maximum clearance to avoid burns. Do not touch if hot.
- The temperature plate surface may be hot.
- Safety goggles should be worn at all times during operation.
- Do not operate near or with explosive substances. Do not use with substances that may generate explosive gases.
- Read this module's Quick Start Guide and Flex manual before operating.
- Unplug when not in use.

## Electrical Warnings

Failure to follow any of these warnings may result in injury, fire, and electric shock.

- Plug into a grounded, Class 1 circuit.
- Only use the power cord provided.
- Do not use a power cord with inadequate current or voltage ratings. Do not replace the power cord unless at the direction of Technical Service.
- Keep the power cord free of obstructions so you can remove it if necessary.
- Do not connect (plug in), disconnect (unplug), or use AC power cables if:
  - The cable is frayed or damaged.
  - Other attached cables, cords, or receptacles are frayed or damaged.
- Using damaged power cords can cause an electric shock hazard resulting in serious injury or damage to the robot.
- Do not replace the AC power cable unless at the direction of Opentrons Support.
- Do not attempt to repair or remove the outer case of the module or its power supply unless directed by Opentrons Support.

## Maintenance

- Power OFF and disconnect any USB or power cables before beginning cleaning.
- No internal parts require cleaning. The modules should never be disassembled for cleaning.
- Gently wipe the module surfaces with a solution of 70% diluted alchohol (ethyl/ethanol, isopropyl or methanol) and distilled water. Do not use 100% alcohol. Air dry.
- Do not use acetone. This may cause damage or dissolve parts of the instrument.
- Do not autoclave.
- See the manual, Cleaning section for detailed cleaning instruction.

## Technical Service

Do not attempt to modify, revise, or repair the instrument. Only authorized Opentrons Support may service the instrument. Unauthorized repair or service may void the warranty and/or cause hazardous conditions.

## Installation

- Read the manual for detailed installation directions before installing any module.
- Ensure the module is properly oriented in the caddy sized for this module.
- Tighten the caddy anchors.
- Firmly connect all cords to the module, as directed in the manual.
- The mains power cord and plug must be easily accessible in the case of a hazard in which the cord and plug need to be pulled to remove power from the product.

## Equipment Operation

Read the manual before operating.

### Power ON/OFF

Plug into the wall outlet. From the front of the caddy, press the left side of the power switch (-) to turn on. A light (left of the switch) will turn on.

To turn off, press the (o) on the right side of the power switch. The light will turn off. Unplug power cord from outlet when not in use.

### Emergency Disconnect

Push down the Emergency Stop Pendant (E-stop) to quickly stop robot motion. The Flex requires a functional, disengaged E-stop to be attached at all times. The E-stop does not cut power.

In case of an emergency, unplug all power cords from the power source. Contact Opentrons Support before restarting the robot.

## Symbol Definitions

|                                                                                   |                                                                                                                           |
|-----------------------------------------------------------------------------------|---------------------------------------------------------------------------------------------------------------------------|
|  | CAUTION, documentation must be consulted                                                                                  |
|  | CAUTION, possibility of electrical shock                                                                                  |
|  | CAUTION, hot surface                                                                                                      |
|  | CAUTION, risk of pinching                                                                                                 |
|  | CAUTION, risk of explosion!                                                                                               |
|  | Manufacturer/ Legal entity                                                                                                |
|  | This product has been evaluated by an OSHA certified Nationally Recognized Testing laboratory (NRTL) in the United States |
|                                                                                   | Conforms to UL Std 61010-1                                                                                                |
|                                                                                   | Certified to CSA Std 61010-1                                                                                              |
|  | European Council Directives conformity marking                                                                            |
|  | United Kingdom conformity marking                                                                                         |

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© OPENTRONS 2025
Opentrons Labworks, Inc.
Opentrons Flex™ (Opentrons Labworks, Inc.) Registered names, trademarks,