

3050 Spruce Street Saint Louis, Missouri 63103 USA Telephone 800-325-5832 • (314) 771-5765 Fax (314) 286-7828 email: techserv@sial.com sigma-aldrich.com

# ProductInformation

Glucose Oxidase from *Aspergillus niger* 

Product Number **G 9010** Storage Temperature 2-8 °C

# **Product Description**

Enzyme Commission (EC) Number: 1.1.3.4 CAS Number: 9001-37-0 Molecular Weight: 160 kDa (gel filtration)<sup>1</sup> pl: 4.2<sup>2</sup> Extinction coefficient:  $E^{1\%} = 16.7 (280 \text{ nm})^3$ Synonyms: GOD,  $\beta$ -D-Glucose: oxygen 1-oxidoreductase

Glucose oxidase from *Aspergillus niger* is a dimer consisting of 2 equal subunits with a molecular weight of 80 kDa each. Each subunit contains one flavin adenine dinucleotide moiety and one mole of iron. The enzyme is a glycoprotein containing approximately 16% neutral sugar and 2% amino sugars.<sup>1</sup> The enzyme also contains 3 cysteine residues and 8 potential sites for N-linked glycosylation.<sup>4</sup>

Glucose oxidase is capable of oxidizing D-aldohexoses, monodeoxy-D-glucoses, and methyl-D-glucoses at varying rates. D-glucose, 2-deoxy-D-glucose, 4-O-methyl-D-glucoses, 6-deoxy-D-glucose, 4-deoxy-D-glucose, 3-deoxy-D-glucose and 3-O-methyl-D-glucose are oxidized at decreasing rates and in the order listed. The pH optimum for glucose oxidase is 5.5, while it has a broad activity range of pH 4 - 7.<sup>2</sup> Glucose oxidase is specific for  $\beta$ -D-glucose with a K<sub>m</sub> of 33 - 110 mM.<sup>5,6</sup>

Glucose oxidase does not require any activators, but it is inhibited by  $Ag^+$ ,  $Hg_{2+}$ ,  $Cu^{2+}$ , phenylmercuric acetate and p-chloromercuribenzoate. It is not inhibited by the nonmetallic SH reagents: N-ethylmaleimide, iodoacetate, and iodoacetamide.<sup>7</sup>

Glucose oxidase can be utilized in the enzymatic determination of D-glucose in solution. As glucose oxidase oxidizes  $\beta$ -D-glucose to D-gluconolactate and hydrogen peroxide, horseradish peroxidase is often used as the coupling enzyme in glucose determinations. Although glucose oxidase is specific for  $\beta$ -D-glucose, solutions of D-glucose can be quantified as  $\alpha$ -D-glucose will mutorotate to  $\beta$ -D-glucose as the  $\beta$ -D-glucose is consumed by the enzymatic reaction.<sup>8</sup>

# **Precautions and Disclaimer**

For Laboratory Use Only. Not for drug, household or other uses.

### **Preparation Instructions**

This enzyme is soluble in 50 mM sodium acetate buffer, pH 5.1 (0.1 ml in 5 ml), yielding a clear solution.

### References

- Tsuge, H. J., et al, Purification, properties, and molecular features of glucose oxidase from *Aspergillus niger*. J. Biochem., **78**, 835-843 (1975).
- Pazur, J. H., and Kleppe, K., The Oxidation of Glucose and Related Compounds by Glucose Oxidase from *Aspergillus niger*. Biochemistry, 3, 578-583 (1964).
- 3. Fasman, G. D., CRC Handbook of Biochemistry and Molecular Biology, CRC Press (Boca Raton, FL), p. 244 (1990) (Product Code C 3671)
- Frederick, K. R., et al., Glucose oxidase from *Aspergillus niger*. Cloning, gene sequence, secretion from *Saccharomyces cerevisiae* and kinetic analysis of a yeast-derived enzyme. J. Biol. Chem., **265**, 3793-3802 (1990).

- Swoboda, B. E. P., and Massey. V., Purification and properties of glucose oxidase from *Aspergillus niger*. J. Biol. Chem., **240**, 2209-2215 (1965)
- Gibson, Q. H., et al., Kinetics and Mechanism of Action of Glucose Oxidase. J. Biol. Chem., 239, 3927-3934 (1964).
- Nakamura, S., and Ogura, Y., Mode of inhibition of glucose oxidase by metal ions. J. Biochem., 64, 439-447 (1968).
- Bergmeyer, H. U., Methods of Enzymatic Analysis, Weinheim (Deerfield Beach, Fla) 3rd ed., 2, 201-202 (1983).

TMG/SAG 4/03

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.