

3050 Spruce Street Saint Louis, Missouri 63103 USA Telephone (800) 325-5832 (314) 771-5765 Fax (314) 286-7828 email: techserv@sial.com sigma-aldrich.com

ProductInformation

Ammonium sulfate

Product Number A 4418
Store at Room Temperature

Product Description

Molecular Formula: (NH₄)₂SO₄ Molecular Weight: 132.1 CAS Number: 7783-20-2

This product is designated as Molecular Biology grade and is suitable for molecular biology applications. It has been analyzed for the presence of nucleases and proteases.

Ammonium sulfate is a widely used reagent in molecular biology and chromatography. Applications include the precipitation and fractionation of proteins, 1 crystallization of proteins 2,3,4 and of protein-nucleic acid complexes. 5 Ammonium sulfate is also widely used in HPLC of proteins, such as in hydrophobic interaction chromatography. 6,7,8

Ammonium sulfate is also utilized in the purification of antibodies. The use of ammonium sulfate in long PCR buffer, in PCR lysis solution, and in second-strand cDNA synthesis has been reported. A protocol for the HPLC analysis of phospholipids using ammonium sulfate has been reported. 12

Precautions and Disclaimer

For Laboratory Use Only. Not for drug, household or other uses.

Preparation Instructions

This product is soluble in water (435 mg/ml, 3.29 M), yielding a clear, colorless solution. Ammonium sulfate saturation charts at 0 $^{\circ}$ and 25 $^{\circ}$ C have been reported. ¹³

References

- Englard, S., and Seifter, S., Precipitation Techniques. Methods Enzymol., 182, 285-300 (1990).
- McPherson, A., Crystallization of macromolecules: general principles. Methods Enzymol., 114, 112-120 (1985).
- Giegé, R., and Ducruix, A., in Crystallization of Nucleic Acids and Proteins, Ducruix, A., and Giegé, R., eds., Oxford University Press (Oxford, UK: 1999), pp. 1-16.
- Tessier, P. M., et al., Self-interaction chromatography: a novel screening method for rational protein crystallization. Acta Crystallogr. D (Biol. Crystallogr.), 58(Pt 10 Pt 1), 1531-1535 (2002).
- Yaremchuk, A., et al., Improved crystals of Thermus thermophilus prolyl-tRNA synthetase complexed with cognate tRNA obtained by crystallization from precipitate. Acta Crystallogr. D (Biol Crystallogr.), 56(Pt 2), 197-199 (2000).
- Mant, C. T., and Hodges, R. S., in High-Performance Liquid Chromatography of Peptides and Proteins, Mant, C. T., and Hodges, R. S., eds., CRC Press (Boca Raton, FL: 1991), pp. 437-450.
- 7. Hyder, S. M., et al., in High-Performance Liquid Chromatography of Peptides and Proteins, Mant, C. T., and Hodges, R. S., eds., CRC Press (Boca Raton, FL: 1991), pp. 451-475.
- 8. Jacob, L. R., in Protein Liquid Chromatography, Kastner, M., ed., Elsevier (Amsterdam: 2000), pp. 235-269.

- 9. Kent, U. M., Purification of antibodies using ammonium sulfate fractionation or gel filtration. Methods Mol. Biol., **115**, 11-18 (1999).
- 10. Page, M., et al., in Basic Protein and Peptide Protocols, Walker, J. M., ed., Humana Press (Totowa, NJ: 1994), pp. 407-432.
- Molecular Cloning: A Laboratory Manual, 3rd ed., Sambrook, J., and Russell, D. W., CSHL Press (Cold Spring Harbor, NY: 2001), pp. 6.22, 8.78, 11.43-11.47.
- 12. Guan, Z., et al., Separation and quantitation of phospholipids and their ether analogues by high-performance liquid chromatography. Anal. Biochem., **297(2)**, 137-143 (2001).
- 13. Green, A. A., and Hughes, W. L., Protein fractionation on the basis of solubility in aqueous solutions of salts and organic solvents. Methods Enzymol., 1, 67-90 (1955).

GCY/AJH 4/05