

3050 Spruce Street, St. Louis, MO 63103 USA
Tel: (800) 521-8956 (314) 771-5765 Fax: (800) 325-5052 (314) 771-5757
email: techservice@sial.com sigma-aldrich.com

Product Information

Guanosine 5'-[β , γ -imido]triphosphate trisodium salt hydrate

Catalog Number **G0635** Storage Temperature –70 °C

CAS RN 148892-91-5 Synonyms: GMP-PNP, Gpp(NH)p, Guanosine 5'-(β - γ -imido)-triphosphate, β : γ -Imidoguanosine 5'-triphosphate, 5'-Guanylyl-imidodiphosphate

Product Description

Guanosine 5'-[β , γ -imido]triphosphate (GMP-PNP) is a synthetically prepared analog of Guanosine 5'-Triphosphate (GTP), where an NH group rather than an O forms the β - γ bridge. This linkage with the NH group in the triphosphate moiety is non-hydrolyzable. This modification makes GMP-PNP an effective substrate for guanylate cyclase, which cleaves the α - β linkage in the triphosphate group. Guanylate cyclase, which is normally membrane-bound, may then be assayed and characterized in the presence of ubiquitous membrane phosphohydrolases.¹⁻³

GMP-PNP is a competitive inhibitor of the ribosome-dependent GTPase, inhibiting the translocation step in protein synthesis. GMP-PNP has also been used to form stabilized complexes with the α -subunit of the signal receptor particle (SRP), in translocation of proteins through the membrane.

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Preparation Instructions

The solubility of this product is routinely tested at 50 mg/mL in water, yielding a clear, colorless solution. 100 mM stock solutions of GMP-PNP can be prepared in 50 mM Tris-HCl, pH 7.5.7

Storage/Stability

Stock solutions, e.g. 10 mg/mL, should be aliquoted and stored at $-70\,^{\circ}$ C. These should be stable for up to 3 months, with ~5% hydrolysis after 6 months. In addition to being temperature-sensitive, GMP-PNP is very unstable in acidic conditions. Solutions at low pH will rapidly hydrolyze.

Solutions prepared in 20 mM Tris, pH 7.8, and stored at -35 °C with repeated thawing and freezing, have been reported to exhibit <3% hydrolysis after several months.⁴ A separate report indicates that 5 mM stock solutions (pH not identified) can be prepared weekly and stored at -20 °C.⁹ One general guideline is to prepare solutions in 50–100 mM buffer in the range of pH 7.0–8.5, i.e. never having a pH <7.

References

- Yount, R.G., Methods Enzymol., 38, 420-427 (1974).
- 2. Pfeuffer, T., and Helmreich, E.J., *J. Biol. Chem.*, **250(3)**, 867-876 (1975).
- 3. Cassel, D., and Selinger, Z., *Proc. Nat. Acad. Sci.*, **75(9)**, 4155-4159 (1978).
- 4. Eckstein, F. et al., *Biochem. Biophys. Res. Commun.*, **45(5)**, 1151-1158 (1971).
- 5. Data for Biochemical Research, 3rd ed. (Dawson, R.M.C., ed.). Clarendon Press (Oxford, UK), pp. 268-269 (1986).
- 6. Rapiejko, P.J., and Gilmore, R., *J. Cell Biol.*, **119(3)**, 493-503 (1992).
- 7. Seedorf, M. et al., *Mol. Cell. Biol.*, **19(2)**, 1547-1557 (1999).
- 8. Sigma data.
- 9. Cooper, B., and Gregerman, R.I., *J. Clin. Invest.*, **57(1)**, 161-168 (1976),

SBC,GCY,MAM 07/19-1