

3050 Spruce Street, St. Louis, MO 63103 USA
Tel: (800) 521-8956 (314) 771-5765 Fax: (800) 325-5052 (314) 771-5757
email: techservice@sial.com sigma-aldrich.com

Product Information

Trizma® base

Catalog Number **T4661**Store at Room Temperature

CAS RN 77-86-1

Synonyms: Tris base, Tris(hydroxyamino)methane, 2-Amino-2-(hydroxymethyl)-1,3-propanediol, Tris(hydroxymethyl)aminomethane, THAM, Trometamol Note: The "Tris" described in this document is **not** the "Tris" used to flame-proof fabric. That compound, Tris(2,3-dibromopropyl)phosphate, has been reported to be a cancer suspect agent.

Molecular Formula: C₄H₁₁NO₃ Molecular Weight: 121.14

pK_a (20 °C):¹ 8.3

Melting Point: 171-172 °C

Product Description

Trizma[®] is the registered trademark of Sigma-Aldrich applied to various compounds of Tris(hydroxymethyl)-aminomethane that are prepared by Sigma-Aldrich.

This product meets all specifications for Catalog Number T1503, and all analytical and packaging requirements per the current United States Pharmacopoeia (USP). This product is not pyrogen-tested. Certificates of Analysis are available with lot-specific information.

Tris(hydroxyamino)methane, or "Tris" for short, is an established basimetric standard and buffer used in biochemistry and molecular biology.

It may be used by itself as a buffer or as a component of mixed buffer formulations.

These different buffer formulations include:

- Tris-EDTA (TE) buffer
- Tris magnesium buffer
- Tris-acetate-EDTA (TAE) buffer
- Tris-borate-EDTA (TBE) buffer
- Tris-buffered saline (TBS)
- Tris-buffered saline with dextrose (TBS-D)
- Tris-glycine buffer
- Tris-phosphate EDTA buffer
- Tris-SDS buffer
- Tris-sucrose
- Tris-Tricine-SDS buffer

When preparing Trizma solutions at a given pH and temperature, it is necessary to choose the proper mixture of Trizma free base and a corresponding Trizma salt to give the desired final pH at the desired temperature.

Trizma has a significant temperature coefficient, which affects the pH of the solution. For a given concentration the following changes are observed:

- From 5 °C to 25 °C, the pH decreases an average of 0.03 pH units per °C.
- From 25 °C to 37 °C, the pH decreases an average of 0.025 pH units per °C.

This product has been used in various studies and application fields, including:

- Transport properties of mucus³
- Enzymatic assays from cell extracts⁴
- Collagen analysis⁵
- Mitochondria preparation⁶

Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Preparation Instructions

This product is soluble in water (666 mg/mL), yielding a clear, colorless solution.

Storage/Stability

Trizma solutions can be autoclaved.

References

- Gomori, G., Preparation of Buffers for Use in Enzyme Studies. *Methods Enzymol.*, 1, 138-146 (1955).
- Sambrook, J., and Russell, D.W., Molecular Cloning: A Laboratory Manual (3rd ed.). Cold Spring Harbor Laboratory Press (Cold Spring Harbor, NY), pp. 5.8, 5.30, 5.43, 5.60, 5.76, 10.25-10.26, 12.75, 12.84, 12.87, 13.52, 16.29-16.31, A1.2-1.3, A1.7-A1.8, A1.17-A1.18, A1.22, A8.42-A8.43 (2001).
- 3. Odziomek, M. et al., Chem. Process Eng., **38(2)**, 217-229 (2017).

- 4. Çağlayan, M. et al., Bio. Protoc., **5(11)**, e1493 (2015).
- Griffiths, M.M. et al., "Collagen-Induced Arthristis in Rats", in *Methods in Molecular Biology: Arthritis* Research – Methods and Protocols, Volume 2 (A.P. Cope, ed.). Humana Press (Totowa, NJ), Chapter 15, pp. 201-214 (2007).
- 6. Guo, R. et al., Cell, 170(6), 1247-1257 (2017).

Trizma is a registered trademark of Sigma-Aldrich Co., LLC.

GCY,MBK,RXR,MAM 12/18-1