1.18751.0001
1.18751.0007

Spectroquant®

COD Cell Test
for photometers of other manufacturers

COD

1. Definition

The COD (chemical oxygen demand) expresses the amount of oxygen originating from potassium dichromate that reacts with the oxidizable substances contained in 1 l of water under the working conditions of the specified procedure.

1 mol $K_2Cr_2O_7$ is equivalent to 1.5 mol O_2
Results are expressed as mg/l COD (= mg/l O_2)

2. Method

The water sample is oxidized with a hot sulfuric solution of potassium dichromate, with silver sulfate as the catalyst. Chloride is masked with mercury sulfate. The concentration of unconsumed yellow $Cr_2O_7^{2-}$ ions is then determined photometrically.

The method corresponds to DIN ISO 15705 and is analogous to EPA 410.4, APHA 5220 D, and ASTM D1252-06B.

3. Measuring range and number of determinations

Measuring range	Number of determinations
0 - 150.0 mg/l COD	25

For programming data for selected photometers see
www.sigmaaldrich.com/photometry.

4. Applications

This test measures organic and inorganic compounds oxidizable by dichromate.

Exceptions: some heterocyclic compounds (e.g. pyridine), quaternary nitrogen compounds, and readily volatile hydrocarbons.

Sample material:

Groundwater and surface water
In-process controls
Wastewater

5. Influence of foreign substances

The determination is not yet interfered with up to a chloride concentration of 2000 mg/l Cl^- .

The tolerable chloride concentration can be raised to 8000 mg/l Cl^- by placing 0.50 g of mercury(II) sulfate in the reaction cell prior to the sample.

6. Reagents and auxiliaries

Please note the warnings on the packaging materials!

Store the pack protected from light!

The test reagents are stable up to the date stated on the pack when stored closed at +15 to +25 °C.

Package contents:

25 reaction cells
1 sheet of round stickers for numbering the cells

Other reagents and accessories:

MQuant® Chloride Test, Cat. No. 1.10079,
measuring range 500 - ≥ 3000 mg/l Cl^-
Mercury(II) sulfate for analysis, Cat. No. 1.04480
Water for chromatography LiChrosolv®, Cat. No. 1.15333
Spectroquant® CombiCheck 10, Cat. No. 1.14676
COD standard solution CRM, 100 mg/l COD, Cat. No. 1.25029

Pipette for a pipetting volume of 2.0 ml
Thermoreactor

7. Preparation

- Analyze immediately after sampling.
- Homogenize the samples.
- Check the chloride content with the MQuant® Chloride Test. Samples containing more than 2000 mg/l Cl^- must be diluted with distilled water **prior to** determining the COD. The tolerable chloride concentration can be raised by adding mercury(II) sulfate (see section 5).

8. Procedure

	Measurement sample	Blank (only 1x per series)	
Suspend the bottom sediment in the reaction cell by swirling.			
Pretreated sample	2.0 ml	-	Carefully allow to run from the pipette down the inside of the tilted reaction cell onto the reagent (Wear eye protection! The cell becomes hot!).
Distilled water ¹⁾	-	2.0 ml	
Tightly attach the screw cap to the cell. In all subsequent steps the cell must be held only by the screw cap! Vigorously mix the contents of the cell. Heat the cell at 148 °C in the preheated thermoreactor for 120 min. Remove the hot cell from the thermoreactor and allow to cool in a test-tube rack. Do not cool with cold water! Wait 10 min, swirl the cell, and return to the rack for complete cooling to room temperature (cooling time at least 30 min). Measure in the photometer.			

¹⁾ It is recommended to use water for chromatography LiChrosolv®, Cat. No. 1.15333.

Notes on the measurement:

- For photometric measurement the cells must be clean. Wipe, if necessary, with a clean dry cloth.
- Measurement of turbid solutions yields false-low readings.
- The measurement value remains stable over a long term.

9. Analytical quality assurance

recommended before each measurement series
To check the photometric measurement system (test reagent, measurement device, handling) and the mode of working, the COD standard solution CRM (see section 6) or Spectroquant® CombiCheck 10 can be used. Besides a **standard solution** with 80 mg/l COD, CombiCheck 10 also contains an **addition solution** for determining sample-dependent interferences (**matrix effects**).

Additional notes see under **www.qa-test-kits.com**.

For quality and batch certificates for Spectroquant® test kits see the website, where you will find all data in production control, that are determined in accordance with ISO 8466-1 and DIN 38402 A51.

10. Note

The test reagents must not be run off with the wastewater!
Information on disposal can be obtained at
www.disposal-test-kits.com.

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