

3050 Spruce Street, St. Louis, MO 63103 USA
Tel: (800) 521-8956 (314) 771-5765 Fax: (800) 325-5052 (314) 771-5757
email: techservice@sial.com sigma-aldrich.com

Product Information

Beta-1,4-galactosyltransferase 1, human recombinant, expressed in HEK 293 cells

Catalog Number **SAE0093** Storage Temperature –20 °C

EC 2.4.1.38

Synonyms: B4GALT1, Beta-1,4-GalTase 1, b4Gal-T1, UDP-Gal:beta-GlcNAc beta-1,4-galactosyltransferase 1, UDP-galactose:beta-N-acetylglucosamine beta-1,4-galactosyltransferase 1, Beta4Gal-T1

Product Description

 $\beta(1\rightarrow 4)$ galactosyltransferase 1 (B4GALT1) is a type II membrane-bound glycoprotein that transfers galactose from uridine diphosphate- α -D-galactose (UDP-galactose) to acceptor sugars, such as *N*-Acetyl-glucosamine (GlcNAc), in a $\beta(1\rightarrow 4)$ linkage. B4GALT1 resides in the Golgi apparatus of higher eukaryotic cells. A major function of B4GALT1 is the addition of $\beta(1\rightarrow 4)$ linked galactose residues to oligosaccharide acceptors with terminal *N*-acetylglucosamine residues. This is a late elongation step in the *N*-glycan processing pathway.

B4GALT1 enzymatic activity is widely distributed in the vertebrate kingdom, in both mammals and non-mammals, including avians³ and amphibians.⁴ B4GALT1 enzymatic activity has also been demonstrated in a subset of plants which diverged from animals an estimated 1 billion years ago.⁵

B4GALT1 interacts with α -lactalbumin (LA), a protein expressed in the mammary gland during lactation, to form the lactose synthase (LS) complex that transfers galactose from UDP- α -D-Gal to glucose, producing the lactose secreted in milk. Defects in B4GALT1 are the cause of congenital disorder of glycosylation type 2D (CDG2D).

Glomerular B4GALT1 expression has been found to be increased in IgA nephropathy. IgA binding and IgA-induced mesangial cell phosphorylation of spleen tyrosine kinase and IL-6 synthesis were inhibited by a panel of $\beta(1\rightarrow 4)$ galactosyltransferase-specific antibodies, which suggests that IgA binds to the catalytic domain of $\beta(1\rightarrow 4)$ galactosyltransferase.⁷

UniProt: P15291

This product is lyophilized from a 0.22 μ m-filtered solution in 25 mM Trizma® HCl, pH 7.9, and 150 mM NaCl. This recombinant human B4GALT1 is expressed in human HEK 293 cells as a glycoprotein with a calculated molecular mass of 40 kDa. The DTT-reduced protein migrates as a ~55 kDa polypeptide on SDS-PAGE because of glycosylation.

This protein is manufactured in human cells, with no serum. The human cells expression system allows human-like glycosylation and folding, and often supports higher specific activity of the protein.

Purity: ≥95% (SDS-PAGE)

Specific activity: ≥2000 units/mg protein

Unit definition: One unit is defined as the amount of enzyme required to transfer 1.0 nanomole of galactose from UDP-Gal to GlcNAc per minute at pH 7.5, 37 °C.

This recombinant B4GALT1 product can be used to study the mode of action of the enzyme, as well as its potential inhibitors. It can also be used as a glycoengineering tool to modify glycoproteins *in vitro*.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

Storage/Stability

Store the lyophilized product at –20 °C. The product is stable for at least 2 years as supplied.

Preparation Instructions

Briefly centrifuge the vial before opening. Reconstitute in water to a concentration of 0.1 mg/mL. **Do not vortex**. This solution can be stored at 2–8 °C for up to 1 week. For extended storage, it is recommended to store the reconstituted B4GALT1 solution in working aliquots at –20 °C.

References

- Qasba, P.K. et al., Curr. Drug Targets, 9(4), 292-309 (2008).
- 2. Kornfeld, R., and Kornfeld, S., *Annu. Rev. Biochem.*, **54**, 631-664 (1985).
- 3. Shaper, N.L. *et al.*, *J. Biol. Chem.*, **272(50)**, 31389-31399 (1997).
- 4. Lo, N-W. et al., Glycobiology, 8(5), 517-526 (1998).
- 5. Powell, J.T., and Brew, K., *Biochem. J.*, **142(2)**, 203-209 (1974).
- 6. Hanßke, B. et al., *J. Clin. Invest.*, **109(6)**, 725-733 (2002).
- 7. Molyneux, K., et al., Kidney Int., **92(6)**, 1458-1468 (2017).

Trizma is a registered trademark of Sigma-Aldrich Co. LLC.

NA,GCY,MAM 08/18-1