



3050 Spruce Street
Saint Louis, Missouri 63103 USA
Telephone 800-325-5832 • (314) 771-5765
Fax (314) 286-7828
email: techserv@sia1.com
sigma-aldrich.com

Product Information

CAMPYLOBACTER SUPPLEMENT VI (BUTZLER)

Product Number **C3597**

Product Description

An antibiotic supplement recommended by the ISO Committee for the selective isolation of the *Campylobacter* species.

Components

Item

Amphotericin B	1.00 mg
Cefoperazone	7.50 mg
Rifampicin	5.00 mg
Colistin Sulphate	5000 units

Precautions and Disclaimer

For laboratory use only. Not for drug, household or other uses.

Preparation Instructions

Rehydrate the contents of one vial aseptically with 5 mls of 50% (v/v) acetone. Mix well and aseptically add to 470 mls of sterile, molten (C 5595) Columbia Blood Agar Base or (C 5720) Columbia Blood Agar Base with 1% Agar. Add 25 mls of defibrinated sheep blood. Mix well and pour into sterile petri plates.

Storage

Store the supplement at 2-8 °C.

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.