User Guide

Millex[®] Vent With Needle (Sterile)

REF SLFGN25VS

Available languages:

English	3	Latviski (Latvian)	21
Français (French)	5	Lietuviškai (Lithuanian)	23
Italiano (Italian)	7	Magyar (Hungarian)	25
Deutsch (German)	9	Česky (Czech)	27
Español (Spanish)	11	Polski (Polish)	29
Português (Portuguese)	13	Русский (Russian)	31
Dansk (Danish)	15	Norsk (Norwegian)	33
Nederlands (Dutch)	17	Slovenčina (Slovak)	35
Svenska (Swedish)	19	繁體中文 (Traditional Chinese)	37

Definition of Symbols Used

Symbol	Definition	Symbol	Definition
	Sterilized using ethylene oxide		Temperature limit
	Non-pyrogenic		Use-by date
	Consult instructions for use		Date of manufacture
	Do not use if package is damaged and consult instructions for use		Manufacturer
	Do not re-sterilize		Caution
	Do not re-use		CE conformity marking
	Catalogue number		Swiss authorized representative
	Batch code		Unique device identifier

Made in Ireland

Merck Millipore Ltd.
Tullagreen, Carrigtwohill,
Co. Cork, Ireland
T45 KD29

an affiliate of Merck KGaA,
Darmstadt, Germany

MDSS CH GmbH
Laurenzenvorstadt 61
5000 Aarau
Switzerland

Millex® Vent With Needle (Sterile)

SLFGN25VS Qty: 25 pack

- 25 mm
- Single use only
- Sterile
- Non-pyrogenic

Intended Purpose

The Millex® Vent With Needle (Sterile) is a medical device intended for use as a gas vent filter for ventilation of medical devices and pharmacy admixture applications. Device is intended for medical and/or laboratory professional use only.

Introduction

This filter equipped with a needle is a bidirectionally supported sterilizing filter for gases that are not compatible with standard membrane filters. It is used for sterile venting of vials and bottles up to 2 L. This sterile, single-use filter is non-pyrogenic and non-toxic. It consists of a 0.2 µm hydrophobic Fluoropore™ membrane filter sealed in a polyvinyl chloride (PVC) housing. Aqueous solutions are blocked from filtration through the hydrophobic membrane of the device. Clinical benefits include the removal of microorganisms, particles, and undissolved powders larger than the membrane's rated pore size from gaseous solutions to limit cross-contamination.

Chemical Compatibility

This filter is compatible with most gaseous solutions. Compatibility of gaseous solutions should be tested before use. Visit [SigmaAldrich.com/MedMillex](https://www.sigmaaldrich.com/MedMillex) for more information.

WARNINGS

- Filter gaseous solutions with a temperature limit of 45 °C (113 °F).
- To ensure sterility, do not use this product if the package is damaged.
- Use aseptic technique to maintain sterility removing product from primary package.
- Take care to avoid needle injury during handling.
- Do not use the same vent filter to filter gaseous solutions in both directions.
- Sudden loss of pressure could indicate failure of the filter.
- Single use only; do not re-use or resterilize.
- This product is not suitable for filtration of aqueous solutions.
- If liquids enter the unit and fill the housing, the hydrophobic filter blocks the passage of air. Replace with new device if necessary.
- Discard appropriately after single use. See "Disposal" below.

Directions for Use

Use aseptic technique.

1. Spike the Millex® Vent With Needle (Sterile) into the vial of the drug (push the needle only partially through, so its opening is exposed in the vial).
2. Fill a syringe with diluent.
3. Add diluent to the vial.
4. Push the Millex® Vent With Needle (Sterile) into the vial up to the hub of the needle.
5. Hold the syringe, vial, and Millex® Vent With Needle (Sterile), and agitate the vial to reconstitute the drug.
6. Invert the vial and draw the drug into the syringe.
7. Push any air out of the syringe, then remove the syringe from the vial.

Specifications and Performance Characteristics

Materials

Membrane	Hydrophobic Fluoropore™ polytetrafluoroethylene (PTFE), type FG
Pore size	0.2 µm
Housing	Polyvinyl chloride (PVC)
Needle	Stainless steel

Dimensions

Inlet to outlet	61 mm (2.4 in.) including needle
Diameter	28 mm (1.1 in.)
Filtration area	4 cm ² (0.62 in ²)

Temperature limit	45 °C (113 °F)
--------------------------	----------------

Maximum inlet pressure	2 bar (29 psi)
-------------------------------	----------------

Sterilization method	Ethylene oxide gas
-----------------------------	--------------------

Connections	Female Luer-Lok™ inlet, male Luer-slip outlet with needle
--------------------	--

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

Disposal

Follow precautions for disposal of items contaminated with potentially infectious or hazardous biological material according to all applicable international, federal, state, and local regulations.

Technical Assistance

Visit the tech service page on our web site at SigmaAldrich.com/techservice.

Any serious incident of this device should be reported to manufacturer and competent authority of country where user is established.

Standard Warranty

The applicable warranty for the products listed in this publication may be found at SigmaAldrich.com/terms.

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Millipore, Millex, MF-Millipore and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.
© 2006-2023 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

The life science business of Merck operates
as MilliporeSigma in the U.S. and Canada.

PR05875w Rev 06/23