



Product Information

PEPSTATIN A

From Microbial Source

Product Number **P 4265**

Storage Temperature: 2-8 °C

CAS #: 26305-03-3

Product Description

Appearance: White powder

Molecular formula: C₃₄H₆₃N₅O₉

Molecular weight: 685.9

Structure:¹ Isovalery-Val-Val-Sta-Ala-Sta
where Sta = statine = (3S,4S)-4-amino-3-
hydroxy-6-methylheptanoic acid

K_i for Pepsin:² approximately 10⁻¹⁰ M.

Pepstatin A is an inhibitor of acid proteases (aspartyl peptidases). It forms a 1:1 complex with proteases such as pepsin,^{1,2} renin,^{1,2} cathepsin D,^{1,2} bovine chymosin,² and protease B (*Aspergillus niger*).³ The inhibitor is highly selective⁴ and does not inhibit thiol proteases, neutral proteases or serine proteases. Solubilized γ -secretase⁵ and retroviral protease⁶ are also inhibited by Pepstatin A. It has been used to characterize proteases from several sources.^{7,8} Pepstatin A is thought to inhibit by a collected-substrate inhibition mechanism.⁹

This inhibitor is often used as a component in a final mixture with other inhibitors (as in Sigma Protease Inhibitor Cocktails). One recommended set of stock solution concentrations is: bestatin (1.7 mM, selective for aminopeptidase), E-64 (0.22 mM, for cysteine proteases), Pepstatin A (2.5 mM, for aspartyl proteases), AEBSF (18 mM, for serine proteases) and disodium EDTA (86 mM, for metalloproteases).¹⁰

Preparation Instructions

Pepstatin A is only sparingly soluble in water.⁴ It is normally dissolved in a solvent such as ethanol, methanol, or DMSO and diluted into buffer. The stock solution should be of a concentration that the solvent used is diluted at least 1000X in the working solution.

It has been dissolved at 10 mg/mL in ethanol with heat. The resulting solution is colorless, but may appear

hazy. To remove haziness, add up to 50 μ l of glacial acetic acid per mL of ethanol.

At 25 mg/mL DMSO Pepstatin A forms a clear, faint yellow solution.

Stock solutions at 1 mg/mL should be stable at least a week at 4 °C. A 1 mM solution in methanol or DMSO should be stable for months at -20 °C. If solutions become more yellow the reagent is hydrolyzing.

An effective working concentration is 1 μ M, stable for at least one day at room temperature.¹⁰ A typical working concentration is 0.5-1.0 μ g/mL.

Storage/Stability

When stored at 2-8 °C this product has a shelf life of 3 years.

References

1. Umezawa, H., *Methods in Enzymology*, **45**, 689 (1976).
2. Marcinišzyn, J., et al., *Adv. Exp. Med. Biol.* **95**, 199 (1977).
3. Takahashi, K., and Chang, W. J., *J. Biochem. (Tokyo)*, **80**, 497 (1976).
4. Dunn, B.M., in *Proteolytic Enzymes: A Practical Approach*, R.J. Beynon and J.S. Bond, eds. (IRL Press, 1989), p. 63.
5. Li, Y.-M., et al., *Proc. Natl. Acad. Sci USA*, **97**, 6138 (2000).
6. Katoh, I., et al., *Nature*, **329**, 654 (1987).
7. Arima, K., et al., *Phytochemistry*, **54**, 559 (2000).
8. Farias, M. E., and Manca de Nadra, M. C., *FEM Microbiol. Lett.* **185**, 263 (2000).
9. Rich, D. H., et al. *Biochemistry*, **24**, 3165 (1985).
10. *Proteolytic Enzymes: A Practical Approach*, R.J. Beynon and J.S. Bond, eds. (IRL Press, 1989), p. 245 (Appendix III).
11. Umezawa, J., et al., *J. Antibiotics*, **23**, 259 (1970).
12. Morishima, H. et al., *J. Antibiotics*, **25**, 551 (1972).

JWM 3/16/01

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.