

3050 Spruce Street Saint Louis, Missouri 63103 USA Telephone (800) 325-5832 (314) 771-5765 Fax (314) 286-7828 email: techserv@sial.com sigma-aldrich.com

ProductInformation

Mineral oil

Product Number **M8410** Storage at Room Temperature

Product Description

CAS[#] 8020-83-5

This material is tested as an overlay in microdrop embryo cultures. (All embryo tested materials are tested using mouse embryos.) Typically BCF1 x BDF1 hybrids are used. The oil is primarily used to prevent evaporation of the medium, particiularly when performing embryo manipulations.¹ The procedure is as follows: 1-cell (BCF1 X BDF1) embryos were cultured in triplicate microdrops of embryo-tested HTF (Human Tubal Fluid) medium supplemented with 0.4% BSA. For a valid assay, at least 70% of 1-cell stage **control** embryos must develop to the blastocyst stage within 96 hours. For product release, at least 80% of 1-cell stage **test** embryos must develop to the blastocyst stage within 96 hours with the tested material. This product is not sterile. Mineral oil is difficult to sterilize. It can only be sterilized by careful filtration of the warmed oil. It is assumed antibiotics will be added to the cell culture medium. Otherwise, Prod. No. M5310 should be used, as it is sterile filtered.

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

References

 Hogan, B., *et al.*, Manipulating the Mouse Embryo, A Laboratory Manual, 2nd Edition, Cold Spring Harbor Press (Cold Spring Harbor, NY: 1994) pp. 142-143. (Product Number Z358568).

LCM,CMH,SAG,MAM 12/05-1

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.