For life science research only. Not for use in diagnostic procedures.



cOmplete Protease Inhibitor Cocktail Tablets provided in a glass vial

Usi Version: 14

Content Version: November 2020

For the complete inhibition of proteases during extractions from animal and plant tissues or cells, yeast, and bacteria

Cat. No. 11 697 498 001 20 tablets in a glass vial

for 50 ml each

Cat. No. 11 836 145 001 3 x 20 tablets in glass vials

for 50 ml each

Store the product at +2 to +8°C.

1.	General Information	3
1.1.	Contents	3
1.2.	Storage and Stability	3
	Storage Conditions (Product)	
	Storage Conditions (Working Solution)	
1.3.	Additional Equipment and Reagent required	3
1.4.	Application	4
2.	How to Use this Product	5
2.1.	Before you Begin	5
	Safety Information	
	Laboratory procedures	
	Waste handling	
	Working Solution Preparation of 25x-concentrated stock solution	
2.2.	Parameters	
	Absorption	
	Working Concentration	5
3.	Results	6
4.	Additional Information on this Product	7
4.1.	Quality Control	7
5.	Supplementary Information	7
5.1.	Conventions	7
5.2.	Changes to previous version	7
5.3.	Ordering Information	7
5.4.	Trademarks	8
5.5.	License Disclaimer	8
5.6.	Regulatory Disclaimer	8
5.7.	Safety Data Sheet	8
5.8.	Contact and Support	8

1. General Information

1.1. Contents

Vial / Bottle	Label	Function / Description	Catalog Number	Content
1	cOmplete, Protease inhibitor	Each tablet is sufficient for a volume of 50 ml solution.	11 697 498 001	1 glass vial, 20 tablets
	cocktail tablets		11 836 145 001	3 glass vials, 20 tablets each

1.2. Storage and Stability

Storage Conditions (Product)

When stored at +2 to +8°C, the product is stable through the expiry date printed on the label.

Vial / Bottle	Label	Storage
1	cOmplete, Protease inhibitor cocktail tablets	Store at +2 to +8°C. Store dry.

Storage Conditions (Working Solution)

Store stock solution at +2 to +8°C for 1 to 2 weeks or -15 to -25°C for at least 12 weeks.

1.3. Additional Equipment and Reagent required

For preparation of stock solution

Double-distilled water or 100 mM phosphate buffer, pH 7.0

For extraction in the acid pH range

Pepstatin*

For determination of proteolytic activity

Universal Protease Substrate, Casein, resorufin-labeled*

1.4. Application

cOmplete Protease inhibitor cocktail tablets are used for the inhibition of a broad spectrum of serine, cysteine, and metalloproteases, as well as calpains. Due to the optimized composition of the tablets, they show excellent inhibition effects and are well suited for the protection of proteins isolated from:

- Mammalian tissue
- Bacteria
- Yeast
- Plants

cOmplete contains both reversible and irreversible protease inhibitors, therefore add cOmplete to all stock buffers and solutions normally protected with protease inhibitors, and not only during the initial purification steps.

Since cOmplete contains EDTA (18.5 mg/tablet yields a 1 mM solution of EDTA in 50 ml), the extraction buffer should not contain divalent cations such as Ca²⁺, Mg²⁺, or Mn²⁺ otherwise the inhibition of the metalloproteases might be incomplete. If the protein of interest will be purified by IMAC (immobilized metal affinity chromatography), for example, poly-His tagged recombinant proteins, EDTA must be eliminated, for example, by dialysis prior to the chromatography.

Alternatively, the product cOmplete, EDTA-free* can be used. These tablets are identical to cOmplete, with the only difference that no EDTA or other chelating agents are present.

2. How to Use this Product

2.1. Before you Begin

Safety Information

Do not eat the tablets.

Laboratory procedures

- Handle all samples as if potentially infectious, using safe laboratory procedures. As the sensitivity and titer of
 potential pathogens in the sample material varies, the operator must optimize pathogen inactivation by the Lysis /
 Binding Buffer or take appropriate measures, according to local safety regulations.
- Do not eat, drink or smoke in the laboratory work area.
- Do not pipette by mouth.
- Wear protective disposable gloves, laboratory coats and eye protection, when handling samples and kit reagents.
- Wash hands thoroughly after handling samples and reagents.

Waste handling

- Discard unused reagents and waste in accordance with country, federal, state, and local regulations.
- Safety Data Sheets (SDS) are available online on dialog.roche.com, or upon request from the local Roche office.

Working Solution

One cOmplete Protease inhibitor tablet is sufficient for the inhibition of the proteolytic activity in 50 ml extraction solution. If very high proteolytic activity is present, use one tablet for 25 ml extraction buffer. Tablets can be added directly to the extraction medium. Alternatively, a 25x-concentrated stock solution can be prepared.

If it is necessary to inhibit proteolytic activity in a smaller volume, use cOmplete, Mini* (1 tablet for 10 ml extraction solution). The composition of the cOmplete Mini tablet is identical to the normal cOmplete tablet, therefore comparable results are achieved with both products.

Preparation of 25x-concentrated stock solution

Dissolve one cOmplete Protease inhibitor tablet in 2 ml double-distilled water or in 2 ml 100 mM phosphate buffer, pH 7.0.

2.2. Parameters

Absorption

A solution of 1 tablet cOmplete inhibitor dissolved in 50 ml water has an absorption of 0.08 at 280 nm.

Working Concentration

1 tablet per 50 ml extraction solution.

3. Results

cOmplete tablets efficiently inhibit serine, cysteine, and metalloproteases in a broad range.

Extract from	Type of protease present
Animal tissues	Serine, cysteine, and metalloproteases
Plants	Serine and cysteine proteases
Bacterial	Serine and metalloproteases

Occasionally, aspartic proteases (acid proteases) can interfere in isolations from animal tissues. These proteases, however, exhibit pronounced activities only at low pH values. If extraction or single isolation steps must be performed at this pH range, add Pepstatin* to inhibit aspartic protease activity as a precaution.

Typical values for the inhibition of different proteases and protease mixtures by cOmplete are shown in the following table. One cOmplete tablet was added per 50 ml incubation solution. Proteolytic activity was determined with the Universal Protease Substrate, Casein, resorufin-labeled*. When extractions or single-step isolations are necessary in the acid pH range, include Pepstatin* along with cOmplete tablets to ensure aspartic (acid) protease inhibition. All experiments were performed at +15 to $+25^{\circ}$ C.

Research communications indicate that acetylcholinesterase is strongly inhibited by cOmplete. Butyrylcholinesterase is inhibited to a lesser extent.

Protease respective protease mixture	Enzyme concentration [mg/ml]	pH value	% Inhibition after immediate addition to the protease	% Inhibition after 1 hour incubation with the protease
Pancreas extract	0.015	7.8	87	99
	0.03	6.5	88	96
Pronase	0.0015	7.8	88	99
	0.003	7.0	90	95
Thermolysin	0.0008	7.8	99	100
Chymotrypsin	0.0015	7.8	97	97
Trypsin	0.0002	7.8	93	89
Papain	1.0	6.5	95	73

The inhibitory power of cOmplete has been demonstrated with many proteases and protease mixtures. In these experiments, drastically higher concentrations of proteases were used compared to the concentration usually present in extracts, see table above.

4. Additional Information on this Product

4.1. Quality Control

For lot-specific certificates of analysis, see section Contact and Support.

5. Supplementary Information

5.1. Conventions

To make information consistent and easier to read, the following text conventions and symbols are used in this document to highlight important information:

Text convention and symbols			
1 Information Note: Additional information about the current topic or procedure.			
⚠ Important Note: Information critical to the success of the current procedure or use of the product.			
1) 2) 3) etc.	Stages in a process that usually occur in the order listed.		
1 2 3 etc.	Steps in a procedure that must be performed in the order listed.		
* (Asterisk)	The Asterisk denotes a product available from Roche Diagnostics.		

5.2. Changes to previous version

Layout changes.

Editorial changes.

Update to include new safety Information to ensure handling according controlled conditions.

5.3. Ordering Information

Product	Pack Size	Cat. No.		
Reagents, kits				
cOmplete, Mini	25 tablets in a glass vial, for 10 ml each	11 836 153 001		
Universal Protease Substrate	40 mg	11 734 334 001		
Pepstatin	2 mg	10 253 286 001		
	10 mg	11 359 053 001		
	50 mg	11 524 488 001		
cOmplete, EDTA-free	20 tablets in a glass vial, for 50 ml each	11 873 580 001		
	3 x 20 tablets in glass vials, for 50 ml each	05 056 489 001		

5.4. Trademarks

All product names and trademarks are the property of their respective owners.

5.5. License Disclaimer

For patent license limitations for individual products please refer to: **List of biochemical reagent products**.

5.6. Regulatory Disclaimer

For life science research only. Not for use in diagnostic procedures.

5.7. Safety Data Sheet

Please follow the instructions in the Safety Data Sheet (SDS).

5.8. Contact and Support

To ask questions, solve problems, suggest enhancements or report new applications, please visit our **Online Technical Support Site**.

To call, write, fax, or email us, visit **sigma-aldrich.com**, and select your home country. Country-specific contact information will be displayed.

